# OPERATING MANUAL

# Ultra-Low Refrigerated Circulators

air-cooled

FP52-SL

FP55-SL

FP90-SL

water-cooled


FPW52-SL

FPW55-SL

FPW90-SL

FPW91-SL

- upgradable with supplementary pump HSP
- upgradable with supplementary heater HST
- special unit with bath opening for internal temperature applications

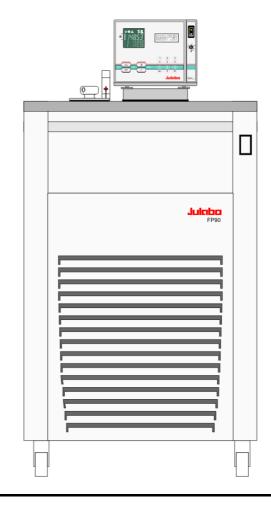



JULABO Labortechnik GmbH 77960 Seelbach / Germany

**+**49 (0) 7823 / 51-0

÷49 (0) 7823 / 24 91

info@julabo.de
www.julabo.de

1.951.2818 03/11

19512818.doc 05.04.11

#### Congratulations!

You have made an excellent choice.

JULABO thanks you for the trust you have placed in us.

This operating manual has been designed to help you gain an understanding of the operation and possible applications of our circulators. For optimal utilization of all functions, we recommend that you thoroughly study this manual prior to beginning operation.

#### The JULABO Quality Management System



Temperature control devices for research and industry are developed, produced, and distributed according to the requirements of ISO 9001:2008. Certificate Registration No. 01 100044846

#### **Unpacking and inspecting**

Unpack the circulator and accessories and inspect them for possible transport damage. Damage should be reported to the responsible carrier, railway, or postal authority, and a damage report should be requested. These instructions must be followed fully for us to guarantee our full support of your claim for protecting against loss from concealed damage. The form required for filing such a claim will be provided by the carrier.

Printed in Germany

Changes without prior notification reserved

**Important:** keep operation manual for future use

# **TABLE OF CONTENTS** Intended use 5 2.5. Setting of temperatures.......34 9.3. Setting the pump pressure ...... 50

9.4.4.Off-Mode	55
9.4.5.ActVar - actuating variable	56
9.4.6.Setting of clock and date	57
9.4.7.Language	57
9.4.8.Unit	
9.4.9.Program type	
9.4.10. Reset	
9.5. Control	
9.5.1. Control internal / external	60
9.5.2.Selftuning	61
9.5.3.Dynamik intern	
9.5.4.CoSpeed - extern	
9.5.5.Control parameters – internal/external	
9.6. Serial Interface	
9.7. ATC Absolute Temperature Calibration, 3-point calibration	65
9.7.1.ATC Fühler - intern / extern	68
9.7.2.ATC Status - yes / no	68
9.7.3.Type 1 -/ 2 -/ 3 point	
9.7.4.Example: 3-point calibration for internal control	
9.8. Limits	
9.8.1.Limits for internal control and external control	72
9.8.2.Limits for external control	73
9.9. Analog inputs/outputs	
9.9.1.Control connector	
9.9.2.Outputs of the connector - REG+E-PROG	79
9.9.3.Input of the connector - REG+E-PROG	
9.9.4.External Stand-by input	
9.9.5.Alarm-output	
10. Troubleshooting guide / Error messages	
11. Electrical connections	
12. Remote control	90
12.1. Setup for remote control	90
12.2. Communication with a PC or a superordinated data system	90
12.3. List of commands	91
12.4. Status messages	
12.5. Error messages	
13. JULABO Service – Online remote diagnosis	
14. Installation of electronic module with analog connectors	
15. Cleaning / repairing the unit	
Anhang: HST Zusatzheizer und HSP Zusatzpumpe	
16. HST booster heater	
16.2. Technical specifications - HST booster heater:	
17. HSP booster pump	
17.1. Pump connection with HSP booster pump	
17.2. Technical specifications - HSP booster pump	

# **Operating manual**

#### 1. Intended use

JULABO circulators have been designed to control the temperature of specific fluids in a bath tank. The units feature pump connections for temperature control of external systems (loop circuit).




JULABO circulators are not suitable for direct temperature control of foods, semiluxury foods and tobacco, or pharmaceutical and medical products. Direct temperature control means unprotected contact of the object with the bath medium (bath fluid).


# 1.1. Description






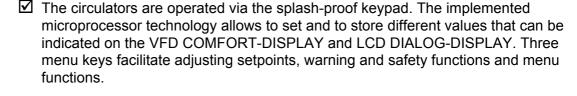




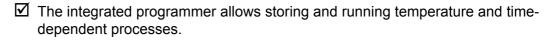



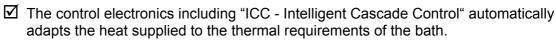



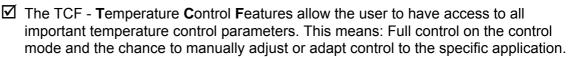



Absolute Temperature Calibration (ATC3) provides high temperature stability in the bath. With the 3-point calibration an offset is adjusted at three temperatures to ensure an accurate temperature pattern at the selected spot in the bath over the full temperature range.



Electrical connections: The serial interface, sw

and set on the LCD DIALOG-DISPLAY.

RS232 RS485 The serial interface, switchable from RS232 to RS485, allows modern process technology without additional interface.

Connection for Pt100 external sensor for external temperature measurement and

Pt100

control. Alarm output for external alarm message or control of JULABO refrigerating baths



or solenoid valve (cooling water). The electronic module (option) provides 3 further analog connections (alarm input, standby input, recorder output, programmer input).



☑ The excess temperature protection conforming to IEC 61010-2-010 is a safety installation independent from the control circuit. This protection can be indicated

SMART **PUMF** 

- ☑ The early warning system for low level signals that bath fluid needs to be refilled before the low level protection conforming to IEC 61010-2-010 causes a complete shut-down of the main functional elements.
- ☑ Intelligent pump system: The pump capacity (electronically adjustable via the motor speed) enables to adapt to varying conditions for internal and external temperature applications.

# 2. Operator responsibility – Safety recommendations

The products of JULABO ensure safe operation when installed, operated, and maintained according to common safety regulations. This section explains the potential dangers that may arise when operating the circulator and also specifies the most important safety precautions to preclude these dangers as far as possible.

- > The operator is responsible for the qualification of the personnel operating the units.
- The personnel operating the units should be regularly instructed about the dangers involved with their job activities as well as measures to avert these dangers.
- Make sure all persons tasked with operating, installing, and maintaining the unit have read and understand the safety information and operating instructions.
- ➤ When using hazardous materials or materials that could become hazardous, the circulator may be operated only by persons who are absolutely familiar with these materials and the circulator. These persons must be fully aware of possible risks.

If you have any questions concerning the operation of your unit or the information in this manual, please contact us!

Contact JULABO Labortechnik GmbH Tel. +49 (0) 7823 / 51-0

Eisenbahnstraße 45 Fax +49 (0) 7823 / 24 91

77960 Seelbach / Germany info@julabo.de www.julabo.de

#### Safety instructions for the operator:

- You have received a product designed for industrial use. Nevertheless, avoid strikes to the housing, vibrations, damage to the operating-element panel (keypad, display), and contamination.
- Make sure the product is checked for proper condition regularly (depending on the conditions of use). Regularly check (at least every 2 years) the proper condition of the mandatory, warning, prohibition and safety labels.
- Make sure that the mains power supply has low impedance to avoid any negative effects on instruments being operated on the same mains.
- This unit is designed for operation in a controlled electromagnetic environment. This means that transmitting devices (e.g., cellular phones) should not be used in the immediate vicinity. Magnetic radiation may affect other devices with components sensitive to magnetic fields (e.g., monitors). We recommend maintaining a minimum distance of 1 m.
- > Permissible ambient temperature: max. 40 °C, min. 5 °C.
- > Permissible relative humidity: 50% (40 °C).
- > Do not store the unit in an aggressive atmosphere.
- > Protect the unit from contamination.
- > Do not expose the unit to sunlight.

#### **Appropriate operation**

Only qualified personnel is authorized to perform configuration, installation, maintenance and repairs of the circulator.

Routine operation can also be carried out by untrained personnel who should however be instructed by trained personnel.

#### Use:

The bath can be filled with flammable materials. Fire hazard!

There might be chemical dangers depending on the bath medium used.

Observe all warnings for the used materials (bath fluids) and the respective instructions (safety data sheets).

Insufficient ventilation may result in the formation of explosive mixtures. Only use the unit in well ventilated areas.

Only use recommended materials (bath fluids). Only use non-acid and non corroding materials.

When using hazardous materials or materials that could become hazardous, **the operator must** affix the enclosed safety labels to the front of the unit so they are highly visible:

When using hazardous materials or materials that could become hazardous, the operator must affix the enclosed safety labels (1 + 2) to the front of the unit so they are highly visible:

Warning label W00: Colors: yellow, black
Danger area. Attention! Observe instructions.
(operating manual, safety data sheet)

Mandatory label M018: Colors: blue, white
Carefully read the user information prior to beginning operation.
Scope: EU

Semi S1-0701 Table A1-2 #9
Carefully read the user information prior to beginning operation.
Scope: USA, NAFTA

Particular care and attention is necessary because of the wide operating range.

There are thermal dangers: Burn, scald, hot steam, hot parts and surfaces that can be touched.



Warning label W26: Colors: yellow, black

Hot surface warning.

(The label is put on by JULABO)

Observe the instructions in the manuals for instruments of a different make that you connect to the circulator, particularly the respective safety recommendations. Also observe the pin assignment of plugs and technical specifications of the products.


# 2.1. Disposal

The circulator contains a back-up battery that supplies voltage to memory chips when the unit is switched off. Do not dispose of the battery with household waste!

Depending on battery regulations in your country, you might be obliged to give back used or defect batteries to gathering places.

The product may be used with oil as bath fluid. These oils fully or partially consist of mineral oil or synthetic oil. For disposal, observe the instructions in the safety data sheets.

This unit contains refrigerants, which at this time are not considered harmful to the ozone layer. However, over the long operating period of the unit, disposal rules may change. Therefore, only qualified personnel should handle the disposal.

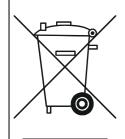

#### Valid in EU countries

Directive 2002/96/EC of the European Parliament and of the Council of 27 January 2003 on waste electrical and electronic equipment (WEEE).

This directive requires electrical and electronic equipment marked with a crossedout trash can to be disposed of separately in an environmentally friendly manner. Contact an authorized waste management company in your country.

Disposal with household waste (unsorted waste) or similar collections of municipal waste is not permitted!

# 2.2. EC Conformity



The products described in the operating instructions conform to the requirements of the following European guidelines:

Directive of the European Parliament and of the Council on the approximation of the laws of the Member States relating to machinery.

EMC guideline with respect to legal harmonization of the member countries concerning electromagnetic compatibility.



# 2.3. Warranty conditions

JULABO Labortechnik GmbH warrants its products against defects in material or in workmanship, when used under appropriate conditions and in accordance with appropriate operating instructions

for a period of ONE YEAR.

Extension of the warranty period – free of charge



With the '1PLUS warranty' the user receives a free of charge extension to the warranty of up to 24 months, limited to a maximum of 10 000 working hours.

To apply for this extended warranty the user must register the unit on the JULABO web site www.julabo.de, indicating the serial no. The extended warranty will apply from the date of JULABO Labortechnik GmbH's original invoice.

JULABO Labortechnik GmbH reserves the right to decide the validity of any warranty claim. In case of faults arising either due to faulty materials or workmanship, parts will be repaired or replaced free of charge, or a new replacement unit will be supplied.

Any other compensation claims are excluded from this guarantee.

# 2.4. Technical specifications

Ultra-Low Refrigerated Circulator		FP52-SL	FPW52-SL
N/ 1: (	0.0	air-cooled	water-cooled
Working temperature range	°C	-60 100	
Special unit with extended	°C	-60 150	
working temperature range			
Temperature stability	°C	±0,05	
Cooling capacity	°C	<u>+20 0 -20</u>	<u>-40 -60</u>
Medium ethanol	kW		0.65 0.1
Cooling compressor		1-stage	
Refrigerant		R404A	
Cooling water flow rate at 20 °C	I/h		ca. 300
Overall dimensions (WxDxH)	cm	59x76x116	
,			
Filling port dia.	cm	7	
Special unit with bath opening	cm	28x23	
(WxL)	OIII	ZONZO	
Bath depth	cm	22	
Filling volume	liters	24	
Weight	kg	156	153
Ambient temperature	°C	5 40	
7 unioniti tomporatare		<b>5</b> III 16	
Cooling water connection			G3/4" male thread
Cooling water connection			US/4 IIIaic IIIIcau
Mains a susan as an att	\//!!	000 440 //00 105 /50 1	000 440 1/00 105 150 1
Mains power connection	V/ Hz	360-440V/3PNPE/50Hz	360-440V/3PNPE/50Hz
400 V/3P/50 Hz Current draw (at 400 V)	٨	19	21
,	A		
Mains power connection 208 V/3P/60 Hz	V/ Hz	190-230V/3PPE/60Hz	190-230V/3PPE/60Hz
	А	25	24
,	•		44
Mains power connection	V/ Hz	207-253V/3PPE/50Hz	
230 V/3P/50 Hz Current draw (at 230 V)	А	25	
Guireili ulaw (at 250 V)	A	25	

Current draw: Refer to notice page 16

Ultra-Low Refrigerated Circulator			FP55-SL air-coole		FPW55-SL water-cooled
Working temperature range	е	°C		-60 100	
Special unit with extended	d	°C		-60 150	
working temperature range	9				
Temperature stability		°C		±0,05	
Cooling capacity		°C		+20 0 -20	<del>-40 -60</del>
Medium ethanol		kW		5.2 4.1 2.2	0.7 0.13
Cooling compressor				2- stage	
Refrigerant				R404A / R23	
Cooling water flow rate at 2	20 °C	l/h			ca. 400
Overall dimensions	(WxDxH)	cm	85x76x1	16	59x76x116
Filling port	dia.	cm		7	
Special unit with bath ope	•	cm		28x23	
Bath depth	(WxL)	cm		23	
Filling volume		liters		27	
			180	21	146
Weight		kg	100	<b>5</b> 40	140
Ambient temperature		°C		5 40	
• " "					
Cooling water connection					G3/4" male thread
Mains power connection		V/ Hz	360-440\	V/3PNPE/50Hz	360-440V/3PNPE/50Hz
	3P/50 Hz		00		00
	at 400 V)	Α	20		20
Mains power connection 208 V/	3P/60 Hz	V/ Hz	190-230\	V/3PPE/60Hz	190-230V/3PPE/60Hz
Current draw (	at 208 V)	Α	28		28

Current draw: Refer to notice page 16

Ultra-Low Refrigerated Circulator		FP90-SL	FPW90-SL
		air-cooled	water-cooled
Working temperature range	°C	-90 100	
Special unit with extended	°C	-90 150	
working temperature range			
Temperature stability	°C	±0,05	
Cooling capacity	°C	<u>+20 0 -20</u>	<del>-40 -60 -80</del>
Medium ethanol	kW	1.8 1.7 1.6	1.35 0.75 0.15
Cooling compressor		2-stage	
Refrigerant		R404A / R23	
Cooling water flow rate at 20 °C	l/h		ca. 250
Overall dimensions (WxDxH)	cm	59x76x116	
, ,			
Filling port dia.	cm	7	
Special unit with bath opening	cm	28x23	
(WxL)			
Bath depth	cm	20	
Filling volume	liters	22	
Weight	kg	195	195
Ambient temperature	°C	5 40	
Cooling water connection			G3/4" male thread
Mains power connection	V/ Hz	360-440V/3PNPE/50Hz	360-440V/3PNPE/50Hz
400 V/3P/50 Hz			
Current draw (at 400 V)	Α	22	19
Mains power connection	V/ Hz	207-253V/3PPE/60Hz	207-253V/3PPE/60Hz
230 V/3P/60 Hz			
Current draw (at 230 V)	Α	29	26
Mains power connection	V/ Hz	207-253V/3PPE/50Hz	207-253V/3PPE/50Hz
230 V/3P/50 Hz			
Current draw (at 230 V)	Α	29	26

Current draw: Refer to notice page 16

Ultra-Low Refrigerated Circulator				
Working temperature range         °C         -91 100           Temperature stability         °C         ±0,2           Cooling capacity         °C         ±20 0 -20 -40 -60 -80 -90 Medium ethanol           KW         5.0 5.2 4.7 4.0 3.5 2.3 0.8 0.2           Cooling compressor         2-stage           Refrigerant         R404A / R508B           Cooling water flow rate at 20 °C         I/h           Ca. 400         -40 0           Overall dimensions         (WXDXH)           Cm         85x76x116           Filling port         dia.           Cm         28x23           Bath depth         cm           Filling volume         liters           Filling volume         liters           Weight         kg           Ambient temperature         °C           Cooling water connection         V/ Hz           400 V/3P/50 Hz           Current draw         (at 400 V)           A 32           Mains power connection         V/ Hz           208 V/3P/60 Hz	Ultra-Low Refrigerated Circulator			· ·
Temperature stability  Cooling capacity  KW  Cooling capacity  KW  Cooling compressor  Refrigerant  Refrigerant  Cooling water flow rate at 20 °C  Coling water flow rate at 20 °C  Cooling water with bath opening (WxL)  Cooling water flow with bath opening (WxL)  Cooling water flow with bath				
Cooling capacity	Working temperature ran	ige	°C	-91 100
Cooling capacity				
Medium ethanol         kW         5.0         5.2         4.7         4.0         3.5         2.3         0.8         0.2           Cooling compressor         2-stage         R404A / R508B         R404A / R508B         R404A / R508B         R400         R4000         R4000         R4000         R4000 <t< td=""><td>Temperature stability</td><td></td><td>°C</td><td>±0,2</td></t<>	Temperature stability		°C	±0,2
Medium ethanol         kW         5.0         5.2         4.7         4.0         3.5         2.3         0.8         0.2           Cooling compressor         2-stage         R404A / R508B         R404A / R508B         R404A / R508B         R400         R4000         R4000         R4000         R4000 <t< td=""><td></td><td></td><td></td><td></td></t<>				
Cooling compressor         2-stage           Refrigerant         R404A / R508B           Cooling water flow rate at 20 °C         I/h         ca. 400           Overall dimensions         (WxDxH)         cm         85x76x116           Filling port         dia.         cm         7           Special unit with bath opening (WxL)         cm         28x23           Bath depth         cm         20           Filling volume         liters         22           Weight         kg         303           Ambient temperature         °C         5 40           Cooling water connection         G3/4" male thread           Mains power connection         V/ Hz         360-440V/3PNPE/50Hz           Current draw         (at 400 V)         A         32           Mains power connection         V/ Hz         207-253V/3PPE/60Hz	Cooling capacity		°C	+20 0 -20 -40 -60 -80 -90
Refrigerant         R404A / R508B           Cooling water flow rate at 20 °C         I/h         ca. 400           Overall dimensions         (WxDxH)         cm         85x76x116           Filling port         dia.         cm         7           Special unit with bath opening (WxL)         cm         28x23           Bath depth         cm         20           Filling volume         liters         22           Weight         kg         303           Ambient temperature         °C         5 40           Cooling water connection         G3/4" male thread           Mains power connection         V/ Hz         360-440V/3PNPE/50Hz           Current draw         (at 400 V)         A         32           Mains power connection         V/ Hz         207-253V/3PPE/60Hz	Medium ethanol		kW	5.0 5.2 4.7 4.0 3.5 2.3 0.8 0.2
Cooling water flow rate at 20 °C         I/h         ca. 400           Overall dimensions         (WxDxH)         cm         85x76x116           Filling port         dia.         cm         7           Special unit with bath opening (WxL)         cm         28x23           Bath depth         cm         20           Filling volume         liters         22           Weight         kg         303           Ambient temperature         °C         5 40           Cooling water connection         G3/4" male thread           Mains power connection         V/ Hz         360-440V/3PNPE/50Hz           Current draw         (at 400 V)         A         32           Mains power connection         V/ Hz         207-253V/3PPE/60Hz	Cooling compressor			2-stage
Overall dimensions (WxDxH) cm 85x76x116  Filling port dia. cm 7  Special unit with bath opening (WxL) cm 28x23  Bath depth cm 20  Filling volume liters 22  Weight kg 303  Ambient temperature °C 5 40  Cooling water connection G3/4" male thread  Mains power connection V/ Hz 360-440V/3PNPE/50Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz  208 V/3P/60 Hz	Refrigerant			R404A / R508B
Overall dimensions (WxDxH) cm 85x76x116  Filling port dia. cm 7  Special unit with bath opening (WxL) cm 28x23  Bath depth cm 20  Filling volume liters 22  Weight kg 303  Ambient temperature °C 5 40  Cooling water connection G3/4" male thread  Mains power connection V/ Hz 360-440V/3PNPE/50Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz  208 V/3P/60 Hz	Cooling water flow rate a	t 20 °C	l/h	ca. 400
Filling port dia. cm 7  Special unit with bath opening (WxL) cm 28x23  Bath depth cm 20  Filling volume liters 22  Weight kg 303  Ambient temperature °C 5 40  Cooling water connection G3/4" male thread  Mains power connection V/ Hz 360-440V/3PNPE/50Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz				
Filling port dia. cm 7  Special unit with bath opening (WxL) cm 28x23  Bath depth cm 20  Filling volume liters 22  Weight kg 303  Ambient temperature °C 5 40  Cooling water connection G3/4" male thread  Mains power connection V/ Hz 360-440V/3PNPE/50Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz	Overall dimensions	(WxDxH)	cm	85x76x116
Special unit with bath opening (WxL) cm 28x23  Bath depth cm 20  Filling volume liters 22  Weight kg 303  Ambient temperature °C 5 40  Cooling water connection G3/4" male thread  Mains power connection V/ Hz 360-440V/3PNPE/50Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz  208 V/3P/60 Hz				
Bath depth cm 20 Filling volume liters 22 Weight kg 303 Ambient temperature °C 5 40  Cooling water connection G3/4" male thread  Mains power connection V/ Hz 360-440V/3PNPE/50Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz	Filling port	dia.	cm	7
Filling volume  Weight  kg 303  Ambient temperature  °C 5 40  Cooling water connection  Mains power connection  V/ Hz 360-440V/3PNPE/50Hz  Current draw  (at 400 V)  Mains power connection  V/ Hz 207-253V/3PPE/60Hz	Special unit with bath opening (WxL)		cm	28x23
Weight kg 303  Ambient temperature °C 5 40  Cooling water connection G3/4" male thread  Mains power connection V/ Hz 360-440V/3PNPE/50Hz  400 V/3P/50 Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz	Bath depth		cm	20
Ambient temperature  °C 5 40  Cooling water connection  G3/4" male thread  Mains power connection  V/ Hz 360-440V/3PNPE/50Hz  400 V/3P/50 Hz  Current draw (at 400 V) A 32  Mains power connection  V/ Hz 207-253V/3PPE/60Hz  208 V/3P/60 Hz	Filling volume		liters	22
Cooling water connection  Mains power connection  V/ Hz  360-440V/3PNPE/50Hz  400 V/3P/50 Hz  Current draw  (at 400 V)  A  32  Mains power connection  V/ Hz  207-253V/3PPE/60Hz  208 V/3P/60 Hz	Weight		kg	303
Cooling water connection  Mains power connection  V/ Hz  360-440V/3PNPE/50Hz  400 V/3P/50 Hz  Current draw  (at 400 V)  A  32  Mains power connection  V/ Hz  207-253V/3PPE/60Hz  208 V/3P/60 Hz	Ambient temperature		°C	5 40
Mains power connection V/ Hz 360-440V/3PNPE/50Hz 400 V/3P/50 Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz 208 V/3P/60 Hz				
Mains power connection V/ Hz 360-440V/3PNPE/50Hz 400 V/3P/50 Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz 208 V/3P/60 Hz	Cooling water connection	١		G3/4" male thread
400 V/3P/50 Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz 208 V/3P/60 Hz	-			
400 V/3P/50 Hz  Current draw (at 400 V) A 32  Mains power connection V/ Hz 207-253V/3PPE/60Hz 208 V/3P/60 Hz	Mains power connection		V/ Hz	360-440V/3PNPE/50Hz
Mains power connection V/ Hz 207-253V/3PPE/60Hz 208 V/3P/60 Hz	•			
208 V/3P/60 Hz	Current draw	(at 400 V)	Α	32
208 V/3P/60 Hz	Mains power connection		V/ Hz	207-253V/3PPE/60Hz
Current draw (at 230 V) A 32		V/3P/60 Hz		
	Current draw	(at 230 V)	Α	32

Current draw: Refer to notice page 16

			SL	
Temperature selection			digital	
·				OC DICDLAY (%C/%E)
via keypad			indication on LCD DIALC	DG-DISPLAY (*C/*F)
remote control via PC			indication on monitor	
Temperature indication			VFD COMFORT-DISPLA	AY (°C/°F)
Resolution		°C	0.01	
ATC3	INT / EXT	°C	±3 / ±9	
Temperature control			ICC - Intelligent Cascade	e Control
Heater wattage (at 230 V	<b>'</b> )		3.0	
Electronically adj. pump of	capacity	stages	1 4	
Flow rate	at 0 bar	l/min	22 26	
Max. pressure	at 0 liters	bar	0.7	
Max. suction	at 0 liters	bar	0.4	
Pump connectors, mal	e thread	mm	M16x1	
Electrical connections:			See page 15	
Ambient temperature		°C	5 40	
Mains power connection	230 V/50-60 Hz	V/ Hz	190 253 / 50-60	
Current draw (	at 208 V / 230 V)	Α	14 / 13	

#### **Electrical connections:**

External alarm device 24-0 V DC / max. 25 mA

Computer interface RS232 or RS485

External Pt100 sensor

Optional for HL, SL

(Order No. 8900100 Electronic module with analog connections)

Programmer input -100 °C to 400 °C = 0 - 10 V or 0 - 20 mA or 4 - 20 mA

Input for the signal of a flow meter or external manipulated variable

Temperature recorder outputs  $0 - 10 \text{ V} (0 \text{ V} = -100 ^{\circ}\text{C}, 10 \text{ V} = 400 ^{\circ}\text{C})$ 

0 - 20 mA (0 mA = -100 °C, 20 mA = 400 °C) 4 - 20 mA (4 mA = -100 °C, 20 mA = 400 °C)

Standby input for external emergency switch-off

Alarm output for external alarm signal

#### Safety installations according to IEC 61010-2-010:

Excess temperature protection adjustable from 0 °C ... 320 °C

Low liquid level protection float switch Classification according to DIN 12876-1 class III

#### Supplementary safety installations

Early warning system for low level float switch

High temperature warning function optical + audible (in intervals)

Low temperature warning function optical + audible (in intervals)

Supervision of working sensor plausibility control

Reciprocal sensor monitoring between

working and safety sensors difference >35 K

Alarm message optical + audible (permanent)
Warning message optical + audible (in intervals)

#### Environmental conditions according to IEC 61 010-1:

Use indoors only.

Altitude up to 2000 m - normal zero.

Ambient temperature: see Technical specifications

Humidity:

Max. relative humidity 80% for temperatures up to +31 °C,

linear decrease down to 50% relative humidity at a temperature of +40 °C

Max. mains voltage fluctuations of ±10% are permissible.

Protection class according to IEC 60 529 IP21

The unit corresponds to Class I

Overvoltage category II Pollution degree 2



#### Caution:

The unit is not suitable for use in explosive environment

EMC requirements according to EN 61326-1

This unit is an ISM device classified in Group 1 (using high frequency for internal purposes), Class A (industrial and commercial range).


# 2.5. Mains power connections



Note concerning current consumption: <u>Mains power connection 3PNPE and 3PPE</u>
The indicated value of current consumption is the maximum useful current of the phase, which also feeds the SL circulator with a 3 kW heater. If the ultra-low refrigerated circulator is operated with the HST additional heater, the 6 kW heating capacity is divided in equal shares (2x3 kW) on the two other phases..

#### Model Mains power connection 3PNPE

FP52-SL	400 V/3PNPE/50 Hz
FPW52-SL	400 V/3PNPE/50 Hz
FP55-SL	400 V/3PNPE/50 Hz
FPW55-SL	400 V/3PNPE/50 Hz
FP90-SL	400 V/3PNPE/50 Hz
FPW90-SL	400 V/3PNPE/50 Hz
FPW91-SL	400 V/3PNPE/50 Hz

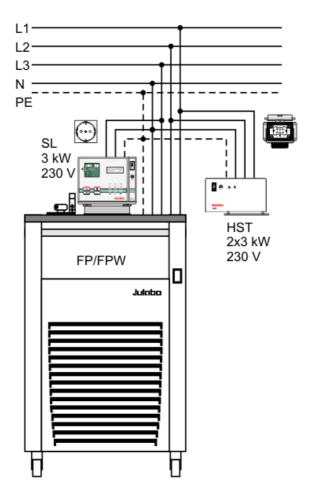


Model	Mains power connection 3PPE
FP52-SL	230 V/3PPE/50 Hz
	208 V/3PPE/60 Hz 208 V/3PPE/60 Hz
	208 V/3PPE/60 Hz 208 V/3PPE/60 Hz
	230 V/3PPE/50 Hz 230 V/3PPE/50 Hz
FP90-SL FPW90-SL	230 V/3PPE/60 Hz 230 V/3PPE/60 Hz

Model Mains power connection 3PPE

FPW91-SL 230 V/3PPE/60 Hz

# Note concerning current consumption:

Each, the SL circulator, the additional HST heater and the cooling unit FPW91 have their own supply mains

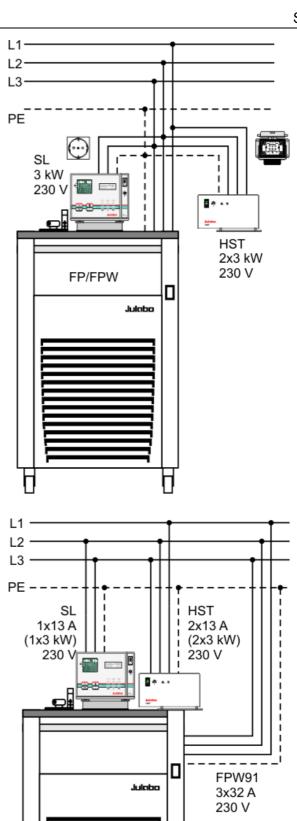

## 2.6. Cooling water connection

Cooling water pressure (IN/OUT) max. 6 bar
Pressure difference (IN - OUT) 3.5 to 6 bar
Cooling water temperature < 20 °C



#### Notice: Cooling water circuit

Risk of oil leaking from the refrigeration system (compressor) of the unit into the cooling water in case of a fault in the cooling water circuit!

Observe the laws and regulations of the water distribution company valid.

Observe the laws and regulations of the water distribution company valid in the location where the unit is operated.



#### Notice:

# Danger of corrosion of heat exchanger due to unsuitable quality of cooling water.

- Due to its high content of lime, hard water is not suitable for cooling and causes scale in the heat exchanger.
- Ferrous water or water containing ferrous particles will cause formation of rust even in heat exchangers made of stainless steel.
- Chlorinated water will cause pitting corrosion in heat exchangers made of stainless steel.
- Due to their corrosive characteristics, distilled water and deionized water are unsuitable and will cause corrosion of the bath.
- Due to its corrosive characteristics, sea water is not suitable.
- Due to its microbiological (bacterial) components, which settle in the heat exchanger, untreated and unpurified river water and water from cooling towers is unsuitable.
- Avoid particulate matter in cooling water.
- Avoid putrid water.

#### Recommended quality of cooling water:

pH	7.5 to 9.0
Sulfate [SO4 2- ]	< 100 ppm
Hydrocarbonate [HCO 3-]/sulfate [SO4 2-]	> 1 ppm
Hardness [Ca 2+, Mg 2+]/[HCO 3-]	> 0.5 °dH
Alkalinity	60 ppm < [HCO 3-] < 300 ppm
Conductivity	< 500 μS/cm
Chloride (Cl -)	< 50 ppm
Phosphate (PO4 3-)	< 2 ppm
Ammonia (NH3)	< 0.5 ppm
Free chlorine	< 0.5 ppm
Trivalent iron ions (Fe 3+)	< 0.5 ppm
Manganese ions (Mn 2+)	< 0.05 ppm
Carbon dioxide (CO2)	< 10 ppm
Hydrogen sulfide (H2S)	< 50 ppm
Content of oxygen	< 0.1 ppm
Algae growth	impermissible
Suspended solids	impermissible

# **Operating instructions**

## 3. Safety notes for the user

## 3.1. Explanation of safety notes



In addition to the safety warnings listed, warnings are posted throughout the operating manual. These warnings are designated by an exclamation mark inside an equilateral triangle. "Warning of a dangerous situation (Attention! Please follow the documentation)."

The danger is classified using a signal word.

Read and follow these important instructions for averting dangers.



#### Warning:

Describes a **possibly** highly dangerous situation. If these instructions are not followed, serious injury and danger to life could result.



#### Caution:

Describes a **possibly** dangerous situation. If this is not avoided, slight or minor injuries could result. A warning of possible property damage may also be contained in the text.



#### Notice:

Describes a **possibly** harmful situation. If this is not avoided, the product or anything in its surroundings can be damaged.

# 3.2. Explanation of other notes



Note!

Draws attention to something special.



Important!

Indicates usage tips and other useful information.

This icon is used in the operating instructions to indicate flashing values or parameters which have to be set or confirmed.

# 3.3. Safety recommendations

Follow the safety instructions to avoid personal injury and property damage. Also, the valid safety instructions for workplaces must be followed.



- Only connect the unit to a power socket with an earthing contact (PE protective earth)!
- The power supply plug serves as a safe disconnecting device from the line and must always be easily accessible.
- Place the unit on an even surface on a base made of nonflammable material.
- Do not stay in the area below the unit.
- Make sure you read and understand all instructions and safety precautions listed in this manual before installing or operating your unit.
- Set the excess temperature safety installation at least 25 °C below the fire point of the bath fluid.
- Observe the limited working temperature range when using plastic bath tanks.
- Never operate the unit without bath fluid in the bath.
- Pay attention to the thermal expansion of bath oil during heating to avoid overflowing of the fluid.
- Prevent water from entering the hot bath oil.
- Do not drain the bath fluid while it is hot! Check the temperature of the bath fluid prior to draining (e.g., by switching the unit on for a short moment).
- Use suitable connecting tubing.
- Avoid sharp bends in the tubing, and maintain a sufficient distance from surrounding walls.
- Make sure that the tubing is securely attached.
- Regularly check the tubing for material defects (e.g., for cracks).
- Never operate damaged or leaking units.
- Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures, or before moving the unit
- Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit.
- Always empty the bath before moving the unit.
- Transport the unit with care.
- Sudden jolts or drops may cause damage in the interior of the unit.
- Observe all warning labels.
- Never remove warning labels.
- Never operate units with damaged mains power cables.
- Repairs are to be carried out only by qualified service personnel.



Some parts of the bath tank and the pump connections may become extremely hot during continuous operation. Therefore, exercise particular caution when touching these parts.



#### Caution:

The temperature controlling i.e. of fluids in a reactor constitutes normal circulator practice.

We do not know which substances are contained within these vessels. Many substances are:

- inflammable, easily ignited or explosive
- hazardous to health
- · environmentally unsafe

i.e.: dangerous

# The user alone is responsible for the handling of these substances! The following questions shall help to recognize possible dangers and to reduce the risks to a minimum.


- Are all tubes and electrical cables connected and installed?
   Note:
  - sharp edges, hot surfaces in operation, moving machine parts, etc.
- Do dangerous steams or gases arise when heating?
   Is an exhaust needed when working?
- What to do when a dangerous substance was spilled on or in the unit?
   Before starting to work, obtain information concerning the substance and determine the method of decontamination.



## Notice: Check the safety installations at least twice a year!

- Excess temperature protection according to IEC 61010-2-010.
   With a screwdriver turn back the adjustable excess temperature protection until the shut-down point (actual temperature).
- Low level protection according to IEC 61010-2-010.
   To check the function of the float, it can be manually lowered with a screwdriver for example.

# 4. Operating controls and functional elements

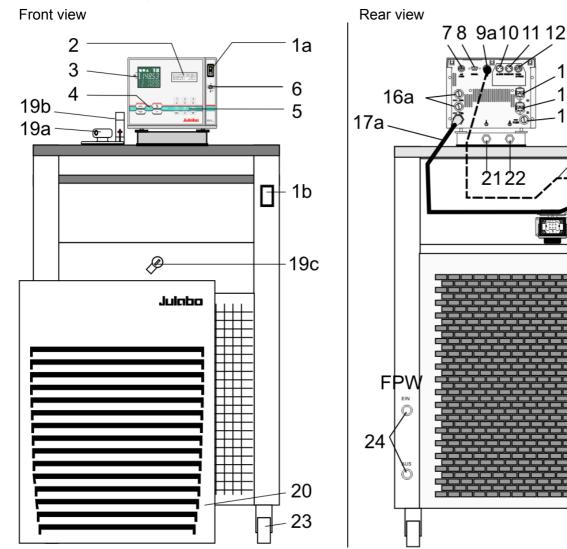

1a

1b



Mains power switch, illuminated for circulator

Mains power switch, illuminated for cooling machine

2



#### **VFD COMFORT-DISPLAY**

Header: Control indicators

Line 1: Actual value internal or external

The display is depending on the selected control mode in the

15

9 9b

18

25

17b

menu > Control < (internal or external).

Line 2: Working temp. setpoint, constantly S xxx.xx

Line 3: Actual value (E = external or I = internal)

Alternating with the display in line 1

Use the keys to indicate further values in line 3

PI Capacity in % - with manipulated variable set to >control<\*

PS Capacity in % - with manipulated variable set to >serial<\* or

#### >eprog<\*

- H Heater capacity in Watts
- U Mains voltage Volts
- F Flow rate in liters/minute (providing EPROG input set to >Flow rate<)

\*refer to >MENU / CONFIGIGURATIO> page 41

2.1	Control indicators in the header:
<u>₩</u> ※ ∧	Heating / Cooling / Alarm /
R	Remote control
2.2	Control indicators in the header:
°C Int	Temperature indication Internal or External actual value
°F Ext	Temperature indication in °C or °F
2.3 🌊 🗒	Display for the adjusted pump pressure stage in the -OFF- mode. Display for the effective pump pressure stage (rotation speed) after start. Four stages, adjustable via the MENU button, in the menu >PUMP<.
3	LCD DIALOG-DISPLAY
Setpoint1: 150.00°C  IntIst : 148.53°C	Line 1: Setpoint and origin of setpoint programming
Power : 80 % Control : intern	(Key / RS232 or RS485 / ext. Pt100 / EProg)
RS232 : 150.00°C	Line 2: Actual value - internal or external,
ExtAct : 148.53°C	identical to line 1 of the VFD-COMFORT-DISPLAY

# ! SAFETY VALUES ! \*SafeTemp : 100.00°C Type Warnin9 \*OverTemp 87.00°C SubTemp 83.00°C SafeSens : 25.00°C

80 %

extern

Power

Control

#### Navigation aid in MENU - Window

Keys / :


Line 3: Heating capacity in %.

Line 4: Control type: internal / external

Selection of menu items / parameters

- Setting in line 3:

Actual value / parameter

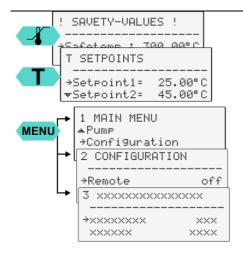

#### Orientation aid in MENU - Window

Line 1 - name with allocation to key



Safety values

T - Setpoint

MENU

1 Main menu Level 1
2 Submenu Level 2
Example: CONFIGURATION

3 Submenu

Level 3

4	Keypad 1	Navigation keys
4.1	ок	1. Key: >OK< Start / Stop (pump / heater )
		2. >OK< in the menu Menu item / select submenu for setting
		Save set value
		Save selected parameter
	~	A beep signals the end of setting
		After the actions Start, Stop and change from VFD Display to standard
		display the key <b>OK</b> is locked for a short time. The above graph "front side" shows an example for standard display.
4.2	4	1. Key: >Return< Stop (pump / heater)
4.2		Special display on VFD Display leave line 3.
		2. >Return< in the menu one menu level down
		Correction function for parameters or values (prior to OK)
		<b>4 (1)</b>
		immediately back to standard display
		οκ 🕑 - Ć 🕑 icon for "keep key pressed down".
4.3		Key: >Up / Down <temperature decrease="" increase="" setpoint<="" td="" –=""></temperature>
		Push key quickly for single steps,
		Keep key pressed for fast change.
		2. >Up/Down< in the menu selection of menu items / parameters
		Menu keys
4.4		Key: start the menu > warning and safety values<
4.5		Key: start the menu >temperature setpoints<
4.6	MENU	Key: display of MENU structure
5	Keypad 2	
	0 9	Numeric keypad: numerals 0 to 9
	⟨ → ⟩	minus / decimal point
	⟨←⟩	Backspace key Setpoint1: 155 °C
		on LCD DIALOG-DISPLAY.
6	160 80 40 40 320 320	Adjustable excess temperature protection according to IEC 61010-2-010

7	ext Pt100	Socket for external measurement and control sensor or external setpoint programming
8	0 0	Interface RS232 / RS485: remote control via personal computer
	SERIAL	
9		Control cable of JULABO refrigerated circulator
9a	*	Socket: output for alarm messages
9b		Socket:
	<b>₩</b>	


Option	ո։ Electronic modւ	ıle Order No. 8 900 100
		The circulator automatically recognizes the connected electronic module.
10		Alarm output (for external alarm signal)
	ALARM	
11		Standby input (for external emergency switch-off)
	STAND-BY	
12		Programmer input and temperature recorder output
	REG+E-PROG	

13	<b>△</b> *	Connector for supplementary pump or solenoid valve 230 V / max. 1.25 A			
	_    <u></u>	No control voltage in the -OFF- condition			
14	M M	Connector solenoid valve			
	<u> </u>	<ol> <li>No control voltage in the -OFF- condition</li> </ol>			
		2. Configurable (Refer to page 74)			
15		Fuses (for connectors 13, 14), T1.25A			
16		Mains fuses, T16A			
17a		Mains power cable with plug for circulator			
17b		Mains power cable with plug cooling machine			
18	(··)	Built-in mains outlet for connection of circulator (230 V / 16 A)			
		Not on - FPW91-SL - 230 V/3PPE/60 Hz			
19	$\bigcirc \ominus \bigcirc \bigcirc \bigcirc$	Drain screw with drain connection			
20		Venting grid, removable			
21					
22		Pump connections Usuction pump Upressure pump			
23		Castor with brake (at the front)			
		Castor without brake (at the back)			
24	OUT IN				
	$\Lambda$	Kühlwasseranschlüsse für FW Kältethermostate			
	<u> </u>				
25		Built-in mains outlet for connection of HST booster heater			
		Not on - FPW91-SL - 230 V/3PPE/60 Hz			
		and Special units with bath opening			


# 5. Preparations

#### 5.1. Installation

Top view:

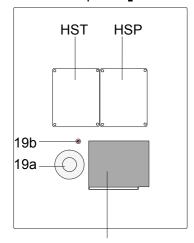

Special unit

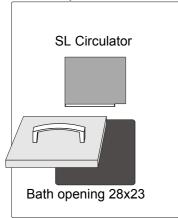





 Place the unit on an even surface on a pad made of non-flammable material. Using the castors move the unit to the intended location.
 For better stability, apply the holding breaks on the front casters.

Upgradable with supplementary pump - HSP see page 102 Upgradable with supplementary heater - HST see page 100

- Cooling machine, pump motor and electronics produce intrinsic heat that is dissipated via the venting openings! Never cover these openings!
- Keep at least 20 cm of open space on the front and rear venting grids.
- The place of installation should be large enough and provide sufficient air ventilation to ensure the room does not warm up excessively because of the heat the instrument rejects to the environment. (Max. permissible ambient temperature: 40 °C).

For a fault (leakage) in the refrigeration system, the standard EN 378 prescribes a certain room space to be available for each kg of refrigerant.

The refrigerant quantity is specified on the type plate.

- > For 0.48 kg of refrigerant R404A, 1 m<sup>3</sup> of space is required.
- > For 0.68 kg of refrigerant R23, 1 m<sup>3</sup> of space is required.
- Do not set up the unit in the immediate vicinity of heat sources and do not expose to sun light
- Before operating the unit after transport, <u>wait about one hour after setting it up.</u> This will allow any oil that has accumulated laterally during transport to flow back down thus ensuring maximum cooling performance of the compressor.

#### Only for water-cooled models:

Ensure circulation of cooling water by connecting the tubing to the cooling water inlet and outlet at the rear of the refrigerated circulator. Cooling water connections: G 3/4"

Important: For cooling water, see page 18

#### 5.2. Bath fluids



#### Caution:

Carefully read the safety data sheet of the bath fluid used, particularly with regard to the fire point!

If a bath fluid with a fire point of ≤65 °C is used, only supervised operation is possible.

Water: The quality of water depends on local conditions.

- Due to the high concentration of lime, hard water is not suitable for temperature control because it leads to calcification in the bath.
- Ferrous water can cause corrosion even on stainless steel.
- Chloric water can cause pitting corrosion.
- Distilled and deionized water is unsuitable. Their special properties cause corrosion in the bath, even in stainless steel.

#### Recommended bath fluids:

Bath fluid	Temperature range
soft/decalcified water	5 °C to 80 °C
mixture water/glycol, mixture 1:1	-20°C to 50°C

#### JULABO bath fluids

JULABO		Thermal	Thermal	Thermal
Description		G	HY	H5
Order Number	10 liters	8 940 124	8 940 104	8 940 106
	5 liters	8 940 125	8 940 105	8 940 107
Temperature range	°C	-30 80	-80 55	-50105
Flash point	°C		78	124
Fire point	°C		80	142
Color		light yellow	clear	clear

JULABO		Thermal	Thermal
Description		H10	H20S
Order Number 10 liters		8 940 114	8 940 108
	5 liters	8 940 115	8 940 109
Temperature range	°C	-20 180	0 220
Flash point	°C	190	230
Fire point °C		216	274
Color		clear	light brown



See website for list of recommended bath fluids.

ATTENTION:

The maximum permissible viscosity is 70 mm<sup>2</sup>/s·



#### Caution:

Fire or other dangers when using bath fluids that are not recommended: Use only non-acidic and non-corroding bath fluids.

JULABO assumes no liability for damage caused by the selection of an unsuitable bath liquid.

Unsuitable bath fluids are fluids which, e.g.,

- are highly viscous (much higher than recommended at the respective working temperature)
- have a low viscosity and have creep characteristics
- have corrosive characteristics or
- tend to crack.
- . No liability for use of other bath fluids!

# 5.3. Temperature application to external systems




Caution:

Securely attach all tubing to prevent slipping.

If the circulator is operated without external system, close the pump connector (22) with the cap nut.


#### Temperature application to external, closed systems

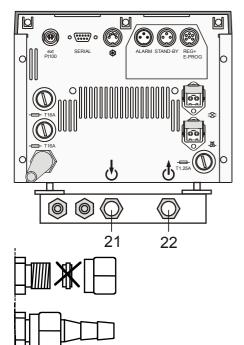

The circulator is used for temperature application to external, closed systems (loop circuit) with simultaneous temperature application in the circulator bath.

# Model combination without HSP booster pump Pump connection:

- Unscrew the M16x1 collar nuts on the pump connectors with a 19 mm (3/4") wrench and remove the sealing disks. Using the collar nuts, screw on the tubing connection fittings (for tubing 8 mm or 12 mm in diameter) delivered with the unit and tighten firmly. (Pressure pump: 22, suction pump: 21)
- Push on the tubing, and secure with tube clamps.
- Attach the tubing to the connectors of the external closed system, e.g., an instrument with a pressure-resistant temperature jacket or a temperature coil, and fasten with tube clamps to prevent slipping.

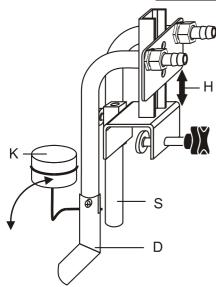

S = Suction pump connection

D = Pressure pump connection

K = Float

H = Height adjustment

The circulator is equipped with both a pressure and suction pump for external temperature application in open systems.


Differing flow rates of the pressure and suction pumps should be compensated. To maintain a constant liquid level, the JULABO "D+S" Level Adapter is recommended for the external bath tank. The flow rate of the pressure pump will be then regulated by a built-in float device. The liquid level may be changed by a height adjustment on the "D+S" Level Adapter.

Accessory: "D+S" Level Adapter Order No. 8 970 410

#### Important:

- The liquid level should be equal in the internal and external baths (absolute height).
- i) If you take out samples (for example Erlenmeyer flasks) from the external bath, turn the circulator off with the Start/Stop key.

#### **Backflow safety device**

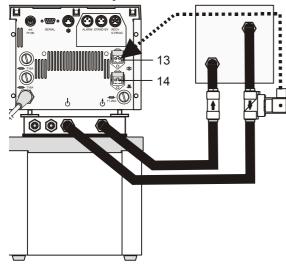

2 INPUTS/OUTPUTS

----
SV-OUT Backflow

\*Chan1 | ActInt

(see page 74)

If the liquid levels in the circulator bath and the external system are at different heights, overflowing must be prevented after the power has been turned off.



#### Flood hazard!

For this reason, solenoid valves for loop circuit or shut-off valves can be integrated in the loop circuit.

- ① Connect the valve to the connector (13).
- (i) If socket (14) is used: In menu >Inputs/Outputs< set the menu item >SV-Out< to >Backflow<.

Order No. Description

8 980 701 Set of solenoid valves

for loop circuit up to +100 °C

8 970 456 Shut-off valve (suitable up to +90 °C) 8 970 457 Shut-off valve (suitable up to +200 °C)

#### 5.3.1. **Tubing**

#### Recommended tubing:

Order No.	Length			Temperature range
8930008	1 m	CR® tubing	8 mm inner dia.	-20 °C to 120 °C
8930012	1 m	CR® tubing	12 mm inner dia.	-20 °C to 120 °C
8930108	1 m	Viton tubing	8 mm inner dia.	-50 °C to 200 °C
8930112	1 m	Viton tubing	12 mm inner dia.	-50 °C to 200 °C
8930410	1 m	Insulation for tubing	8 mm inner dia.	-50 °C to 100 °C
8930412	1 m	Insulation for tubing	12 mm inner dia.	-50 °C to 100 °C
8 930 209	0.5 m	Metal tubing, triple in	nsulated, M16x1	-100 °C to +350 °C
8 930 210	1.0 m			
8 930 211	1.5 m			
8 930 214	3.0 m			
8 930 220	0.5 m	Metal tubing, insulate	ed, M16x1	-50 °C to +200 °C
8 930 221	1.0 m		•	
8 930 222	1.5 m			
8 930 223	3.0 m			



#### Warning: Tubing:

At high working temperatures the tubing used for temperature application and cooling water supply represents a danger source.

A damaged tubing line may cause hot bath fluid to be pumped out within a short time. This may result in:

- Burning of skin
- Difficulties in breathing due to hot atmosphere

#### Safety recommendations


- Employ suitable connecting tubing.
- Make sure that the tubing is securely attached.
- Avoid sharp bends in the tubing, and maintain a sufficient distance from surrounding walls.
- Regularly check the tubing for material defects (e.g. for cracks).
- Preventive maintenance: Replace the tubing from time to time.


# 5.4. Filling / draining



#### Notice:

- Pay attention to the thermal expansion of bath oil during heating to avoid overflowing of the liquid.
- Do not drain the bath fluid while it is hot!
  Recommendation: Temperature range 5 °C to 40 °C
  Check the temperature of the bath fluid prior to draining (by switching the unit on for a short moment, for example).
- Store and dispose the used bath fluid according to the laws for environmental protection.

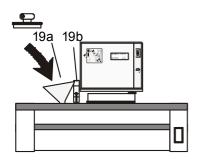

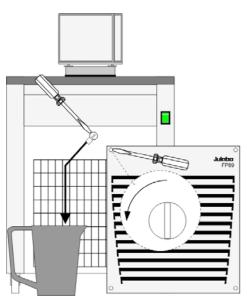

#### **Filling**

Take care that no liquid enters the interior of the circulator. Remove the bath cover and, if necessary, use a funnel. Observe the level indication (19b).

- Recommended maximum filling level with water as bath fluid: 30 mm below the tank rim
- Recommended maximum filling level with bath oils:
   40 mm below the tank rim

#### Special unit with bath opening 28x23 cm

- (i) After filling, immerse the samples in the bath, or place the lid on the bath if the opening is not to be used.
- (i) The circulator has an early warning system for low level (see description on page 38) that may be triggered when changing samples in the bath.

#### Draining:

- The drain tap is located behind the venting grid on the front of the unit, which can be taken off by removing 4 screws.
- Slide a short piece of tube onto the drain connection and hold it into a container.
- Loosen the drain screw a few turns and drain the unit completely.
- Tighten the drain tap.

#### Penetration of ice flakes/water into bath oil

At working temperatures below 0 °C humidity from the atmosphere may penetrate into bath oil through condensation. This may subsequently result in the formation of ice flakes, preventing the pump from proper operation.

For this reason, the bottom of the bath tank provides a groove sloping down towards the drain tap. Thus water having penetrated into the bath oil gathers in the groove immediately behind the drain tap.

Shut down the refrigerated circulator at a bath temperature of +10 °C. Wait for a while and open the drain tap to drain the water.


# 6. Operating procedures

#### 6.1. Power connection



#### Caution:

- Only connect the unit to a power socket with earthing contact (PE protective earth)!
- The power supply plug serves as safe disconnecting device from the line and must be always easily accessible.
- Never operate equipment with damaged mains power cables.
- Regularly check the mains power cables for material defects (e.g. for cracks).
- We disclaim all liability for damage caused by incorrect line voltages!

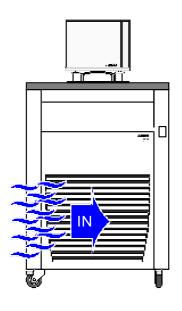


Check to make sure that the line voltage matches the supply voltage specified on the identification plate.

- Connect the circulator with mains power cable (17a) to the mains outlet (18).
- Connect the control cable (9) to the \$\pi\$ sockets (9a, 9b).
- Connect the refrigerated circulator with mains power cable (17b) to the mains socket.
- (i) Note the direction of rotation of the built in fan! For proper cooling of the compressor, the air must be drawn in from the front of the unit.
- For the supplementary heater see chapter 16. HST booster heater
- For the supplementary pump see chapter
   17. HSP booster pump

# 6.2. Switching on / Selecting the language

#### 6.2.1. Switching on the circulator



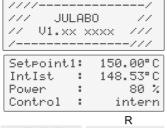



#### Switching on:

- Turn on the mains power switch (1a).
- (i) The unit performs a self-test.

  Then the software version (example: V 1.xx-xxxx) appears. The display "OFF" or "R OFF" indicates the unit is ready to operate.
- The circulator enters the operating mode activated before switching the circulator off:

**keypad control mode** (manual operation) or **remote control mode** (operation via personal computer).


#### 2 KONFIGURATION ≜Zeit/Datum →Sprache deutsch ▼Einheit °C

#### 2 CONFIGURATION ▲Time/Date →Language english ▼Unit °C

#### Selecting the language:

There are two options for the language of the LCD DIALOG-DISPLAY: German or English. Select the desired language in the menu >Configuration< under the submenu >Language/Sprache<.

Press the respective keys in the following order:

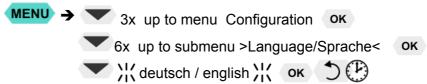

#### 6.2.2. Switching on the cooling machine



Switching on:

• Switch on the cooling machine using the switch (1b).



## (i) Control of the cooling machine

With the mains switch (1b) turned on, the circulator automatically switches the cooling machine off and on.

- It is switched off if:
  - the current working temperature is increased by > 30 °C (cooling is not required).
  - the heater operates at full power (> 800 W) for longer than 5 minutes.
- It is switched on if:
  - cooling is necessary to maintain the bath temperature.

    After switch-off, the cooling machine automatically switches on only after a delay of 5 minutes to protect the cooling compressor.
- (1b) To save energy, turn off the cooling machine with the mains switch (1b) whenever cooling is not required.

# 6.3. Start - Stop



#### Start:

Press OK key.
 The actual bath temperature is displayed on the VFD COMFORT-DISPLAY. The circulating pump starts with a slight delay.

#### Stop:

Press OK key.
 or
 Keep Description

The VFD COMFORT-DISPLAY indicates the message "OFF".

#### **Setting of temperatures** 7.

key to call up the menu for temperature selection.

3 different working temperatures can be adjusted. Their values are freely selectable within the operating temperature range.


(i) The temperatures can be set in start or stop mode.

#### Setting of working temperature in the





Select SETPOINT 1 or 2 or 3 using the key or \_\_\_\_\_\_.

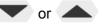

- 3. Confirm by pressing the **OK** key.
- The circulator uses the new working temperature value for temperature control.

#### **Example:**

#### Adjustment/modification of the pre-setting of "SETPOINT 3"

(i) If the active Setpoint is changed, the new value is immediately used for the control of the working temperature. The heater control indicator flashes.

Setpoint3: Change the value from 70.00°C to 80.00°C. See standard display line 1

Use the numeric keypad to enter 8 and 5. Then press enter **OK** to store the value.



Notice: Refer to SETPOINT MAX / MIN in chapter "LIMITS"

# 8.

# Safety installations, warning functions



Check the safety installations at least twice a year! Refer to page 21.


Settings for the excess temperature protection > SafeTemp< and for the warning functions > OverTemp< and > SubTemp< are made in a menu which is called up by pressing the key

Menu item > Type< allows choosing between a warning and an alarm cut-off for the menu items > OverTemp< and > SubTemp<.

# ▼Setpoint2 = 37.00°C Setpoint3 = 70.00°C

Factory settings:

T SETPOINTS

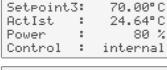

⇒Setpoint1 = 25.00°C

85.**∭**0°C Setpoint3: ActIst 24.64°C 80 % Power . Control internal




#### Warning:



The excess temperature protection **must** be set at least 25 °C below the fire point of the bath fluid used!

In case of wrong setting there is a fire hazard!

We disclaim all liability for damage caused by wrong settings!

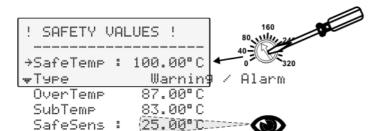

#### SafeTemp:

Setting range: 20 °C ... 320 °C

(i) Rough setting can be effected by using the temperature scale.

#### Settings:

• Press the key and by pressing the key or select the menu item : and confirm the by pressing the key or .

#### For setting proceed as follows.

SafeTemp:

Set the new cut-out value within 30 seconds with a screw driver.
 Exact display and setting on LCD display
 Recommendation:
 Set over temperature-protective setting >SafeTemp< 5 °C to 10.</li>

Set over temperature-protective setting >SafeTemp< 5 °C to 10 °C above the set point of the working temperature.

Type:

Switch-over from warning to shutdown function

Change the parameter by pressing the key pressing the key ok or
 or pressing the key if the parameter is to be retained.

Possible parameters: warning or alarm

OverTemp:

Set value on numeric keypad and confirm by pressing

SubTemp:

Example: 83.00 °C

Press the keys 8 and 3 and confirm by pressing OK

SafeSens:

Indicated is the temperature value of the safety sensor.

#### SafeTemperature:

Setting range: 20 °C ... 320 °C

Indicated is the cut-out value set with a screwdriver on the excess temperature protection device.

This safety installation is independent of the control circuit. When the temperature of the bath fluid has reached the safety temperature, a complete shutdown of the heater and pump is effected.

The alarm is indicated by optical and audible signals (continuous tone).

The following error message appears on the VFD COMFORT-DISPLAY:







Press **OK** for help on the LCD **↓** 

Temperature exceeds the adjusted safety-temperature. Check adjustment!

① Depending on the setting of >ControlType - internal or external< the actual temperature values for both sensors are shown/indicated on both displays.

Examples: I 94.06/1\(\text{17}\) and Ext. : xx.xx°C


## **Type:** >Warning< or >Alarm<

For the two menu items >OverTemp< and >SubTemp< choose between a warning message being signaled or a complete shutdown of the main functional elements such as heater and circulating pump being effected.

#### **OverTemp**erature: **SubTemp**erature:

If for a sensitive temperature application task adherence to a working temperature value >Setpoint< is to be supervised, then set high and low temperature warning values.

In the example below, the >Setpoint< of 85 °C is surrounded by the values >OverTemp< 87 °C and >SubTemp< 83 °C. The electronics immediately registers when the actual temperature attains a temperature out of the limits and it follows a reaction according to what is set in the menu item >-Type<.

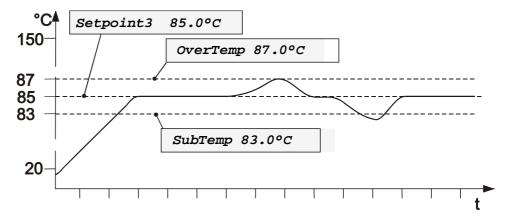

(i) The warning functions are only triggered when the actual bath temperature, after start from the "OFF" or "rOFF" mode, lies within the set limits for 3 seconds.

#### Setting >WARNING


A mere warning function with optical and audible warning signal (interval tone) A message appears on the VFD COMFORT-DISPLAY:





#### The LCD –DISPLAY shows the message:

Setpoint1: 150.00°C IntAct : 148.53°C excess temperature! -check limits-

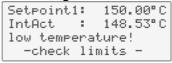

#### Setting >ALARM

Temperature limit with shutdown of heater and circulating pump. An audible alarm sounds (continuous tone) and a message appears on the VFD COMFORT-DISPLAY:



**- [] F F -**RL RRM CO JE O Y

OverTemp

SubTemp

#### The LCD -DISPLAY shows the message



A L A R M !
Ext. : xx.xx°C
Low Temperature!

Press **OK** for help on the LCD

Actual temperature obove high temperature value.
Check limits!

Actual temperature below low temperature value. Check limits!

#### 8.0.1. Early warning system, low level protection



DBGM: 20306059.8

This low level protection is independent of the control circuit and is divided into two sections:

1. Switch in stage 1 recognizes a critical fluid level . An audible warning (interval tone) sounds and a message appears on the VFD COMFORT-DISPLAY.




Setpoint1: 150.00°C IntAct : 148.53°C low level! -refill medium-

Refill bath fluid!

2. Switch in stage 2 recognizes a low fluid level 🙃.

If stage 2 of the low level protection device (according to IEC 61010-2-010) is triggered, a complete shutdown of the heater and circulating pump is effected.

A continuous alarm tone sounds and a message >ALARM< >CODE 01< appears on the VFD COMFORT-DISPLAY.

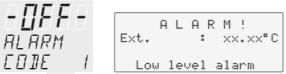

Press OK for help on the LCD ♥

Low liquid level or float brocken. Check float and fill medium!

Turn off the unit with the mains switch, refill bath fluid and turn the unit on again!

3. Float

5

- 4. Circulating pump
- 5. Heater



#### Warning:

For refill always use the same bath fluid type that is already in the bath. Bath oils must not contain any water contaminants and should be pre-heated to the actual bath temperature! Explosion hazard at higher temperatures!

# 9. Menu functions

MENU

The term "menu functions" refers to adjustments such as



1 MAIN MENU

→Start Program ▼Manage Programs

Pump
Configuration
Control
Interface serial
ATCalibration
Limits
Inputs/Outputs

#### Menu level 1

 Start program

Page 41

This menu will start a previously set program.



1 MAIN MENU ▲Start Program →Manage Programs ▼Pump Administration and creation of programs

Page 46



1 MAIN MENU Manage Programs →Pump ▼Configuration Electronically adjustable pump capacity

Page 50

The capacity of the circulating pump is set by adjusting the motor speed.



1 MAIN MENU ◆Pump →Confi9uration ▼Control Configuration of the unit

Page 51

Page 58

Remote – on / off (remote control via RS232) Setpoint ext - Switch over external setpoint setting

Autostart - AUTOSTART on / off

Off-Mode – Pump on / off

ACTVAR - Switch over input variable


Time / Date – setting time and date

Language – english / deutsch

Unit – Temperature values can be displayed in °C or °F.

Program Type – Time or Gradient (grad)

Reset – factory settings

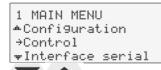

Control characteristics and parameters

ControlType - Control internal or external

Selftuning

Dynamic - internal

CoSpeed – external

Control parameters - XP-, TN-, TV- internal

Control parameters - XP-, TN-, TV- XPU-, external

#### Menu level 1

# Adjustable interface parameters BAUDRATE, PARITY, HANDSHAKE

Page 64



→ATCalibration •Limits ATC - Absolute Temperature Calibration, Sensor calibration internal sensor, Sensor calibration external sensor

3-point calibration

Page 65

Page 72

1 MAIN MENU ♠ATCalibration →Limits ♥Inputs/Outputs Limitations of temperature and capacity

Setpoint Max / Min - Maximum and minimum setpoint

Heating Max – Set maximum heating

Cooling Max - Set maximum cooling

Internal Max / Min – Limitation of the temperature range

Upper band / Lower band – Band limit

1 MAIN MENU ▲Limits →Inputs/Outputs Analog inputs/outputs

Recorder output - CHANNEL 1, 2, 3

EPROG – External programmer input

Ext-StBy - Stand-By input

ALARM - output

Page 74

## 9.1. Start of a program



The start menu of the integrated programmer allows one of six previously stored temperature programs to be called up and started. The profiles are started manually or via the integrated timer.

### (i) Requirements:

- 1. Create a program. (refer to next chapter)
- 2. Return to the Start-MENU and confirm the desired setting of each MENU item with the key OK
- 3. Set a start time (>TIME< >DATE< >YEAR<) if the program is to be started by the internal timer.

>Program Number< Select the program that should be started.

0 ... 5

> at Step< Select the Step at which the program should be

started. 0 ... 59

> Runs < Select the number of program repetitions. 1 ... 99


> Program End< Status at program end .


StdBy, PSetp, Setp 1, Setp 2, Setp 3

> RUN < Start time (now / time)

① Press the 3 key if a parameter is to be retained. (prior to OK)



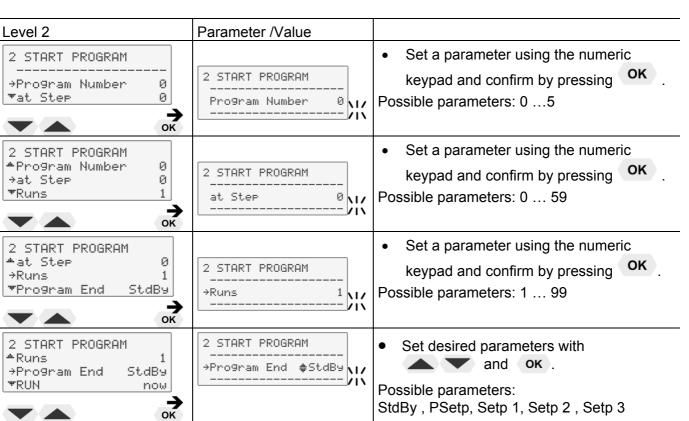

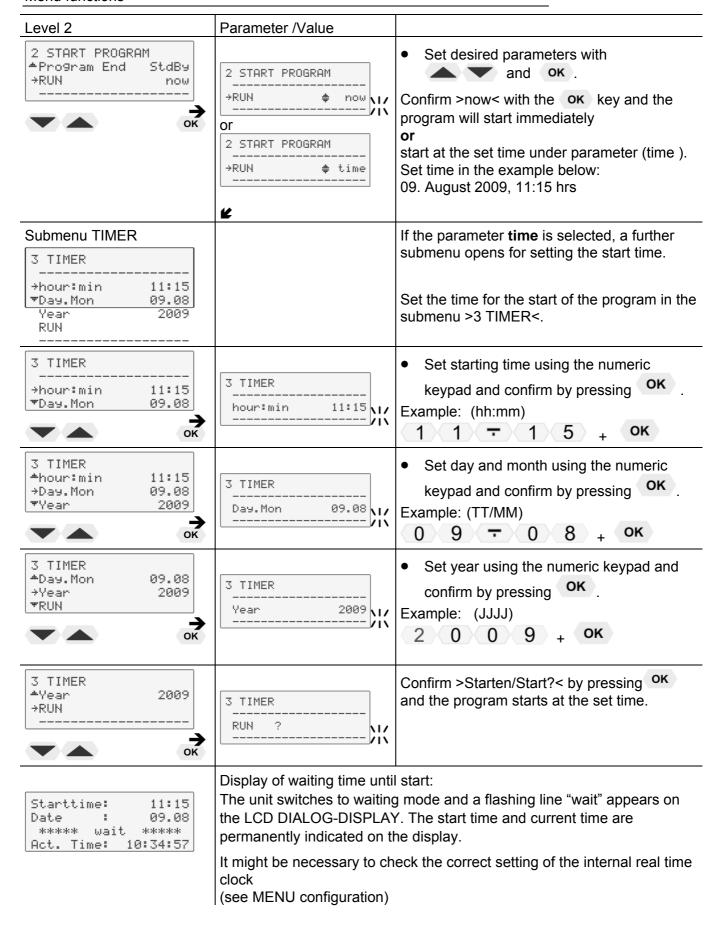

#### 9.1.1. Status at the end of a program

Setpoint	.3:	80.00°C
IntAct	:	80.00°C
Power	:	3 %
Control	:	internal

T SETPOINT	S	
<b>-</b> Setpoint1	=	25.00°C
⇒Setpoint2	=	58.00°C
₩Setpoint3	=	85.00°C
Setpoint2:		58.00°C
IntAct :		60.00°C
Power :		0 %
Control :		intern



#### StdBy, PSetp, Setp 1, Setp 2, Setp 3 Program end

This parameter is used to determine if the unit switches to the -OFFcondition at the end of a profile or if temperature control is continued. and the working temperature setpoint that is to be used if temperature control is continued.

**StdBy** – If the end of the profile is reached, the circulator switches to the -OFF- condition. The VFD display indicates the working temperature setpoint that was displayed in line 2 before the program was started (example: S 25.00 °C).

**PSetp** –(Setpoint of the **ProG**rammer) At the end of the program, the setpoint of the last program step is written to the position of setpoint 3 (example: 80.00 °C) and indicated in line 1 of the LCD display. The circulator operates and controls to this temperature until the key **OK** is pressed or a new profile is started.

#### Setp 1, Setp 2, Setp 3 – ("Setp" short for Setpoint)

Before starting the programmer, a temperature value is set for one of the three setpoints in the **T** menu.

(example: setpoint 2 58.00 °C)

The circulator operates and controls to this temperature at the end of the profile until the key **OK** is pressed or a new program is started.

#### 9.1.2. Indication after successful start

The started programmer displays the currently calculated setpoint in line 1. The value changes until the final temperature of the section is reached (temperature ramp).

1<sup>st</sup> line: Setpoint of the programmer

2<sup>nd</sup> line: Actual temperature value

> for **internal control** = IntAct: XXX.XX for external control = ExtAct: xxx.xx

3<sup>rd</sup> line: Selected profile and the current section (step).

4<sup>th</sup> line: With the keys, the following information can be

displayed in line 4.

- Remaining time of the current section (standard display)
- Total remaining time: profile x number of repetitions (All: Days/hours:minutes:seconds)

#### Particularities:

1. If a total time of more than 999 days is calculated, only the hours are displayed by the unit.

(All: hours:minutes:seconds)

2. If the time period of a section is edited after the start, then the remaining total time can no longer be displayed.

### Examples:

Set	:	XX.XX°C
IntAct	:	XX.XX°C
ProgNr.	:2	Step: 2
Remain	:	03:32:25

Set	#		XX.X	ΚΧ¤	C
IntAct	:		XX.X	XX"	С
ProgNr.	i	2	Step:		2
911	=	t.t.t./	hh:mr	n: <	<

#### 9.1.3. Editing after Start



The programmer allows a running program to be edited.


#### Initiation and exit of the editing function:

- After the start enter the submenu "3 EDITOR".by pressing the key
  - See chapter "program administration, creation" page 46
- Exit the editing modus anytime by pressing the key .


#### Particularity for changes in a section that is currently in process.

- (i) If the target temperature is changed, the slope of the ramp is automatically calculated based on the remaining time of the section.
- (i) If the time period is changed, the slope of the ramp is automatically calculated based on the new total time of the section.
- (i) The "remaining total time" (profile x number of repetitions) as described on page 43 can no longer be displayed.

#### 9.1.4. Interrupting a program

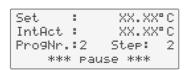

- Press the hand time period set for the corresponding section are
  - thus stopped at the values presently achieved. The instrument is put on hold and the message "pause" flashes on the LCD DIALOG DISPLAY.
- (i) A program can be interrupted or restarted by an external shut-off.

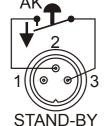

## Important:

For this, in the >Inputs/Outputs< menu, the menu item >Ext. StBy< must be set to >active< and in the >Configuration< menu, >Autostart< needs to be set

(i) Setpoint control and timer are interrupted by breaking the contact "AK".

The instrument is put on hold and the message "Stand-By" flashes on the LCD DIALOG-DISPLAY.



Notice: This is not an actual emergency switch-off.



H

\*\*\*\*\* StandBy \*\*\*\*\*

Set

IntAct :

ProgNr.:2

#### Warning:

XX.XX°C

XX.XX°C

Step:

Following a power interruption, it would be possible in this condition for the instrument to restart automatically. The safety and warning functions of the instrument should always be used to their fullest capacity.

See **Warning** page 55.

#### 9.1.5. Interruption after a power failure

It is possible to operate the programmer safely in case of a power failure.


#### Important:

For this, >Autostart< needs to be set in the >Configuration< menu.

- i If the AUTOSTART function is activated, the programmer starts again at a point approx. 20 seconds before the interruption took place. However, an uncontrolled change of the bath temperature has occurred.
- (i) Power failure when the instrument is on hold if start is effected via the built-in timer:


If the starting time is not yet reached, the programmer is put on hold again.

If the starting time is exceeded, the programmer starts immediately.

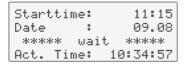

#### 9.1.6. Termination of a program

Set : XX.XX°C IntAct: XX.XX°C Pro9Nr.:2 Step: 2 Remain: 03:32:25

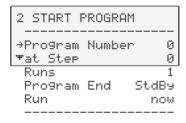

(1) A program can be terminated by pressing (2). The programmer switches back to the Start menu.

Further procedures are defined in the MENU >START PROGRAM<, menu item >Program End <.

Example: >Program End - StdBy<



The VFD COMFORT-DISPLAY displays "OFF".

#### Interruption after a power failure

In case of a power failure, the reaction of the circulator is the same as when switched off and on again with the mains switch.

The VFD COMFORT-DISPLAY displays "OFF".

## 9.2. Program administration, creation

2 MANAGE	PROGRAMS
 >Edit	PrgNr.
<b>™</b> Delete Print	Pr9Nr. Pr9Nr.

The integrated programmer permits fast and easy programming of setpoint temperature sequences. This temperature sequence is called program. A program is composed of individual sections (Step). The sections are defined by duration (t) or Gradient (°/t) and target temperature. The target temperature is the setpoint (SP), which is achieved at the end of a section. The programmer calculates the temperature ramp from the difference in time and temperature.

#### **Create programs**


A flashing segment indicates that a number needs to be entered. Under the submenu "Edit", enter a program number. Six programs may be stored (numbers 0 to 5).

When the program is running, only sections with complete information for the target temperature and time period are considered.

It makes sense to skip section numbers in the profile in order to use them later for corrections to the profile.

#### Important:

If a time of 00:00 is set for a profile, the profile is continued with the next section only after the setpoint temperature (±0.2 °C) is reached.

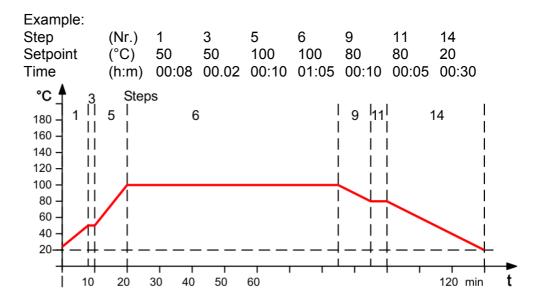

### **Editing**





1 MAIN MENU ▲Start Program →Manage Programs ▼Pump

ок 🖊

PrgNr: Set the program you wish to create.  $0 \dots 5$ 

SP: Setpoint

TI: Time [hh:mm] or GRD: Gradient  $[^{\circ}/t] = [^{\circ}/min]$ 

Change type of program in MENU >configuration<,

menu item >Programmtyp/program type< .

ST: Step 0 ... 59

	(i) Pr	ess the 5 key if a parameter is to be retained. (prior to OK)
Level 2		
2 MANAGE PROGRAMS		Example: Program (PrgNr.) 2 is to have a target temperature with a set point (SW) of 100.00 °C in section t (ABS) 3. Duration (ZI) 1 hour.
2 MANAGE PROGRAMS		Select the program number.
Edit # Pr9Nr.2	Ж	Set the desired >PrgNr.< by pressing     and ok .
Level 3		For the selected section >ST< a setpoint >SP< and the duration >TI< or the gradient >GRD< are entered in the right half of the input window.
3 EDITOR >PrgNr.2	\ \ \ \	Input window for >ST<, >SP< and >TI< / >GRD<
SP:C 100.00C TI:T 1.00T ST: 2 # 3		1. Set the desired Step >ST < by pressing and ok .
or 3 EDITOR >PrgNr.2		Set set point >SP< by using the numeric keypad and confirm by pressing OK .
SP:CC GRD:%% ST: # 0		3. Set the time >TI< or a gradient >GRD< using the numeric keypad and confirm by pressing OK .
(m)		By confirming time or GRD the values on the left side of the input window are moved.
3 EDITOR >Pr9Nr.2 SP: 100.00CC TI: 1.00TT ST: 3 # 4	\\\ \\\	<ul> <li>The values for the next section are entered in the right half of the input window.</li> <li>See above 1. / 2. / 3.</li> </ul>

#### **Delete**

It is possible to delete an entire program or consecutive sections

Set the program you wish to delete Delete PrgNr:

0 ... 5

Delete from Step X to Step X - (0 ... 59)





1 MAIN MENU ∸Start Program ⇒Mana9e Pro9rams ₩Pump

Press the be retained. (prior to OK)





Level 2	Submenu >Delete<
2 MANAGE PROGRAMS  Edit PrgNr.  Delete PrgNr.  Print PrgNr.	Set menu item >Delete PrgNr.< by pressing and ok .
2 MANAGE PROGRAMS	• Set desired PrgNr with  and  ok .
Delete # PrgNr.0	
Level 3	Input window: delete sections.
3 DELETE >Pr9Nr. 0	
→from Step 0 ▼to Step 59 delete	• Select menu items > → from Step< and >→ to Step< one after the other by pressing and oκ.
	Select the desired section for every menu item by using the
3 DELETE >Pr9Nr. 0	numeric keypad and confirm by pressing <b>ok</b> .  Example: >from 0< >to 59<
from Step 0	,
3 DELETE >Pr9Nr. 0	
to Step 59	
	• Select menu item >delete< by pressing ▼ and OK
3 DELETE >Pr9Nr. 0	• Confirm >Delete?< by pressing OK
delete?	
3 DELETE >Pr9Nr. 0	① The delete procedure is confirm by indication of the
! delete !	message >delete< on the LCD display.

### **Print**





1 MAIN MENU ∸Start Pro9ram →Mana9e Pro9rams ₩Pump

Every program can be printed via the serial interface for control and documentation purposes.



Level 2	Submenu >Print<
2 MANAGE PROGRAMS ◆Delete Pr9Nr. →Print Pr9Nr.	Set menu item >Print PrgNr.< by pressing and ok.
2 MANAGE PROGRAMS Print # Pr9Nr.4	Set desired >PrgNr.< by pressing     and ok     Example: 4
Level 3	
3 PRINT Pr9Nr.4  →Print	• Confirm → > Print< by pressing OK .
3 PRINT Pr9Nr.4 Print?	Confirm >Print?< by pressing οκ .
3 PRINT Pr9Nr.4 ! print!  3 PRINT Pr9Nr.4 done!	Printing and end of printing are confirmed by successive messages on the LCD display as shown on the left.


#### Setting the pump pressure 9.3.





1 MAIN MENU ▲Manage Programs ⇒Pump ⇔Confi9uration OK T

The capacity of the circulating pump is set by adjusting the motor speed

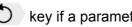

① Press the bkey if a parameter is to be retained. (prior to OK)

Level 2		Submenu >Pump<
2 PUMP 		The set pump stage is displayed. Example: 1  • Confirm by pressing OK .
2 PUMP 	\ / / \	Set required pump stage by pressing     and ok     Example: 3

### Example:

Adjusted

Effective



Adjustable pump capacity stage 1 ... 4

Illuminated display: For pump pressure


- i Indication of the adjusted pump pressure stage in the –OFF- mode.
- i Indication of the effective pump pressure stage (rotational speed) after start.


To protect the pump motor, the rotational speed changes with the load applied, for example, depending on the viscosity of the bath fluid at different working temperatures.

22 ... 26 I/min Flow rate:


Pump capacity stage		1	2	3	4
Pump pressure	[bar]	0.4	0.5	0.6	0.7
Suction pump	[bar]	0.2	0.26	0.33	0.4
Total capacity in a loop circuit	[bar]	0.6	0.76	0.93	1.1

#### 9.4. Configuration

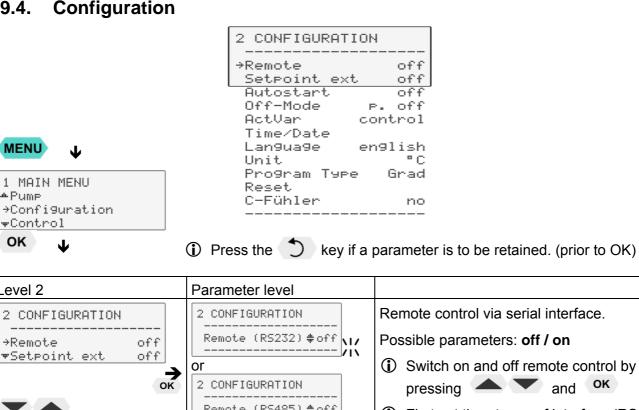

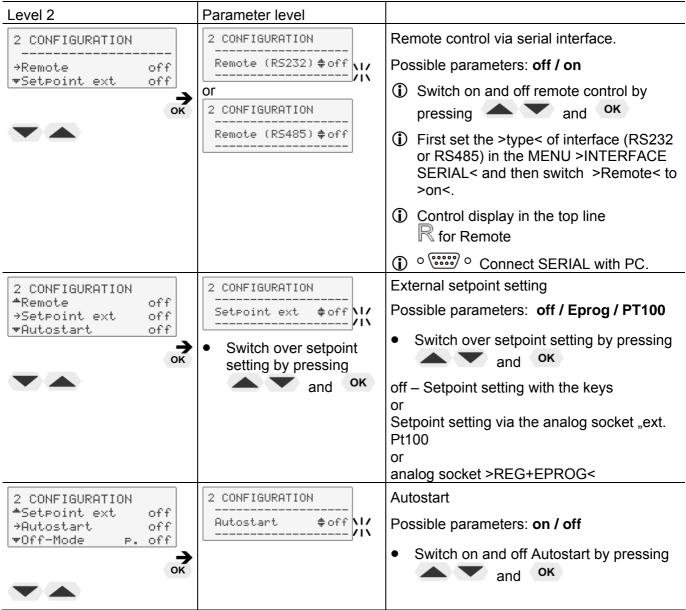

Level 2	Parameter level	
2 CONFIGURATION #Autostart off +Off-Mode p. off #ActVar control OK	2 CONFIGURATION	Pump-Mode Possible parameters: P. on / P. off  • Switch on and off >Off-Mode by pressing and ok  P.on continuous operation of circulating pump  P.off circulating pump is linked to Start/Stop
2 CONFIGURATION  *Off-Mode p. off  →ActVar control  *Time/Date	2 CONFIGURATION	ActVar - actuating variable  Possible parameters: control, Eprog, serial  i Set desired parameters with and ok.  i Programming of variables for the parameters > serial < or > EPROG < is only accepted, if the unit is in Start mode
2 CONFIGURATION  ActVar control  Time/Date Language english		Time, Date, Year see next page
2 CONFIGURATION #Time/Date +Language english #Unit °C	2 KONFIGURATION  Language \$english	Language Possible parameters: deutsch / english  • Set desired parameters with  and ok .
2 CONFIGURATION *Language english *Unit °C *Program type time  OK	2 KONFIGURATIONUnit	Unit  Possible parameters: °C / °F  • Set desired parameters with  and ok .
2 CONFIGURATION  *Unit °C  *Program type time  *Reset  OK	2 KONFIGURATION	Program type  Possible parameters: time / grad  • Set desired parameters with and ok .

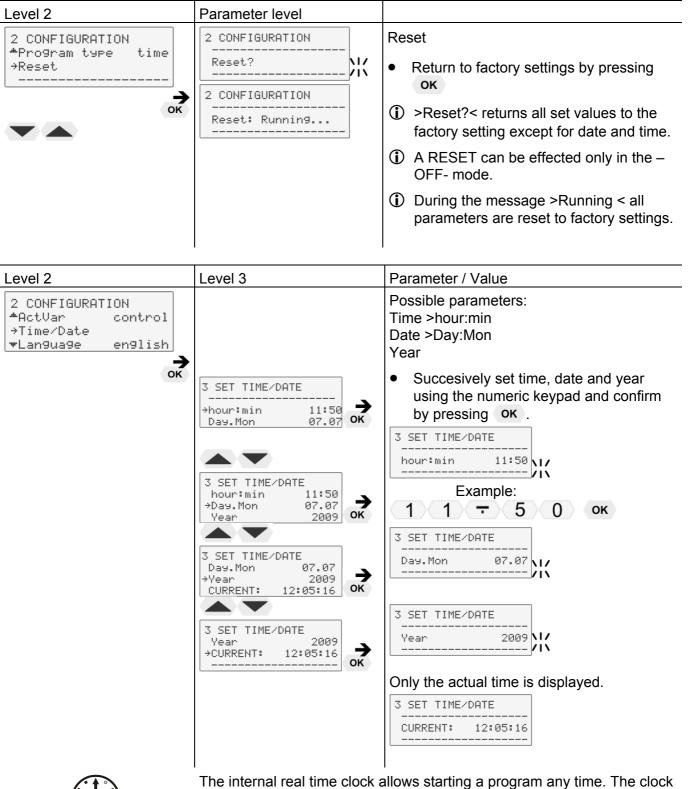



is set to the local mean time (MEZ) at the factory.

- (i) If the unit is operated in a different time zone, the clock can be adjusted in this menu.
- (i) Change summer/winter time in this menu

#### 9.4.1. Remote control via the serial interface

#### Factory setting: off

,	0	
RS232	:	25.00°C
IntAct	:	27.00°C
Power	:	0 %
Control	:	intern
RS485	:	25.00°C
IntAct		27.00°C
Power	:	0 %
Control	:	intern

Setpoint is set via the serial RS232/RS485 interface through a PC or superordinated data system.

In the header of the VFD COMFORT-DISPLAY, an "**R**" illuminates. It indicates that remote control mode is active.

← The selected setting is shown on the LCD DIALOG-DISPLAY.

### o o SERIAL


**IMPORTANT:** additional measures for remote control

- (i) RS232 or RS485:
- First set >type< of interface in the MAIN MENU > Interface serial < the switch >Remote< to >on<.
- (i) Connect the circulator to the PC using an interface cable.
- (i) Check the interface parameters of both interfaces (circulator and PC) and make sure they match.

(See 12.1. Setup for remote control page 90)

#### 9.4.2. Keypad control or setpoint setting via the analog input

#### Factory setting: off

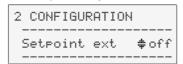

152.00°C 50.64°C

100 %

The selected mode is indicated on standard display.

Setpoint1:

IntAct :

Power :

Control	:	intern
Pt100 IntAct Power	:	85.00°C 83.00°C 60 %
Control	:	external
Epro9 IntAct	:	50.00°C
Power	:	80 %

Control : internal

2 CONFIGURATION

→Remote off

Setpoint ext Epro9

Autostart off
Off-Mode p. off

ActVar Epro9

Time/Date

In addition to the serial interface via remote control the circulator offers the possibility to adjust the setpoint via analog interface >ext. Pt100< or >REG+E-PROG<.

#### Possible parameters:

**off** - Setpoint is set via the keypad or via the integrated programmer. (factory setting

**Pt100 -** Setpoint setting via the analog socket "ext. Pt100" using an external temperature sensor or an appropriate voltage/current source.

**Eprog** - Can only be adjusted when an electronic module with analog connections is used (option).

Setpoint setting via the analog interface REG+E-PROG connection with an external voltage or current source or a programmer.

#### Important:

- (i) Connect the external voltage or current source or a programmer to the circulator via the socket REG+E-PROG (see page 80).

  In the menu >MENU Inputs/Outputs< set the parameter >Erog-INPUT< and the input variables >Eprog-signal< (see page 80).
- The E-Prog input can only be used either under menu item >Setpoint ext < or under menu item >ActVar<</p>

#### 9.4.3. Autostart



#### Possible parameters:

on - Autostart on

off - Autostart off (Factory setting)

#### Note:

The circulator has been configured and delivered by JULABO according to N.A.M.U.R. recommendations. This means for the start mode that the unit must enter a safe operating state after a power failure (non-automatic start mode). This safe operating state is indicated by "**OFF**" or "**R OFF**" on the VFD-COMFORT-DISPLAY. A complete shutdown of the main functional elements such as the heater and circulating pump is effected simultaneously.

The values set on the circulator remain stored, and the unit is returned to operation by pressing the start/stop key (in manual control mode). In remote control mode, the values need to be resent by the PC via the interface.

Should such a safety standard not be required, the AUTOSTART function (automatic start mode) may be activated, thus allowing the unit to be started directly by pressing the mains power switch or using a timer.

The AUTOSTART function can be used only if the setpoint is set via >keyboard<, >eprog<, or >ext. Pt100<.



#### Warning

For supervised or unsupervised operation with the "AUTOSTART" function avoid any hazardous situation to persons or property

Take care to fully observe the safety and warning functions of the circulator.

#### 9.4.4. Off-Mode



#### **OFF-Mode**


Normally the circulating pump is switched via the start/stop signal. However, if circulation should be maintained even in the -off- condition, the parameter >pump on< needs to be set.

#### Possible parameters:

P.on Circulating pump in continuous operation

P.off Circulating pump linked to Start/Stop (factory setting)


#### 9.4.5. ActVar - actuating variable

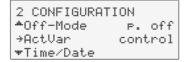

The variable (ACTuating VARiable) corresponds to the extent to which the heater or cooling unit of the circulator is controlled. Heat or cold is applied to the bath according to this variable. If this happens with the control electronics of the circulator, called > control < in this particular case, the bath temperature is exactly heated and maintained constant at the adjusted setpoint.

(i) Programming of variables for the parameters > SERIAL < or > EPROG < is only accepted, if the unit is in Start mode.



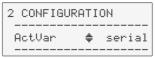



#### Possible parameters:

control – The internal control electronics of the circulator controls the heater and the connected cooling unit. Self-tuning is possible.. (factory setting)

**serial** — The heater or the connected cooling unit receives the control signal via the serial interface. Self-tuning is not possible.

**Eprog -** The heater or the connected cooling unit receives the control signal via the E-Prog input. Self-tuning is not possible.

- Setting requires electronic module.

#### Important:

Set the parameter >Eprog< and the input variables also in the menu >Inputs/Outputs< (see page 80).

#### 

#### Note:

The E-Prog input can only be used either under menu item >Setpoint ext< or under menu item > ActVar <.



Time/Date

#### Warning:

The working temperature range of the circulator is determined during configuration. If set to >Control<, this range cannot be exceeded.

If set to > Serial < and > Eprog <, heat or cold is applied to the bath without control. The permissible maximum temperature can be exceeded. The user has to take adequate precautions for temperature control.

Materials, such as gaskets or insulations for example, may be damaged or destroyed, if the permissible maximum temperature is exceeded.

The safety and warning functions > < of the instrument must always be used to their fullest capacity.

### 9.4.6. Setting of clock and date



2 CONFIGURATION \*ActVar control →Time/Date \*Lan9ua9e en9lish The internal real time clock allows starting a program any time. The clock is set to the local mean time (MEZ) at the factory.

- i If the unit is operated in a different time zone, the clock can be adjusted in this menu.
- (i) Change summer/winter time in this menu


#### 9.4.7. Language

2 CONFIGURATION \*Time/Date →Language english \*Unit °C There are two options for the language of the LCD DIALOG-DISPLAY: German and English.

#### Possible parameters:

#### deutsch / english

#### 9.4.8. Unit

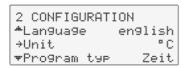

Temperature values can be displayed in °C or °F.


With this change, all adjusted temperature values are converted and displayed in the new unit.

#### Possible parameters:

°C (factory setting)

۰F

#### 9.4.9. Program type

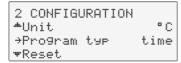


A program can be created with the parameters time or gradient.

#### Possible parameters:

time (factory setting)

**Grd** (Gradient) [°/t] = [°/min]

#### 9.4.10. Reset

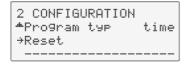


A Reset will return all values to factory setting except for date and time.

(i) A Reset can be effected in the >OFF< mode only.

#### 9.5. Control

The circulator is designed for internal and external temperature control. This can be switched in this submenu. Only the set of parameters corresponding to the particular setting is displayed.

#### Control internal

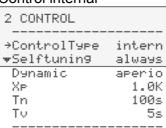

#### Control external


2 CONTROL	
 →ControlType	extern
♥Selftunin9	always
CoSpeed	0.0
XP	0.7K
Tn	720s
Τv	55s
XPu	2.9K







- 1 MAIN MENU ▲Confi9uration →Control ▼Interface serial
- OK J

- (i) For external temperature control and measurement connect a Pt100 external sensor to the socket at the rear of the circulator.
- Press the key if a parameter is to be retained. (prior to OK)

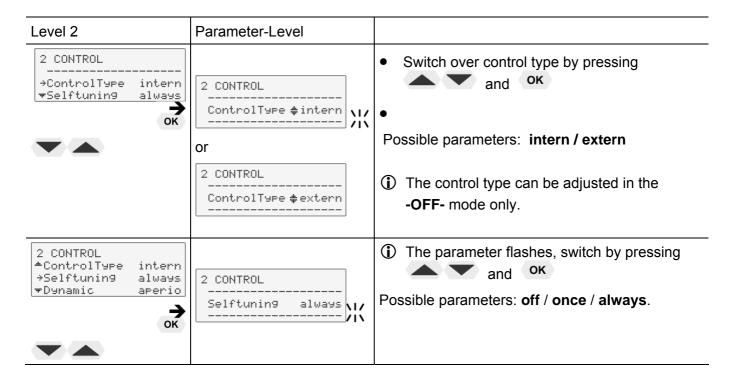

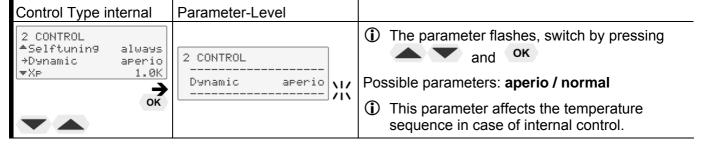

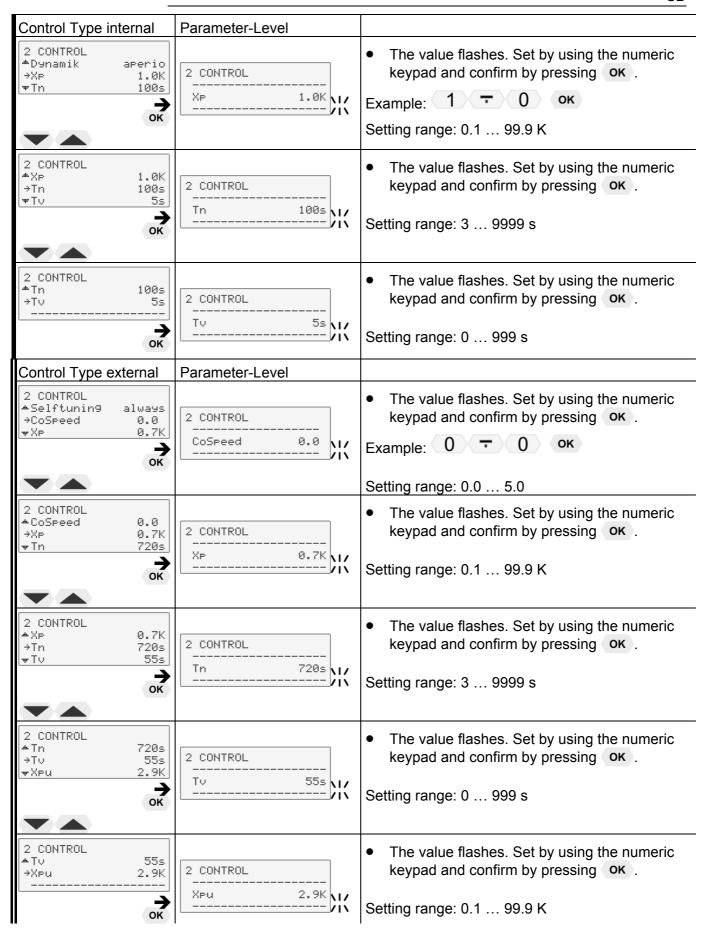

#### 9.5.1. Control internal / external



(i) The control type can be adjusted in the **-OFF-** mode only.

Possible parameters:

**intern** internal temperature control

**extern** external temperature control with external Pt100 sensor



#### **IMPORTANT:**

Additional measures for external temperature control.

- (i) Connect a Pt100 sensor to the socket on the rear of the circulator.
- (i) Sensor calibration of the external Pt100 sensor is performed in the >ATCalibration< menu.
- (i) Suggested adjustments for external temperature control: Upper band / Lower band and IntMax / IntMin see Section >Limits<.
- (i) External temperature control does not allow the setpoint to be set via the "ext. Pt100" socket.



#### Notice:

Place the external sensor into the temperature-controlled medium and securely fix the sensor.



#### Accessory: Pt100 external sensor

Order No.	Description	Material	Cable
8981003	200x6 mm Ø,	stainless steel	1.5 m
8981005	200x6 mm Ø,	glass	1.5 m
8981006	20x2 mm Ø,	stainless steel	1.5 m
8981010	300x6 mm Ø,	stainless steel	1.5 m
8981015	300x6 mm Ø,	stainless steel / PTFE coated	3 m
8981013	600x6 mm Ø,	stainless steel / PTFE coated	3 m
8981016	900x6 mm Ø,	stainless steel / PTFE coated	3 m
8981014	1200x6 mm Ø,	stainless steel / PTFE coated	3 m
8981103	Extension cable fo	r Pt100 sensor	3.5 m
8981020	M+R in-line Pt100	sensor	

The M+R in-line Pt100 sensor is a flow sensor and can be installed loop circuit

#### 9.5.2. Selftuning

When performing a selftuning for the controlled system (temperature application system), the control parameters Xp, Tn and Tv are automatically determined and stored.

#### Possible parameters:

#### off - no selftuning

The control parameters ascertained during the last identification are used for control purposes.

#### once - single selftuning (factory setting)

The instrument performs a single selftuning of the controlled system after each start with the ok key or after receiving a start command via the interface.


### always - continual selftuning

The instrument performs a selftuning of the controlled system whenever a new setpoint is to be reached.

Use this setting only when the temperature application system changes permanently.

#### 9.5.3. Dynamik intern

This parameter affects the temperature sequence only in case of internal control.

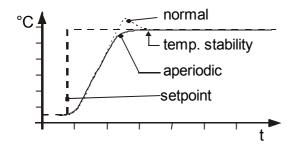

Factory setting: APER (aperiodic)

### Possible parameters:

normal temp. stability aperiodic temperature ramp

**normal** Allows for reaching the setpoint faster – with setpoint change or ramp function – but overshooting of up to 5 % is possible.

aperio Ramp function: the increase of temperature occurs temporally offset and achieves the target temperature without overshooting.

Setpoint change: The temperature increases at the same rate, the target temperature is achieved without overshooting.


(i) With both settings constant temperature is achieved after approximately the same time.

#### 9.5.4. CoSpeed - extern

#### **CoSpeed:**

This parameter affects the temperature pattern only in case of **external** control.

Possible parameters: 0 ... 5

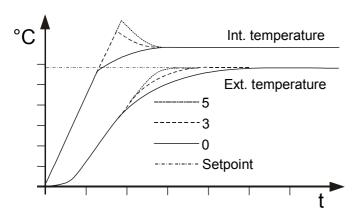

During selftuning, the control parameters Xp, Tn, and Tv of a controlled system are automatically determined and stored. The time required for tuning may vary depending on the controlled system. This controller design allows protection of sensitive objects requiring temperature control. As soon as a co-speed factor is set, it is considered in calculating the control parameters. As shown in the diagram, tuning times are shorter the higher the co-speed factor is, but overshoot can occur in the internal system.

#### 9.5.5. Control parameters – internal/external

#### Control parameters – internal/external

The control parameters preset at the factory are in most cases adequate to achieve an optimal temperature pattern in the samples requiring temperature control.

Each parameter may be manually set via the keypad if necessary to allow optimal control performance.

#### Setting range: internal / external 0.1 ... 99.9 K

#### Proportional range >Xp<

The proportional range is the temperature range below the setpoint temperature value in which the control circuit reduces the heating power from 100% to 0%.

### Setting range: internal / external 3 ...9999 s

#### Resetting time >Tn< (Integral component)

Compensation of the remaining control deviation due to proportional control. An insufficient resetting time may cause instabilities to occur. An excessive resetting time will unnecessarily prolong compensation of the control difference.

#### Setting range: internal / external 0 ... 999 s

#### Lead time >Tv< (Differential component)


The differential component reduces the control settling time. An insufficient lead time will prolong the time required to compensate for disturbance effects and cause high overshoot during run-up. An excessive lead time could cause instabilities (oscillations) to occur.

# Setting range: 0.1 ... 99.9 K

#### Proportional range >Xpu<


The proportional range Xpu of the cascaded controller is only needed for external control.


## **Optimization instructions for the PID control parameters:**

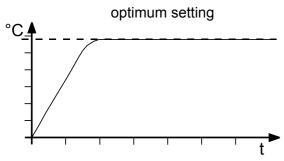



The heat-up curve reveals inappropriate control settings.

## Inappropriate settings may produce the following heat-up curves:

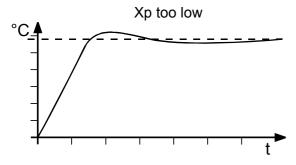

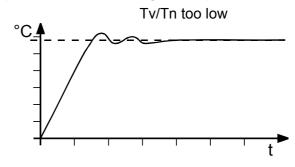

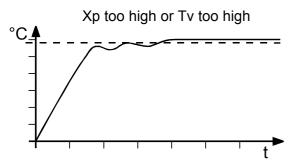


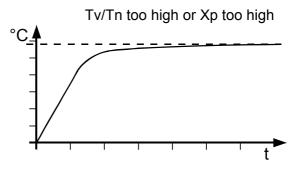

#### 9.6. Serial Interface

Factory settings:
4800 Baud
even
hardware handshake

For communication between the circulator and a PC or a superordinated process system, the interface parameters of both units must be identical. The adjustment is usually carried out one time only in the >Interface serial< menu.

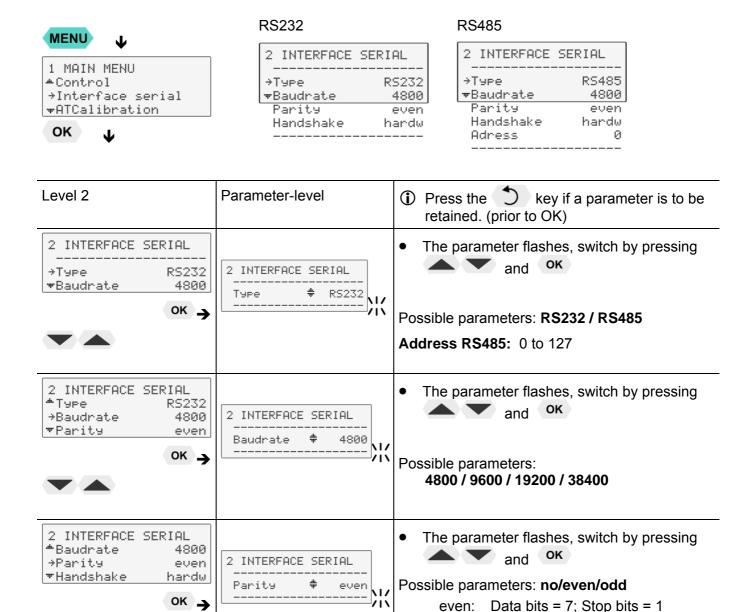

odd:

and οκ

Possible parameters: hardw / softw

Data bits = 7; Stop bits = 1 Data bits = 8; Stop bits = 1

The parameter flashes, switch by pressing

Xon/Xoff-protocol (Software handshake)
Protocol RTS/CTS (Hardware handshake)

2 INTERFACE SERIAL

even

hardw

ok -

2 SERIELLES INTERF.

Handshake # hardw

**NI**/

**^**Parity

→Handshake

#### **ATC Absolute Temperature Calibration, 3-point calibration** 9.7.

ATC is used to compensate for a temperature difference between the circulator and a defined measuring point in the bath tank that may develop for physical reasons.

ATC

internal					
2 ATCalibration					
⇒Sensor	intern				
₩Status	no				
Type	3-point				
TmpValue1	: xx.xx°C				
CalValue1	: xx.xx°C				
TmpValue2	: xx.xx°C				
CalValue2	: xx.xx°C				
TmpValue3	: xx.xx°C				
CalValue3	: xx.xx°C				

2 ATCalibration				
⇒Sensor		exterr		
₩Status		no		
Type		3-point		
TmpValuei	H	××.××°C		
CalValuei	H	××.××°C		
TmpValue2	H	××.××°C		
CalValue2	H	xx.xx°C		


external


CalValue3 : xx.xx°C

TmpValue3 : xx.xx°C

### Example:

1-point calibration

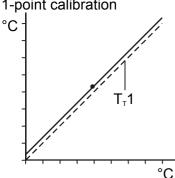

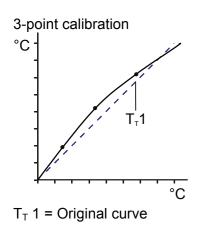


#### Principle:

M

For ATC calibration, in steady state the bath temperature at the location of the temperature sensor (CT) is determined at the respective adjusted working temperature. This value is then set on the circulator in the menu >ATCalibration< under menu item > CTEMP X <.

B (INT)

This can be a 1-point, 2-point or 3-point calibration.

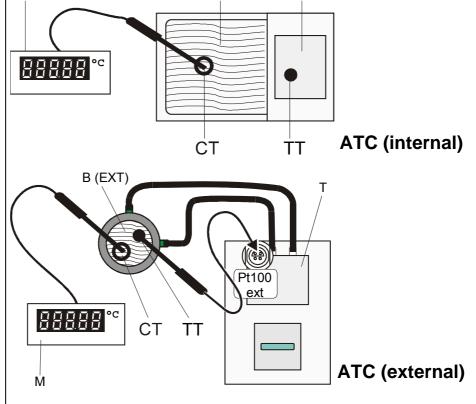


M = Temperature measuring instrument with temperature sensor


B = Bath tank (INTernal or EXTernal)

T = circulator

CT = Temperature on measuring point

TT = Temperature on circulator

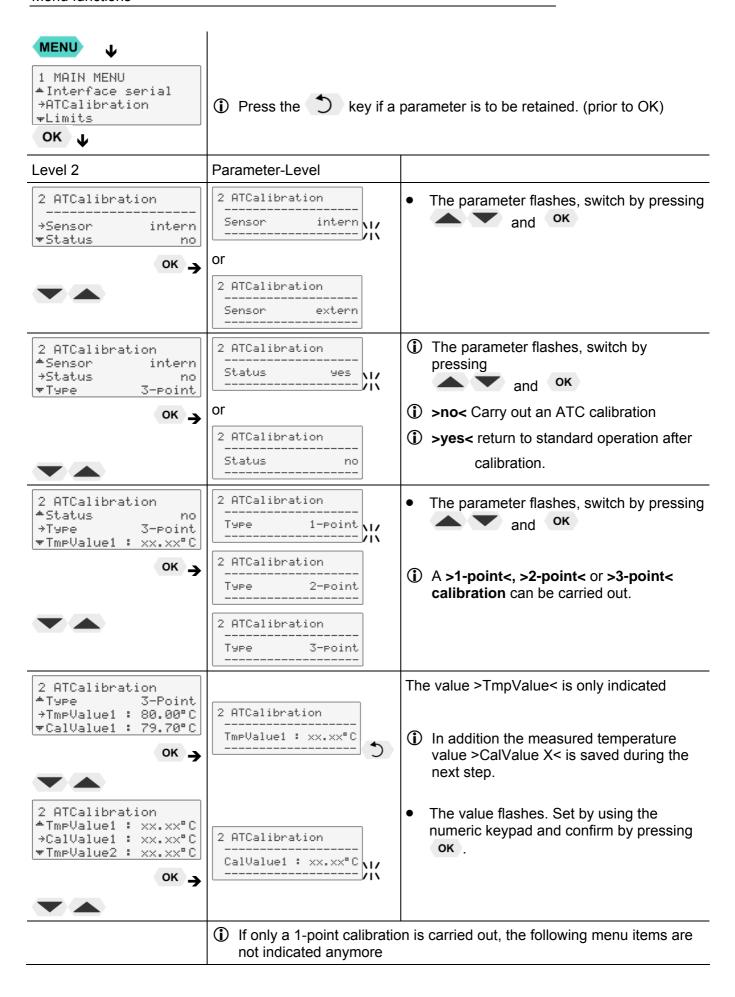

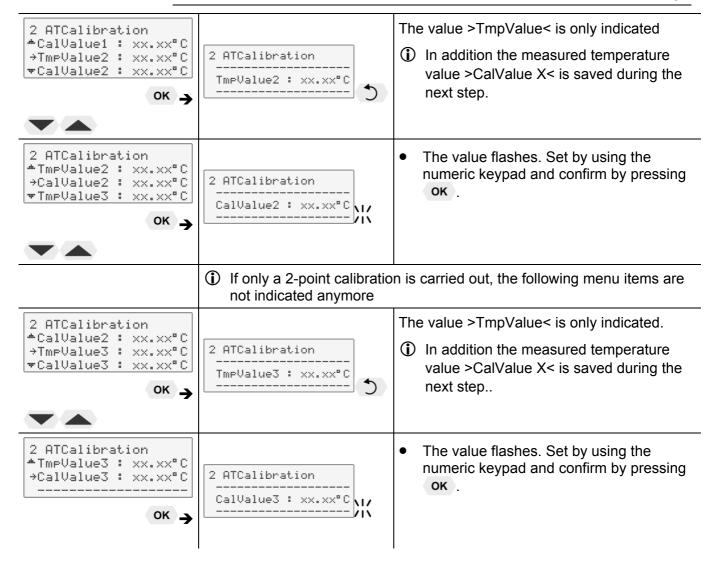

#### 9.7.1. ATC Fühler - intern / extern



In the first submenu the ATC function is set for the >intern< internal or the >extern< external temperature sensor.

Calibration can be carried out for the internal temperature sensor and for the external temperature sensor connected to the socket "ext. Pt100". The circulator is able to save both parameter sets. However only the one which has been set under menu item > sensor < is displayed.

#### 9.7.2. ATC Status - yes / no



In the second submenu the ATC function for the temperature sensor selected above is activated >yes< or deactivated >no<.

>yes< (factory setting) The controller of the circulator uses the original curve of the temperature sensor or the new curve measured during the ATC calibration.


Important: Set to >NO< during the calibration process

>no< An ATC calibration is to be carried out.

**Important:** Set to **>yes<** after calibration.

(i) In the > Status < >yes< the ATC calibration always affects the current working temperature; also the one set via interface.

#### 9.7.3. Type 1 -/ 2 -/ 3 point

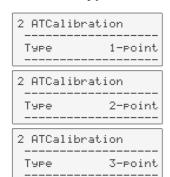


A >1-point<, >2-point< or >3-point< calibration can be carried out.

First geometrically define the location for calibration (measuring point CT), then determine the temperature values of the calibration points. The type of calibrations also determines the number of the following pairs of values indicated on the LCD DIALOG-DISPLAY.

#### Pairs of values:

**TmpValue X:** Circulator temperature 1 or 2 or 3 (actual value TT)

The actual temperature of the bath is simultaneously saved with the "calibration value" >CalValue< and can be indicated for control purposes (value does not flash).

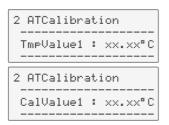

CalValue X: Calibration temperature 1 or 2 or 3 (actual value CT)
The "calibration value" is determined with a temperature measuring device and saved under menu item >CalValue<.


(value flashes (1))

#### 9.7.4. Example: 3-point calibration for internal control

In the temperature range from 80 °C to 160 °C the calibration curve of the temperature sensor (TT) is to be adjusted to the actual temperatures at measuring point (CT).

#### 1. Set circulator to internal control:

The type of control can be set only in the –OFF- mode.

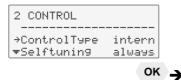




80.00°C

120.00 °C

160.00 °C

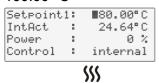

2. Set working temperature setpoint

See standard display line 1

• Example 80.00 °C

Press the keys 8 and 0 and confirm by pressing

The bath is heated up.
 Wait for approx. 5 minutes until the temperature is constant.

Pt100

#### 3. Reading of temperature measuring device

Read the value of measuring point CT on the device and enter under menu item > CalValue X< by using the keypad.


- ÷ CalValue 1< (79.70 °C)
- → CalValue 3< (159.30 °C)</p>




MENU

1 MAIN MENU

#### 4. Calibration

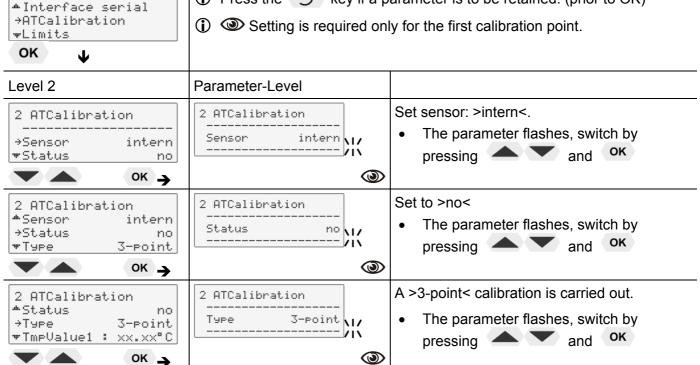



#### Menu functions

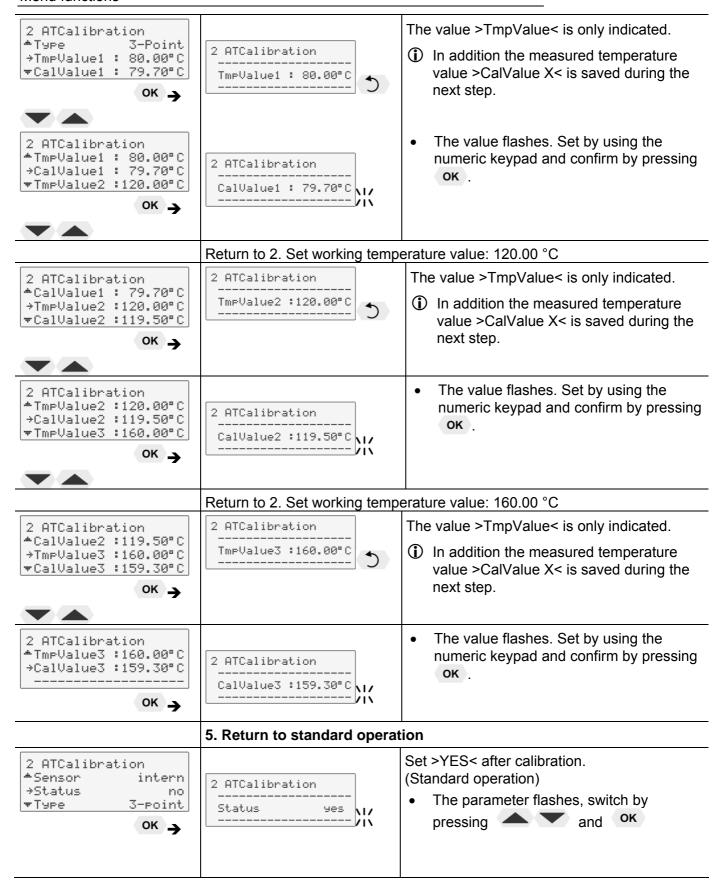


#### 9.8. Limits

The >Limits< menu allows the minimum and maximum values to be set for all important setting ranges and capacity variables.

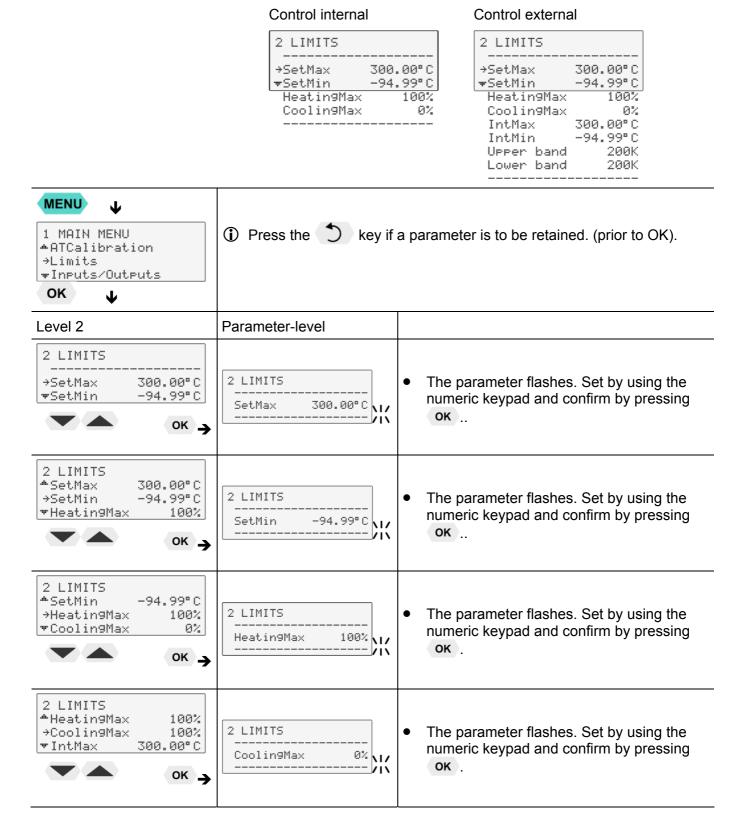

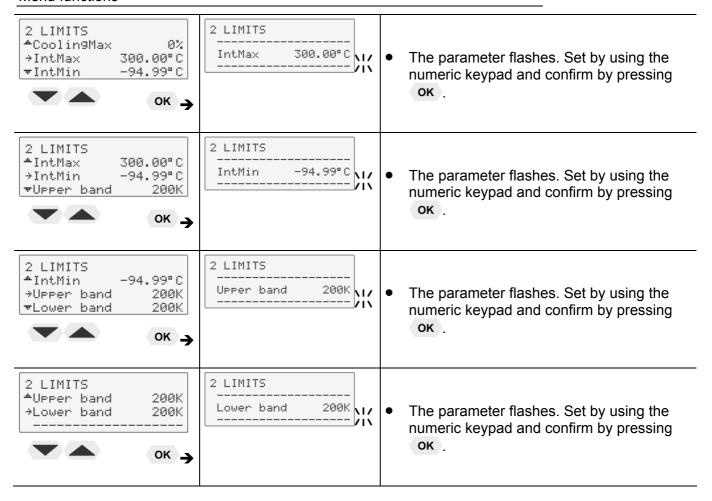

#### 9.8.1. Limits for internal control and external control

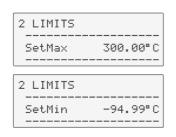

**Set Max / Min–** Maximum and minimum setpoint Restriction of the adjustable temperature range.


The limitation of the operating temperature range effects the temperature setting in the menu with the key .

Only setting of working temperatures which lie within the determined limits is possible

Existing settings for Setpoint 1, -2, -3, as well as those for >OverTemp< and > SubTemp <, are automatically deferred into the limit range.

Setting range: -94,90 °C ... +300,0 °C

### Set maximum heating / cooling

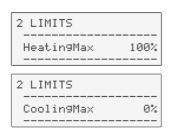


The heating and cooling capacity of the unit are adjustable. 100 % corresponds to the technical specification of the equipment.

#### Setting range:


**Heating Max** – 0 to 100 % in 1 % steps **Cooling Max** – 0 to 100 % in 1 % steps

#### 9.8.2. Limits for external control



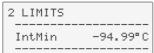



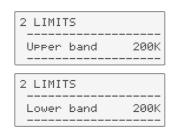

#### **INTERN MAX / MIN**

Restriction for the temperature range of the internal bath.

Setting range: -94,9 °C ... +300,0 °C

The limits IntMax and IntMin are only active in external control. IntMax and IntMin determine fixed limits for the temperature within the internal bath. The temperature controller cannot exceed these limits even if it would be necessary for achieving the temperature in an external system. Therefore it is possible that the external setpoint cannot be achieved.

Sense of limit setting:

- ☑ Protects the bath fluid from overheating.
- ✓ Prevents an undesired alarm shutdown by the excess temperature protection >ALARM CODE 14<.</p>
  Set the value of > Int Max at least 5 °C below the value of >SafeTemp<.</p>
- ☑ Protects the pump motor from high viscosity of the bath fluid at low temperatures.
- ☑ For refrigerated circulators. Freezing protection when using water as bath fluid.


#### Upper band: and Lower band:

The band limitation is active during external control. Varied, practiceoriented setting are feasible for heat-up and cool-down phases.

Setting range: 0 °C ... 200 °C

**Upper band:** and **Lower band** allow for the limitation of the difference between the temperatures in the internal bath and the external system to any maximum value for the heat-up and cool-down phase.


During the heat-up phase this difference value is always added to the actual external temperature. During the cool-down phase the difference value is subtracted.

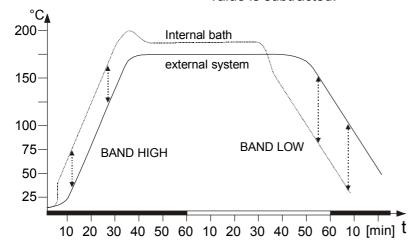

Sense of a band limitation:

- ☑ Protection of objects and samples by gentle temperature control
- Protection of e.g. glass reactors from thermal shock.


# 9.9. Analog inputs/outputs

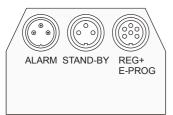


in order to use the analog inputs and outputs, the circulator must be equipped with the electronic module, which is available as option.

Order No. 8 900 100 Electronic module

This submenu enables the input and output values to be set for the programmer input and the temperature recorder outputs of the REG+E-PROG socket.

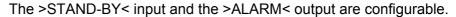

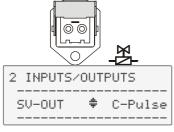

← If the electronic module is not installed, this menu can be used only to switch the control of the solenoid valve (SV).

2 INPUTS/OUTPUTS			
⇒SV-OUT	C-Pulse		
♥Chan1	ActInt		
Chan2	Power		
Chan3	Setpoint		
EPROG	Setroint		
Ext.StB9	inactiv		
AlarmOut	StdBy		

SV-Output Control connector (14) 230 V max. 1.25 A

Channel 1 voltage output for recorder (V)

Channel 2 voltage output for recorder (V))


Channel 3 current output for recorder (mA)


EPROG external programmer input


External standby Standby input


(for external emergency switch-off)

Alarm output Output for external alarm signal

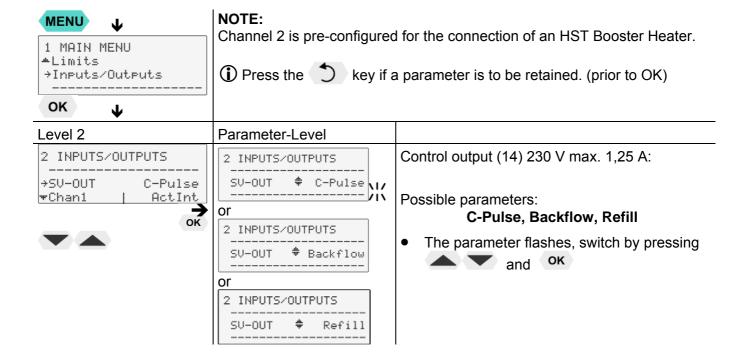

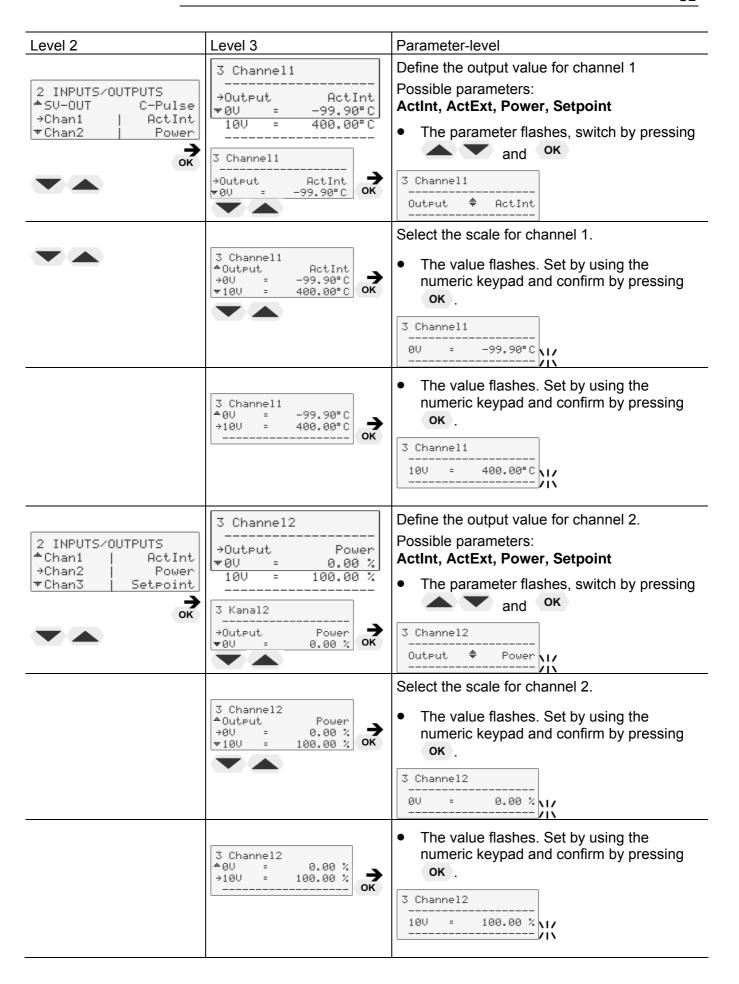

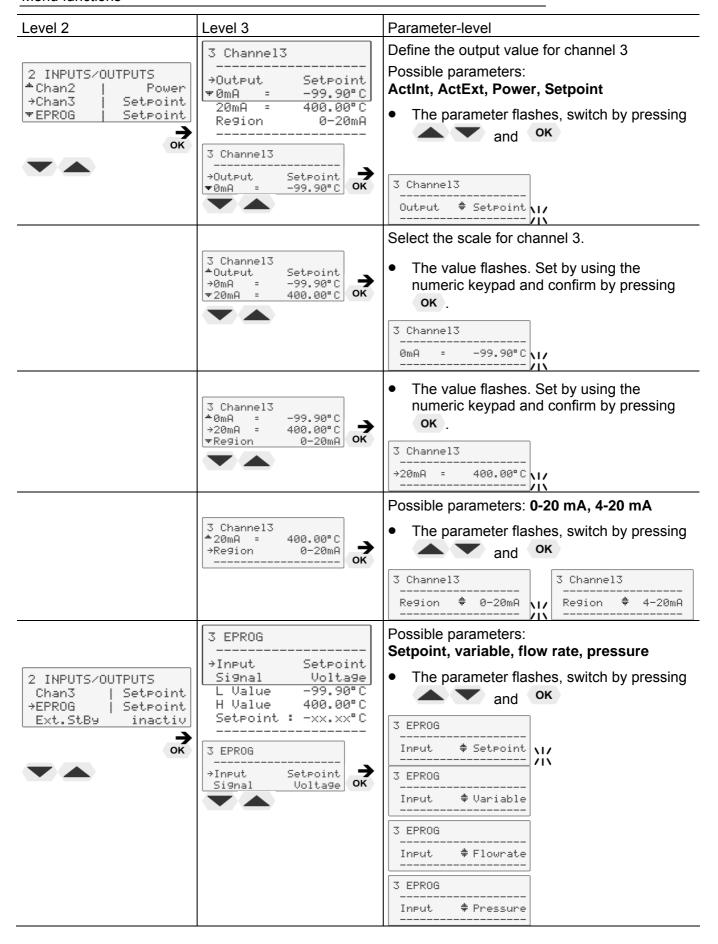

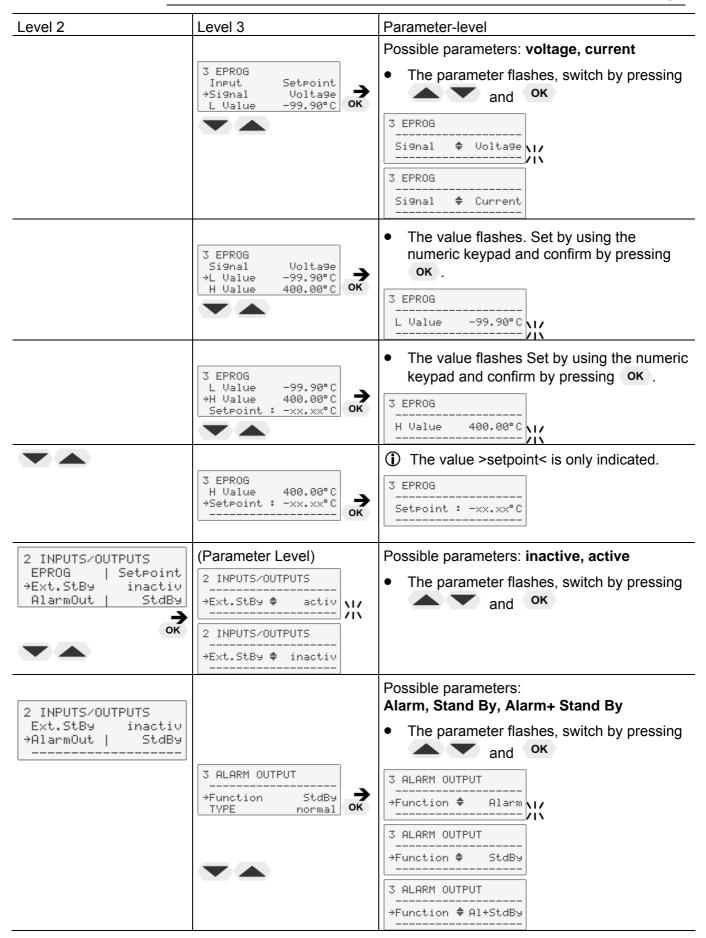

Level 2	Level 3	Parameter-level	
	3 ALARM OUTPUT Function StdB9 →TYPE normal	Possible parameters: <b>normal, inverse</b> • The parameter flashes, switch by pressing  and OK	

#### 9.9.1. Control connector



## Control connector (14):

SV-Out >Backflow< or >C-Pulse< or >Refill<

The control output (14) can be adapted to solenoid valves for different tasks.


>Backflow< In the OFF condition, the control output (14) is not powered.

See - Backflow safety device- on page 29

- >C-Pulse< The control output (14) can be used to emit a cooling pulse. For the pulsing period, the output is powered. (only used with Heating Circulators)
- >Refill< Control of a Automatic Refill Device (Accessories)

#### 9.9.2. Outputs of the connector - REG+E-PROG

#### **Outputs of the connector**

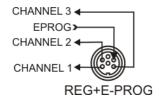

1. First define the desired output value for channels 1 to 3:

ActInt internal actual temperature value (bath temperature)
ActExt external actual temperature value (external sensor)

**Power** periodic or intermittent heating or cooling

**Setpoint** active setpoint temperature

(setpoint 1, 2, 3/integr. programmer/external programmer)

#### 2. Then select the scale for channels 1 to 3:

## Voltage outputs channels 1 and 2


Assign the voltage values of 0 V to the lowest and 10 V to the highest necessary temperature or power rating required as an output value (°C/%).

## **Current output channel 3**

Examples:

Assign the current values 0 mA or 4 mA to the lowest and 20 mA to the highest temperature or power rating required as an output value (°C/%).

3. The current output (channel 3) offers 2 ranges for selection: 0 to 20 mA or 4 to 20 mA

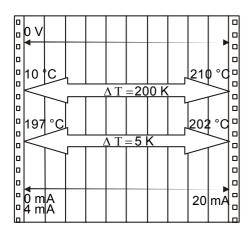

3 Channel1 \*Output ActInt >0V = 10.00°C ▼10V = 210.00°C

lowest temperature value: 10 °C highest temperature value 210 °C Fig. shows 200 °C scaled to paper width

slope: 50 mV/°C

lowest temperature value: 197 °C highest temperature value: 202 °C Fig. shows 5 °C scaled to paper width

slope: 2000 mV/°C

#### 9.9.3. Input of the connector - REG+E-PROG

## Menu >Configuration<

2 CONFIGURATION

\*Remote off

\*Setroint ext Erro9

1. \*Autostart off

2 CONFIGURATION

\*Off-Mode p. off

ActVar Eprog

2. \*Time/Date

#### Menu >Inputs/Outputs<

2 INPUTS/OUTPUTS
Chan3 | Setpoint
→EPROG | Flowrate
3. Ext.StB9 inactiv

2 CONFIGURATION

→Remote off

Setpoint ext Epro9
Autostart off
Off-Mode p. off
ActVar Epro9
Time/Date

#### **E-PROG - Input**

Setting is necessary if

- the Setpoint is to be set via an external voltage or current source or programmer
  - For this, in the menu >Configuration<, first set the menu item >Setpoint< to >Eprog<.
- 2. the heater variable should be controlled via an external control pulse.

For this, in the menu >Configuration<, set the menu item > ActVar.< to >Eprog<.

- 3. the signal of an external flow meter should be registered.
- The E-Prog input can only be used either under menu item >Setpoint< or under menu item > Actvar <.

  If the input is occupied neither by >Setpoint< or > ActVar <, the signal of a flow meter can be connected.
- i First set menu >Configuration<, then menu >Inputs/Outputs<.
- Connect the external voltage or current source or programmer to the REG+E-PROG socket (12) of the circulator.

#### Selecting the signal:

The programmer (E-PROG) input of the circulator can be matched to the output signal of the external voltage or current source.

Voltage voltage input Current current input

#### Example:

3 EPROG	
→Input Si9nal	Setpoint Current
L Value H Value	0.00°C
Setpoint :	50.00°C

"L Value" - Setting the LOW value: (See below  $\bigcirc$ )

First adjust and set the lowest voltage or current on the external voltage or current source (e.g., 0 V or 0 mA). Then after approx. 30 seconds, enter the corresponding temperature value (e.g., 20.00°C).on the circulator by pressing the appropriate keys on the keypad and press **OK** to set.

"H Value" - Setting the HIGH value: (See below  $\bigcirc$ )

First adjust and set the highest voltage or current on the external voltage or current source (e.g., 10 V or 20 mA). Then after approx. 30 seconds, enter the corresponding temperature value (e.g., 300 °C).on the circulator by pressing the

# Example:

(i) Set the external voltage or current source output for the equivalent of a 50 °C temperature setpoint.


appropriate keys on the keypad and press **OK** to set.

The value adjusted and set on the external programmer is displayed in line 4 of the LCD DIALOG-DISPLAY for inspection.

Setpoint1	ŧ	50.00°C
IntAct	i	24.64°C
Power	i	0 %
Control	ŧ	intern

After returning the LCD display to standard display by pressing the key , the temperature value adjusted and set on the external voltage or current source is displayed in line 1 (Example: Eprog 50.00 °C).

This EPROG input enables the use of different voltage and current values as program parameters.

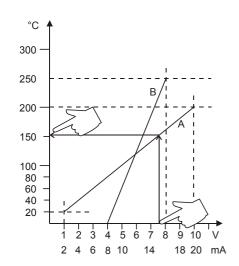

- "L Value" Setting the Low value: (See below 
   )
   1) Adjust and set the lowest desired value on the voltage or current source (Example A: 1 V).

   Wait approximately 30 seconds.
  - 2) Assign a lower temperature threshold value to this adjusted voltage/current value by pressing the appropriate keys on the keypad of the instrument (Example A: 20 °C) and set by pressing **OK**
- "H Value" Setting the High value: (See below 2)
  1) Adjust and set the highest desired value on the voltage or current source (Example A: 10 V).
  Wait approximately 30 seconds.
  2) Assign an upper temperature threshold value to this adjusted voltage/current value by pressing the appropriate keys on the keypad of the instrument
- (i) Example B in the diagram illustrates that the end point values are freely selectable (e.g., 8 mA and 16 mA).

(Example A: 200 °C) and set by pressing OK .

# 3 EPROG →Input Setpoint Signal Voltage L Value 20.00°C H Value 000.00°C Setpoint : 152.00°C

Setpoint1: 152.00°C IntAct : 50.64°C Power : 100 % Control : intern

# Example out of diagram A:

Adjusting the voltage source for an output of 7.6 V!

Line 5 of the LCD DIALOG-DISPLAY shows the externally set setpoint value. The instrument calculates this value from the slope of the two specified end points (in example A, 7.6 V corresponds to an external setpoint temperature of 152.0 °C).

After returning the LCD display to standard display by pressing , this value is displayed in line 1

(Example: EPROG 152.00 °C).



#### Notice:

If this adjustment is not correctly performed at two different points, the setpoint setting will be incorrect.



## Important:

The usable temperature range between > L Value < and > H Value < is limited to the configured working temperature range of the circulator or unit combination. For the working temperature range, see technical specifications.

#### 9.9.4. External Stand-by input

2 INPUTS/OUTPUTS EPROG | Setpoint →Ext.StBy inactiv AlarmOut | StdBy


#### Ext. StBy (External STAND-BY input)

Stand-By input for external switch-off.

#### Possible parameters:

inactive - Stand-By input is ignored

active - Stand-By input is active

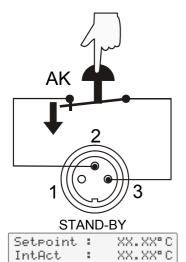

#### Activate the standby input:

- 1. Under menu item >Ext. StBy<, set the parameter to >active<.
- 2. Connect an external contact "AK" (e.g., for external switch-off) or an alarm contact of the superordinated system.

If the connection between pin 2 and pin 3 is interrupted by opening the contact "AK", a complete shutdown of the circulating pump and heater is effected, and the unit enters the condition "**E OFF**".

As long as the contact remains open, line 4 of the LCD DIALOG-DISPLAY flashes and displays the message "Stand-By".

If the contact is reclosed, the instrument returns to the standby state and " **E OFF** " is displayed. Press **OK** to start.



Stand-By


Power

0 %

# (i) Additional tips for using the STAND-BY input:

The standby function can be used in conjunction with the AUTOSTART feature.

- 1. If the Autostart function is NOT turned ON, the standby input is used as described above.
- 2. If the Autostart function is enabled, the instrument will revert back to the original method of entering the setpoint (i.e., keypad, RS232, analog input, etc.).

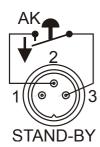

Set : XX.XX°C IntAct: XX.XX°C ProgNr.:2 Step: 2 \*\*\*\*\* StandBy \*\*\*\*\*\* Entering the setpoint with the keypad. e.g.,

As described above, a complete shutdown is accompanied by the "Stand-By" display and the OFF state. The programmable controller starts again when the contact is reclosed. The temperature of the bath fluid changed during the Stand-By state.

#### Entering the setpoint with the programmer.


The display "StandBy" appears. The setpoint value and the time are both held at the current value. The temperature of the bath fluid will be held constant at this temperature. The programmer continues once the contact is reclosed.

 $\Lambda$ 

Notice: this is not an actual emergency switch-off.

#### 9.9.5. Alarm-output



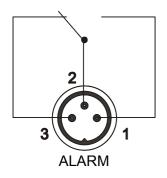

#### Alarm output (10)


(for external alarm signal)

This socket is a potential-free change-over contact. With the adjustments in the menu >Inputs/Outputs<, all operating conditions can be signaled without having to change the pin assignments.



Meaning of the terms under menu item >Function<:

#### The circulator is in condition

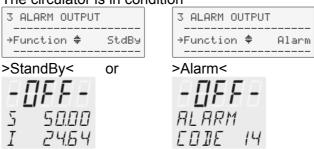

For >Type: normal<, pins 2 and 3 are connected in any case according to the selected >Function<.

For >Type: inverse<, pins 2 and 1 are connected in any case according to the selected >Function<.

Switching capacity max. 30 W/40 VA Switching voltage max. 125 V~/– Switching current max. 1 A

# 10. Troubleshooting guide / Error messages



## Alarm with complete shutdown:

If one of the following failures occur a complete, all-pole shutdown of the heater and circulating pump is effected.



"🃤" lights up and a continuous signal sounds.

The code for the cause of alarm is indicated on the VFD COMFORT-DISPLAY.

## Alarm without shutdown:

The code for the cause of alarm is indicated on the VFD COMFORT-DISPLAY. The warning signal sounds in regular intervals. The messages appear every 10 seconds.



- Press the key OK to stop the signal
- Press ok to indicate the help text in the LCD DIALOG-DISPLAY when alarm messages appear.

# ALARM COJE O I

The circulator is operated without bath fluid, or the liquid level is insufficient. Replenish the bath tank with the bath fluid.

Low liquid level or float brocken. Check float and fill medium! Tube breakage has occurred (insufficient filling level due to excessive bath fluid pumped out). Replace the tubing and replenish the bath tank with the bath fluid.

The float is defect (e. g., because damaged in transit). Repair by authorized JULABO service personnel.

# ALARM COJE OZ

During the self-test after switch-on a short-circuit is registered between pin 2 and pin 4 of the control cable or the control cable is interrupted during operation.

Control cable of the cooling machine is defective.

Ckeck connection!

Reconnect the cable or eliminate the short-circuit.

# WARNING COUE O3

## ALARM COJE O3

 Excess temperature warning or Excess temperature alarm

Actual temperature obove high temperature value. Check limits!

Warn-Type: Set to >Warning< or >Alarm<

WARNING EDJE DY

ALARM COJE OY  Low temperature warning or Low temperature alarm.

Actual temperature below low temperature value. Check limits!

Warn-Type: Set to >Warning< or >Alarm<

# ALARM COJE OS

The cable of the working temperature sensor is disconnected or short-circuited.

Internal working sensor is shorted or interrupted. Call service!

ALARM COUL OB  Sensor difference b. working- and safety- temp. Check pump stage and viscosity!	Defect of the working or excess temperature sensor.  The working temperature and excess temperature sensors report a temperature difference of more than 35 K.
ALARM EDJE D7  Internal hardware	Other errors (I <sup>2</sup> C-BUS errors)
error.  Call service!	
RLARM EDIE 12	Error in A/D converter
Error in A/D converter of the measuring system. Call service!	
ALARM	Excess temperature sensor is defective.
EDIE 14	The safety temperature is below the working temperature setpoint. Set the
Temperature exceeds the adjusted safety-temperature. Check adjustment!	safety temperature to a higher value.
ALARM EDJE 15	External control selected, but external Pt100 sensor is not connected or is defective.
External sensor is not connected or brocken. Check the external sensor!	
WARNING COJE 20	Cooling of the condenser is impaired. Clean the air-cooled condenser.  For a water-cooled condenser, check the flow rate and temperature of the cooling water.
WARNING EDJE 21	The first-stage compressor does not work.  After a short cooling interval, the compressor motor will be automatically reconnected and the message "CODE 21" will no longer appear.
WARNING	The second-stage compressor does not work.
COJE 22	Cooling compressor overload protection  The motor of the cooling compressor is equipped with an overload protector, which will be activated in case of excessive temperature in the capsule or excessive current draw.
	Causes of motor shut-off: - poor air circulation - short distance to walls - dirt accumulated on condenser
	<ul><li>high ambient temperature</li><li>compressor switched off and on at short intervals</li></ul>
WARNING COUE 23	Excess temperature in the first-stage compressor.
WARNING EDJE 24	Excess temperature in the second-stage compressor.

*WARNING* EDDE 25 Short-circuit in the control cable for the cooling machine during self-test.

RL RRM EDJE 33

Cable of the excess temperature sensor disconnected or short-circuited.

Internal safetytemperature sensor Call service!

shorted/interrupted.

RLRRMEDIE38

ГППЕ ЧП

Ext. Pt100 sensor input without signal, but setpoint programming set to external Pt100.

Ext. sensor input without signal but setpoint programming set to ext. Pt100!

*WRRNING* 

The early warning system for low level signals a critical fluid level. Replenish the bath tank with bath fluid.



After eliminating the malfunction, turn the mains power switch off and on again to cancel the alarm state.

If the unit cannot be returned to operation, contact an authorized JULABO service center.

CONFIGURATION **ERROR** CONFIRM BY PRESSING : <OK> Special message "CONFIGURATION ERROR"

The configuration of the circulator does not correspond to its present use.

Press **OK** to automatically perform a one-time modification of the configuration.

Then contact an authorized JULABO service center.

#### Disturbances that are not indicated.

The electronic pump motor is protected against overload by an electronic current limiter. If the viscosity of the bath fluid is or becomes too high, the motor stops running.



#### Fuses:

The mains fuses on the rear of the unit may easily be replaced as shown at

Circulator: Fine fuses (15)- T 1.25 A, 250 V~, D 5 x 20 mm Circulator: Fine fuses (16a)- T 16 A, 250 V~, D5 x 20 mm



#### Warning:

Before exchanging the fuses, turn off the mains power switch and disconnect the power plug from the mains socket!

Only use fine fuses with a nominal value as specified.

#### Example:

Manufacturer	Supplier	Туре	Order-No.
Schurter	Schurter	G-fuse link SPT T16A 5x20mm	No. 0001.2516
Wickmann	Wickmann	G-fuse link T1,25A 5x20 mm	No. 19195

## 11. Electrical connections



## Notice:


Use shielded cables only.

The shield of the connecting cable is electrically connected to the plug housing.

The unit ensures safe operation if connecting cables with a maximum length of 3 m are used. The use of longer cables does not affect proper performance of the unit, however external interferences may have a negative impact on safe operation (e.g. cellular phones).




#### Socket for external Pt100 sensor

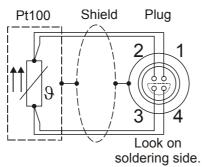

# Pin assignment:

Signal
l+
U+
U-
I-

The shield of the connecting cable is electrically connected to the plug housing and the sensor tube.

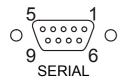

#### RS232/RS485 serial interface

This port can be used to connect a computer with an RS232 or RS485 cable for remote control of the circulator.

#### Pin assignments RS232:

Pin 2	RxD	Receive Data
Pin 3	TxD	Transmit Data
Pin 5	0 V	Signal GND
Pin 7	RTS	Request to send
Pin 8	CTS	Clear to send

Pin 1; 4; 6, 9 Reserved - do not use!

#### Pin assignments RS485:

Pin 3 B
Pin 5 0 V Signal GND
Pin 6 +5 V (max. 50 mA)
Pin 8 A

Pin 1; 2; 4; 7; 9 Reserved - do not use!



## Notice: RS485 serial interface

The max. load current between pin 5 and pin 6 is 50 mA.


## RS232 interface cable

Circulator (9-pol)		PC (9-pol)
Pin 2 RxD	$\Leftrightarrow$	Pin 3 TxD
Pin 3 TxD	$\Leftrightarrow$	Pin 2 RxD
Pin 5 GND	$\Leftrightarrow$	Pin 5 GND
Pin 7 RTS	$\Leftrightarrow$	Pin 8 CTS
Pin 8 CTS	$\Leftrightarrow$	Pin 7 RTS

Accessories:	Order No.	Description
	8 980 073	RS232 interface cable 9-pol./9-pol. , 2,5 m
	8 900 110	USB interface adapter cable




The connector may only be used for control of a JULABO refrigerated circulator or JULABO MVS Solenoid valve controller for cooling water.

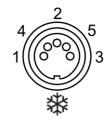

## Pin assignment:

Pin	Signal (only with attached JULABO equipment)
1	+24 V (I max. current 25 mA)
2	0 V
3	Alarm relay
4	Reserved - do not use!
5	Cooling pulse

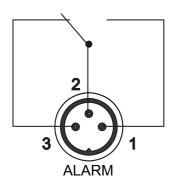


## **Alarm output**

(for external alarm signal)

This potential-free change-over contact is activated in case of an alarm when pins 2 and 3 are connected.

Switching capacity	max.	30 W / 40 VA
Switching voltage	max.	125 V~/−
Switching current	max.	1 A

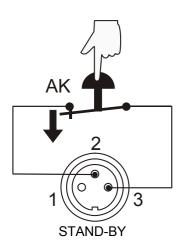

## Stand-by input

(for external switch-off)

Pin assignment:	<u>Pin</u>	Signal
	1	not connected
	2	5 V / DC
	3	0 V





<u>Pin</u>		Signal		
1 Voltage output	Channel 1	0 10 V		
2 Voltage output	Channel 2	0 10 V		
3 GND for outputs	0 V			
4 Programmer input	EPROG	0 to 10 V / 0 to 20 mA		
5 Current output	Channel 3	0 to 20 mA / 4 to 20 mA		
6 GND for Progammer 0 V				





# Control connectors (13, 14)

The control output (13) is not powered in the OFF condition.


An external pump motor can be connected.

Output voltage: 230 V~ / max. 1.25 A

The control output (14) can be adapted to solenoid valves for different tasks.

## 12. Remote control

# 12.1. Setup for remote control

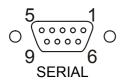

1. Check the interface parameters for both interfaces (on the circulator and PC) and make sure they match.

In the >Interface< menu, set the menu item >Type< to >RS232< or >RS485<.



- 2. In the >Configuration< menu, set the menu item >Setpoint< to >RS232< or >RS485<.
- 3. Connect both units with an interface cable.



Like all parameters that can be entered through the keypad, interface parameters are stored in the memory even after the circulator is turned off.

# 12.2. Communication with a PC or a superordinated data system

If the circulator is put into remote control mode via the configuration level, the VFD COMFORT-DISPLAY will read "R -OFF-" = REMOTE STOP. The circulator is now operated via the computer.

In general, the computer (master) sends commands to the circulator (slave). The circulator sends data (including error messages) only when the computer sends a query.



In remote control mode, the start command and all values to be set must be resent by the PC via the interface in case of a power interruption.

AUTOSTART is not possible.

A transfer sequence consists of:

- address (RS485 interface only)
- command
- space (⇔; Hex: 20)
- parameter (decimal separation with a period)
- end of file (↓; Hex: 0D)
- The response (data string) after an in command is always followed by a line feed (LF, Hex: 0A).

The commands are divided into in or out commands.

in commands: retrieve parametersout commands: set parameters

## Important times for a command transmission:



To ensure secure data transfer, the time gap between two commands should be at least 250 ms.

The circulator automatically responds to an **in** command with a data string followed by an LF (Line Feed). The next command should be sent only after 10 ms.



The **out** commands are valid only in remote control mode.

Examples:

When the RS485 interface is used, the three-digit instrument address precedes each command.

(example: address Ad32 = A032)

Command to set the working temperature >Setpoint1< to 55.5 °C

out\_sp\_00 ⇔ 55.5↓ A032\_out\_sp\_00 ⇔ 55.5↓

Command to retrieve the working temperature >Setpoint1<

in\_sp\_00↓ A032\_in\_sp\_00↓

Response from the circulator:

55.5↓ LF A032\_55.5↓ LF



## Notice: out commands

Temperature values can be displayed on the circulator in °C or °F.

The settings sent via the interface must also be made in the selected unit according to this setting.

## 12.3. List of commands

out commands: Setting parameters or temperature values.

Command	Parameter	Response of circulator
out_mode_01	0	Use working temperature >Setpoint1<
out_mode_01	1	Use working temperature >Setpoint2<
out_mode_01	2	Use working temperature >Setpoint3<
out_mode_02	0	Selftuning "off". Temperature control using the stored parameters.
out_mode_02	1	Selftuning "once" Single selftuning of the controlled system after the next start.
out_mode_02	2	Selftuning "always"  Continual selftuning of the controlled system whenever a new setpoint is to be reached.
out_mode_03	0	Set external programmer input to voltage.  Voltage 0 V 10 V
out_mode_03	1	Set external programmer input to current.  Current 0 mA 20 mA
out_mode_04	0	Temperature control of internal bath.
out_mode_04	1	External control with Pt100 sensor.
out_mode_05	0	Stop the unit = R –OFF

## Remote control

Command	Parameter	Response of circulator	
out_mode_05	1	Start the unit.	
out_mode_08	0	Set the control dynamics - aperiodic	
out_mode_08	1	Set the control dynamics - standard	
out_sp_00	XXX.XX	Set working temperature. "Setpoint 1"	
out_sp_01	XXX.XX	Set working temperature. "Setpoint 2"	
out_sp_02	XXX.XX	Set working temperature. "Setpoint 3"	
out_sp_03	xxx.xx	Set high temperature warning limit "OverTemp"	
out_sp_04	XXX.XX	Set low temperature warning limit "SubTemp"	
out_sp_06	XXX.XX	Set manipulated variable for heater via serial inte -100 +100 [%]	erface
out_sp_07	х	Set the pump pressure stage. (1 4)	
out_par_04	X.X	CoSpeed for external control 0 5.0.	
out_par_06	xxx	Xp control parameter of the internal controller.	0.1 99.9
out_par_07	xxx	Tn control parameter of the internal controller.	3 9999
out_par_08	XXX	Tv control parameter of the internal controller.	0 999
out_par_09	xxx	Xp control parameter of the cascade controller.	0.1 99.9
out_par_10	XXX	Proportional share of the cascade controller.	1 99.9
out_par_11	XXX	Tn control parameter of the cascade controller.	3 9999
out_par_12	XXX	Tv control parameter of the cascade controller.	0 999
out_par_13	XXX	Maximum internal temperature of the cascade co	ntroller.
out_par_14	XXX	Minimum internal temperature of the cascade con	ntroller.
out_par_15	xxx	Band limit (upper) 0 200 °C	
out_par_16	XXX	Band limit (lower) 0 200 °C	
out_hil_00	-xxx	Set the desired maximum cooling power (0% to 1 Note: Enter the value with a preceding negative so This setting is meaningful only for FP cooling maximum.	sign!
out_hil_01	XXX	Set the desired maximum heating power (10% to	100%).

in commands: Asking for parameters or temperature values to be displayed.

Command	Parameter	Response of circulator
version	None	Version number of the software (V X.xx)
status	none	Status message, error message (see page 94)

Command	Parameter	Response of circulator
in_pv_00	none	Current bath temperature.
in_pv_01	none	Heating power being used (%).
in_pv_02	none	Temperature value registered by the external Pt100 sensor.
in_pv_03	none	Temperature value registered by the safety sensor.
in_pv_04	none	Setpoint temperature ("SafeTemp") of the excess temperature protection
in_sp_00	none	Working temperature "Setpoint 1"
in_sp_01	none	Working temperature "Setpoint 2"
in_sp_02	none	Working temperature "Setpoint 3"
in_sp_03	none	High temperature warning limit "OverTemp"
in_sp_04	none	Low temperature warning limit "SubTemp"
in_sp_05	none	Setpoint temperature of the external programmer (REG+E-PROG).
in_sp_06	none	Temperature indication in °C or °F
in_sp_07	none	Adjusted pump stage in the -OFF- condition.     Pump stage corresponding to the effective rotational speed after start.
in_sp_08	none	Value of a flow rate sensor connected to the E-Prog input
in_sp_09	none	Value of a pressure sensor connected to the E-Prog input
in_par_00	none	Temperature difference between working sensor and safety sensor
in_par_01	none	Te - Time constant of the external bath.
in_par_02	none	Si - Internal slope
in_par_03	none	Ti - Time constant of the internal bath.
in_par_04	none	CoSpeed - Band limit (max. difference between the temperatures in the internal bath and external system).
in_par_05	none	Factor pk/ph0: Ratio of max. cooling capacity to max. heating capacity
in_par_06	none	Xp control parameter of the internal controller.
in_par_07	none	Tn control parameter of the internal controller.
in_par_08	none	Tv control parameter of the internal controller.
in_par_09	none	Xp control parameter of the cascade controller.
in_par_10	none	Proportional share of the cascade controller.
in_par_11	none	Tn control parameter of the cascade controller.
in_par_12	none	Tv control parameter of the cascade controller.
in_par_13	none	Adjusted maximum internal temperature of the cascade controller.
in_par_14	none	Adjusted minimum internal temperature of the cascade controller.
in_par_15	none	Band limit (upper)
in_par_16	none	Band limit (lower)

Command	Parameter	Response of circulator
in_mode_01	none	Selected setpoint:
		0 = Setpoint 1
		1 = Setpoint 2
		2 = Setpoint 3
in_mode_02	none	Selftuning type:
		0 = Selftuning "off"
		1 = Selftuning "once"
		2 = Selftuning "always"
in_mode_03	none	Type of external programmer input:
		0 = Voltage 0 V to 10 V
		1 = Current 0 mA to 20 mA
in_mode_04	none	Internal/external temperature control:
		0 = Temperature control with internal sensor.
·		1 = Temperature control with external Pt100 sensor.
in_mode_05	none	Circulator in stop/start condition:
		0 = stop
		1 = start
in_mode_08	none	Adjusted control dynamics
		0 = aperiodic
		1 = standard
in_hil_00	none	Max. cooling power (%).
in_hil_01	none	Max. heating power (%).

# 12.4. Status messages

_	Status messages	Description
-	00 MANUAL STOP	Circulator in "OFF" state.
	01 MANUAL START	Circulator in keypad control mode.
	02 REMOTE STOP	Circulator in "r OFF" state.
	03 REMOTE START	Circulator in remote control mode.

# 12.5. Error messages

Error messages	Description
-01 LOW LEVEL ALARM	Low liquid level alarm
-02 REFRIGERATOR ALARM	Control cable of the refrigerated circulator or MVS solenoid valve controller short-circuited or disconnected.
-03 EXCESS TEMPERATURE WARNING	High temperature warning
-04 LOW TEMPERATURE WARNING	Low temperature warning.

Error messages	Description
-05 WORKING SENSOR ALARM	Working temperature sensor short-circuited or disconnected.
-06 SENSOR DIFFERENCE ALARM	Sensor difference alarm. Working temperature and safety sensors report a temperature difference of more than 35 K.
-07 I <sup>2</sup> C-BUS ERROR	Internal error when reading or writing the I <sup>2</sup> C bus.
-08 INVALID COMMAND	Invalid command.
-09 COMMAND NOT ALLOWED IN CURRENT OPERATING MODE	Invalid command in current operating mode.
-10 VALUE TOO SMALL	Entered value too small.
-11 VALUE TOO LARGE	Entered value too large.
-12 TEMPERATURE MEASUREMENT ALARM	Error in A/D converter.
-13 WARNING : VALUE EXCEEDS TEMPERATURE LIMITS	Value lies outside the adjusted range for the high and low temperature warning limits. But the value is stored.
-14 EXCESS TEMPERATURE PROTECTOR ALARM	Excess temperature protector alarm
-15 EXTERNAL SENSOR ALARM	External control selected, but external Pt100 sensor not connected.
-20 WARNING: CLEAN CONDENSOR OR CHECK COOLING WATER CIRCUIT OF REFRIGERATOR	Cooling of the condenser is impaired. Clean air-cooled condenser. Check the flow rate and cooling water temperature for a water-cooled condenser.
-21 WARNING: COMPRESSOR STAGE 1 DOES NOT WORK	First-stage compressor does not work.
-22 WARNING: COMPRESSOR STAGE 2 DOES NOT WORK	Second-stage compressor does not work.
-23 WARNING: HIGH TEMPERATURE ON COMPRESSOR STAGE 1	Excess temperature at first-stage compressor.
-24 WARNING: HIGH TEMPERATURE ON COMPRESSOR STAGE 2	Excess temperature at second-stage compressor.
-25 REFRIGERATOR WARNING	Error in the cooling machine.
-26 WARNING: STAND-BY PLUG IS MISSING	External standby contact is open.
-30 CONFIGURATION ERROR: CONFIRM BY PRESSING <ok> ON CIRCULATOR</ok>	The configuration of the circulator does not correspond to its present use.  Press OK to automatically perform a one-time modification of the configuration.
-33 SAFETY SENSOR ALARM	Excess temperature sensor short-circuited or disconnected.
-38 EXTERNAL SENSOR SETPOINT PROGRAMMING ALARM	Ext. Pt100 sensor input without signal and setpoint programming set to external Pt100.
-40 NIVEAU LEVEL WARNUNG	Low liquid level warning in the internal reservoir.
·	


# 13. JULABO Service – Online remote diagnosis

JULABO circulators of the HighTech series are equipped with a black box. This box is implemented in the controller and records all significant data for the last 30 minutes.

In case of a failure, this data can be read out from the unit by using special software. This software is available as a **free** download from <a href="www.julabo.de">www.julabo.de</a> \ EasyBlackBox.

Installation is easy and is performed step by step.

Please observe the instructions.

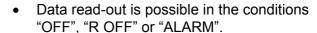


- Connect the circulator to the computer using an interface cable.
- Start the EasyBlackBox program.
   The program asks for the port used (COM1, .....) and the baud rate of the unit.
   You do not have this information on hand?
   Simply try it out!

The program continues to send the request until the correct settings are made.

- Data is read out and shown on the monitor divided into the sections
  - >Einstellungen/Settings<,
  - >Alarmspeicher/Alarms stored<,
  - >Blackbox<
  - ← see example
- After pressing >Speichern/Save<, a text file is created. The program suggests a filename ->C:\model description and barcode no.<.</li>
   Modifications are possible.
- E-mail this file to <u>service@julabo.de</u>, JULABO's service department. JULABO is thus able to provide rapid support.



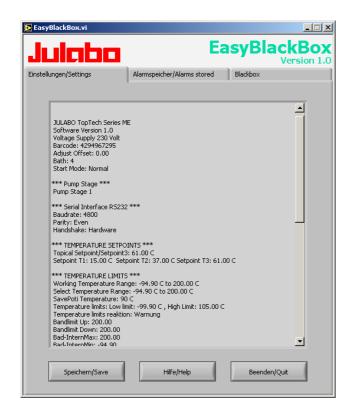


# 14. Installation of electronic module with analog connectors



#### Caution:

Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures, or before moving the unit. Configuration, installation, maintenance, and repairs on the circulator may be performed only by qualified personnel.

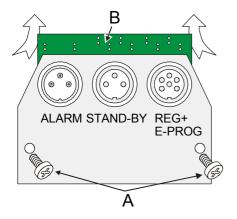


- Turn off the unit with the mains switch and disconnect the power plug.
- Remove the screws (A) and remove the dummy plate.
- Carefully insert the upper edge (B) into the rear panel from below, align the electronic module at the top and bottom, and connect the 15-pin connector by applying light pressure.
- Attach the electronic module using the screws (A).
- The circulator is ready for use. It automatically recognizes the installed electronic module.

# 15. Cleaning / repairing the unit



#### Caution:

- Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit.
- Prevent humidity from entering into the circulator.
- Electrical connections and any other work must be performed by qualified personnel only.

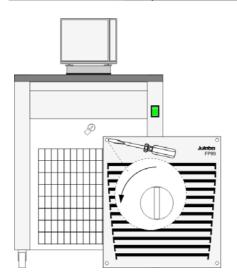

In dusty environments, the refrigerating machine's condenser should be cleaned by vacuuming every 1 to 2 months. The condenser is located behind the venting grid on the front of the unit, which can be taken off by removing the screws.

- Switch off the unit, disconnect mains power cable.
- Remove the venting grid.
- Clean the ribbed condenser with a vacuum cleaner.
- · Replace the venting grid.
- Switch on the unit.

#### Cleaning:

For cleaning the bath tank and the immersed parts of the circulator, use low surface tension water (e.g., soap suds).

Clean the outside of the unit using a wet cloth and low surface tension water.

The circulator is designed for continuous operation under normal conditions. Periodic maintenance is not required.

The tank should be filled only with a bath fluid recommended by JULABO. To avoid contamination, it is essential to change the bath fluid from time to time.

#### Repairs

Before asking for a service technician or returning a JULABO instrument for repair, please contact an authorized JULABO service station.


When returning the unit:

- Clean the unit in order to avoid any harm to the service personnel.
- Attach a short fault description.
- During transport the unit has to stand upright. Mark the packing correspondingly.
- When returning a unit, take care of careful and adequate packing.
- JULABO is not responsible for damages that might occur from insufficient packing.



JULABO reserves the right to carry out technical modifications with repairs for providing improved performance of a unit.

# Anhang: HST Zusatzheizer und HSP Zusatzpumpe

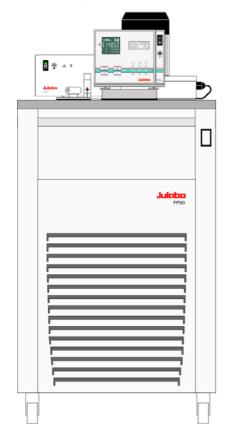

#### **Upgradable with supplementary heater - HST**

For control a HST booster heater a electronic module is required.

Order No. Description

8810012 HST booster heater 6 kW

8900100 Electronic module

Technical specifications on page 101

Upgradable with supplementary pump - HSP

Order No. Description

8810015 HSP booster pump, 30 l/min. –3 bar max.


Technical specifications on page 103



#### Caution:

To ensure full operating safety, the HST booster heater and HSP booster pump must be securely fixed to the refrigerated bath using the screws supplied with the unit!

#### Top view:

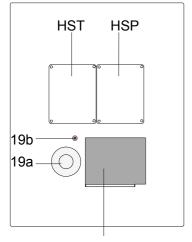

SL

circulator

#### Mounting the optional devices:


On the basic version of the ultra-low refrigerated circulator, the mounting openings for the HSP booster pump and the HST booster heater are closed with stainless steel lids.


- If the optional devices are retrofitted, remove the four screws on each of the stainless steel lids.
- Place the booster pump and/or the booster heater in the respective opening and fix it using the screws supplied with the unit.


#### Connecting the power cable and the control cable:

- For the circulator basic version see page 22 and page 32
- For the HST booster heater see page 100
- For the HSP booster pump see page 102

#### 16. **HST** booster heater

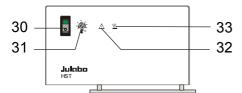

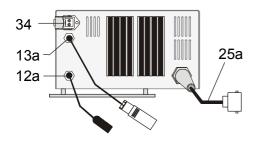

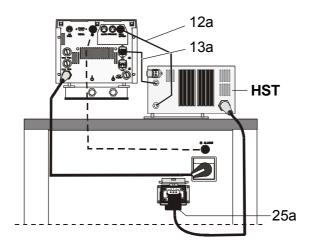

Factory setting

12a

31

Connecting cable to the REG+EPROG socket (12) of the circulator.

(i) Heating pulse transmission of the HST booster heater.

#### NOTE:

3 Kanal2 ⇒Aus9an9 Leistun9 0.00 % 10U 100.00 %

Channel 2 is pre-configured for connecting a HST booster heater (see page 74).

If the HST booster heater is retrofitted, please check the setting for Channel 2 because it might have been configured for a different use in the meantime.

Connecting cable to the control connector (13) of the circulator. 13a

(i) Power supply for the control electronics of the HST booster heater.

Connecting cable for the HST booster heater. 25a

> Built-in mains outlet for connection of HST booster heater Not on -FPW91-SL - 230 V/3PPE/60 Hz (see page 17)

i Power supply for the heater.

Mains power switch, illuminated 30

i Turn the power supply (and the HST booster heater) on and off.

Adjustable excess temperature protection on the HST (see page 101)



(i) A safety device independent from the circulator. If the actual bath temperature exceeds the adjusted safety temperature value at the safety potentiometer, the supplementary heater is switched off permanently.

There is an optical and audible alarm indication with a permanent signal sound. The error is not reported to the circulator!

32	Λ	Indicator light: Alarm indication (see 31)
33	<u>sss</u>	Indicator light: Heating  i Blinks in time with the heating pulse/booster heater, resp.
34		Control connector 230 V / max 1 25 A for HSP booster pump

connector 230 V / max. 1.25 A for HSP booster pump

(i) If the ultra-low refrigerated circulator is provided with HST booster heater and HSP booster pump, this connector replaces the unusable control connector (14) on the rear side of the circulator. The function is independent from the switch position of the mains power switch (30) at the HST.

# 16.1. Excess temperature protection on the HST booster heater



## Warning:



The excess temperature protection >SafeTemp< should be set at least 25 °C below the fire point of the bath fluid used.

In the event of wrong setting there is a fire hazard!

We disclaim all liability for damage caused by wrong settings!



A safety device not depending on the circulator. If the actual bath temperature exceeds the adjusted safety temperature value at the safety potentiometer, the supplementary heater is switched off permanently. There is an optical and audible alarm indication with a permanent signal sound.

Adjust the safety potentiometer to the value for the high temperature protection (see above) by means of a screwdriver and the scale which is printed on. Example above:  $50\,^{\circ}\text{C}$ 

#### Recommendation::

When the working temperature in the bath is reached, the safety potentiometer should be turned back slowly to the switched-off point by means of a screwdriver. Afterwards use the scale to adjust the value by 10 °C higher. Press the mains power switch of the HST off and on again to cancel the alarm

# Setting range:

0 °C to 220 °C state.

# 16.2. Technical specifications - HST booster heater:

Heater capacity kW 6

Mains power connection V / Hz 230/50-60/2phases

(See chapter "Mains power connection" page 16)

Current draw (at 230 V) A 16 / Phase

Overall dimensions HST (WxDxH) mm 210x200x160 (from assembly level on)


Weight kg 5.5

Safety Installations

Excess temperature protection adjustable from 0 to 220 °C

Alarm indication optical + audible (continuous tone)

# 17. HSP booster pump

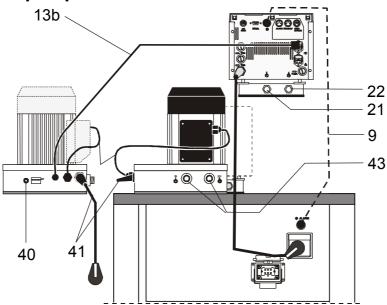

13b Connecting cable to the control connector (13) of the circulator.

Feed (OUT), Return (IN)

**(i)** Power supply for the control electronics of the HSP booster pump.

40	*****	Overflow connector  Only if the upper shaft seal of the pump motor fails, the penetrating bath fluid is let off at this overflow and thus the destruction of the motor is prevented.
41		Mains cable with plug of the HSP booster pump.     Connect the HSP booster pump with mains power cable to the mains socket.
21		Important::
22		If an HSP booster pump is used, the pump connections (21, 22) cannot be used any more They have to be closed tightly with union nuts with inserted blind sieves.



## Important notice:

With the installation of the HSP booster pump, the pump capacity of the ultra-low refrigerated circulator increases.

Pump connector for external temperature application - M16x1

Before starting the HSP booster pump, please check the insert capacity of the employed test tubes.

For this, see chapter tubing on page **Fehler! Textmarke nicht definiert.** in the operating manual.

Cooling capacity: On ultra-low refrigerated circulators, the use of the supplementary pump reduces the cooling capacity by approx. 400 W.

43


# 17.1. Pump connection with HSP booster pump



Caution: Securely attach all tubing to prevent slipping.

#### Important::

If an HSP booster pump is used, the pump connections (21, 22) cannot be used any more They have to be closed tightly with union nuts with inserted blind sieves.

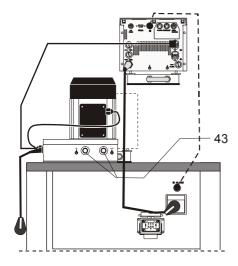

Pump connectors (43) M16x1:

- Unscrew the M 16 x 1 collar nuts on the pump connectors (43) (feed the return to with a 19 mm (3/4") wrench and remove the sealing disks.
- Attach the tubing to the connectors of the external open system and fasten with tube clamps to prevent slipping.

# 17.2. Technical specifications - HSP booster pump

Pressure, max.	Psi / bar at 0 liters	55 / 3.8	43.5 / 3.0
Discharge, max.	lpm at 0 bar	36	30
with tubing connecti	ons mm ID	13.5	13.5
Mains power connection ±10 %	V/ Hz	230/60	230/50
Current input (at 230 V)	Α	4	4
Total power consumption	W	920	920
Weight	kg	15.5	15.5

## → Cooling capacity using a HSP booster pump:

If a HSP booster pump is used, the ultra-low refrigerated circulators do not reach the specified lowest bath temperature because the pump dissipates frictional heat of approx. 400 W to the bath fluid.