

# Case Study

## JULABO CF30, CF41, FL300

### Performance evaluation of JULABO models CF30, CF41, and FL300 with DrySyn COOL reaction systems

Case Study  
03-2008  
Page 1 from 7



#### Objective

The objective of each test was to evaluate the performance of various JULABO units with ASYNT's DrySyn COOL reaction system.

#### Test description

The experiment was conducted once each with an insulated and uninsulated DrySyn COOL reaction vessel. The DrySynCOOL module was equipped with a 1000 ml round-bottom flask filled one-third (330 ml) with ethanol. The times required to achieve the temperature control unit's lowest rated temperature inside the unit itself and inside the round-bottom flask were recorded.

#### Test conditions

Units	CF31 Cryo Compact Circulator CF41 Cryo Compact Circulator FL300 Recirculating Cooler
Mains voltage	230 Volt / 50 Hz
Ambient temperature	20 °C (room temperature)
Medium	Ethanol
Special unit configuration	Temperature control via JULABO EasyTemp



#### Results

The results show that the tested JULABO models are capable of achieving temperatures below -33 °C inside the 1000 ml flask in DrySyn COOL.

#### Compilation of actual measurement results

MODEL	INSULATION	LOWEST TEMP: IN MODEL / TIME	LOWEST TEMP: IN ROUND-BOTTOM / TIME
CF31 with DrySyn COOL	insulated	-30 °C / after 2h 20min	-25.2 °C / after 2h 45min
CF31 with DrySyn COOL	not insulated	-21 °C / after 2h 30min	-15.6 °C / after 3h 45min
CF41 with DrySyn COOL	insulated	-40 °C / after 2h 20min	-33.4 °C / after 3h 30min
CF41 with DrySyn COOL	not insulated	-40 °C / after 2h 40min	-27.9 °C / after 4h 00min
FL300 with DrySyn COOL	insulated	-20 °C / after 50min	-16.8 °C / after 1h 45min
FL300 with DrySyn COOL	not insulated	-20 °C / after 1h 10min	-11.3 °C / after 2h 10min

# Case Study

## JULABO CF30, CF41, FL300



Case Study  
03-2008  
Page 2 from 7



### Test 1

#### CF31 Asynth DrySyn COOL, insulated

Lowest temp. internally: achieved after::	-30 °C 2h 20min
Lowest temp. externally: achieved after:	-25.2 °C 2h 45min

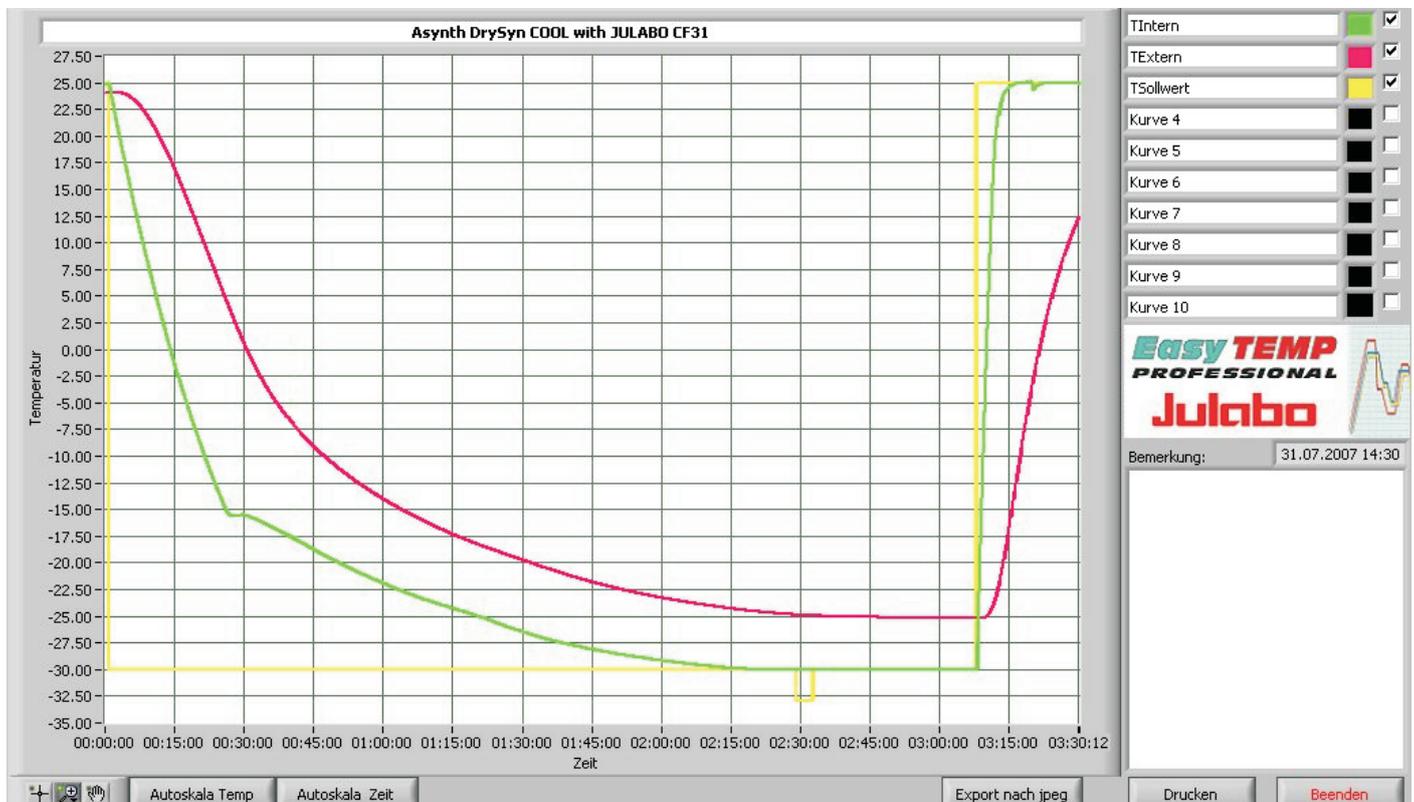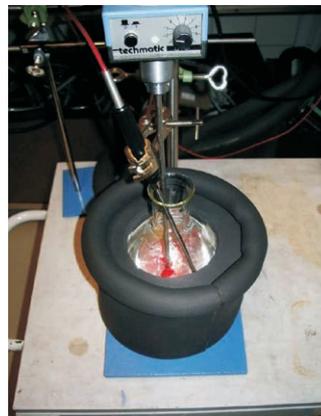

# Case Study

## JULABO CF30, CF41, FL300


Case Study  
03-2008  
Page 3 from 7



### Test 2

#### CF31 Asynth DrySyn COOL, not insulated

Lowest temp. internally: achieved after:	-30 °C Temperature not achieved, Lowest achieved temperature -21.0 °C after 2h 30min
Lowest temp. externally: achieved after:	-15.6 °C 3h 45min

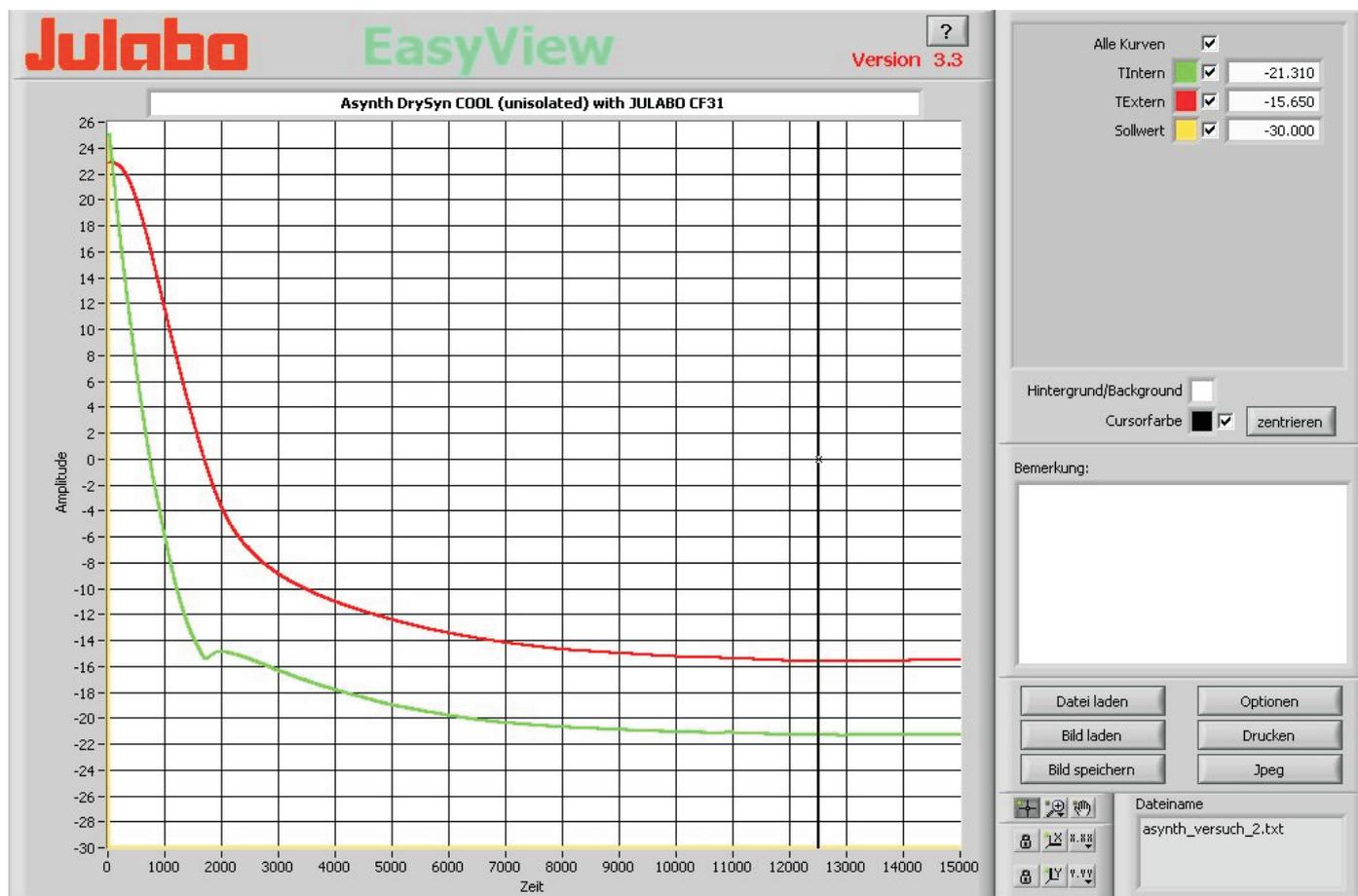

# Case Study

## JULABO CF30, CF41, FL300


Case Study  
03-2008  
Page 4 from 7



### Test 3

#### CF41 Asynt DrySyn COOL, insulated

Lowest temp. internally: -40 °C  
achieved after: 2h 20min  
Lowest temp. externally: -33.4 °C  
achieved after: 3h 30min

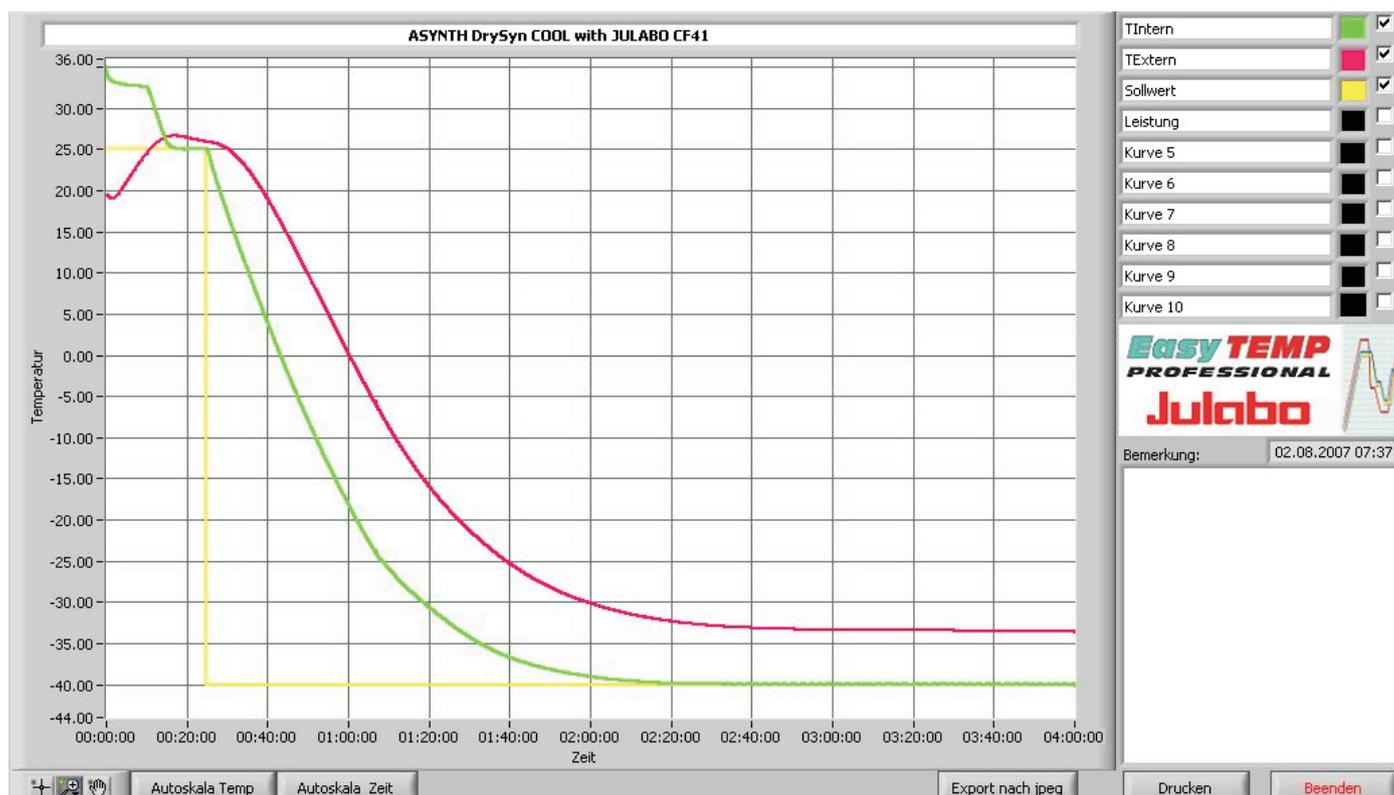

# Case Study

## JULABO CF30, CF41, FL300



Case Study  
03-2008  
Page 5 from 7



### Test 4

#### CF41 Asynt DrySyn COOL, not insulated

Lowest temp. internally:	-40 °C
achieved after:	2h 40min
Lowest temp. externally:	-27.9 °C
achieved after:	4h 00min

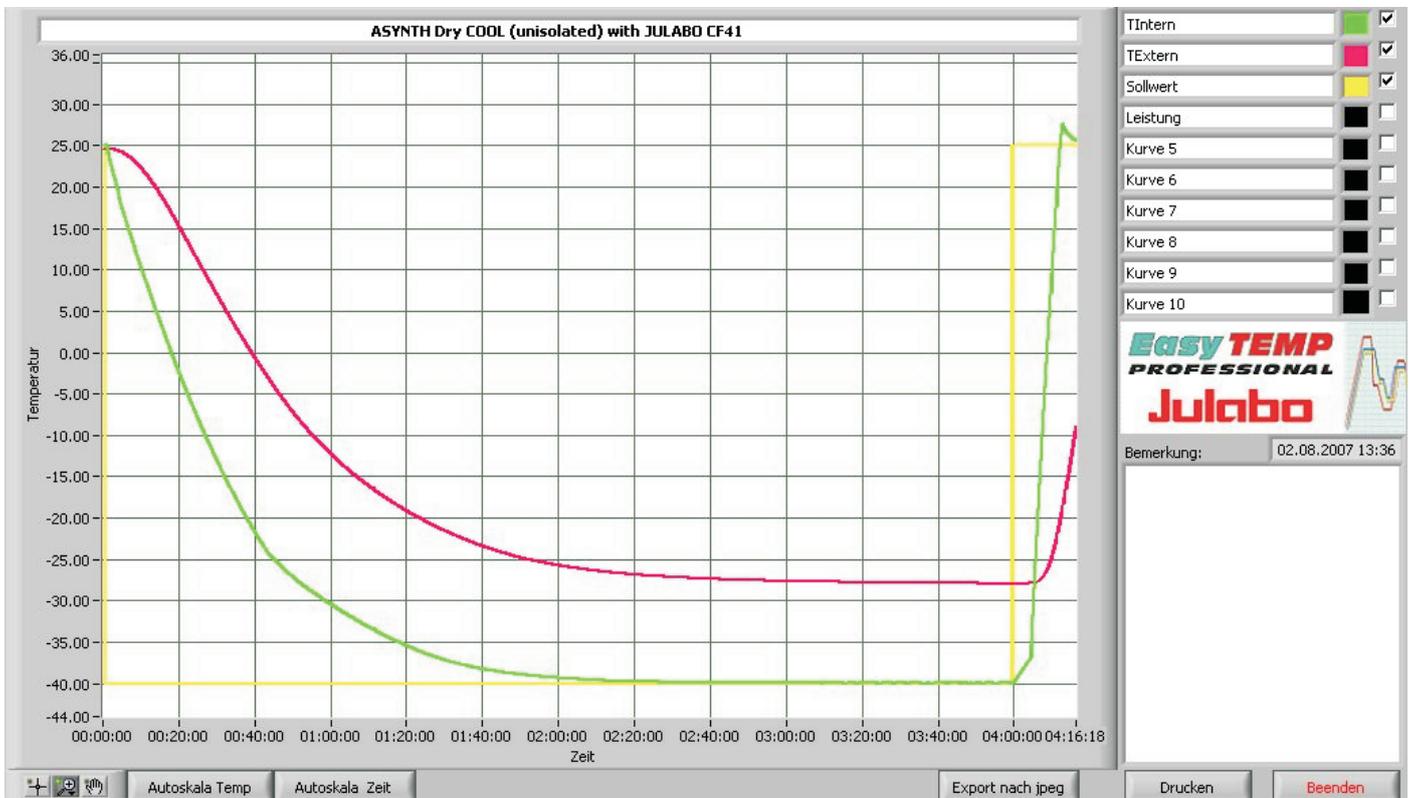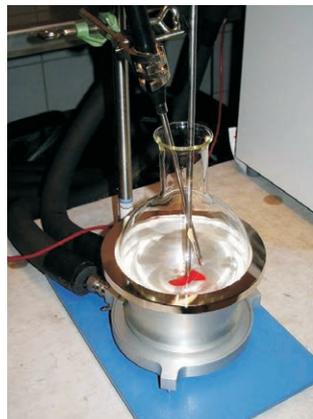

# Case Study

## JULABO CF30, CF41, FL300


Case Study  
03-2008  
Page 6 from 7



### Test 5

#### FL300 Asynth DrySyn COOL, insulated

Lowest temp. internally:	-20 °C
achieved after:	50min
Lowest temp. externally:	-16.8 °C
achieved after:	1h 45min

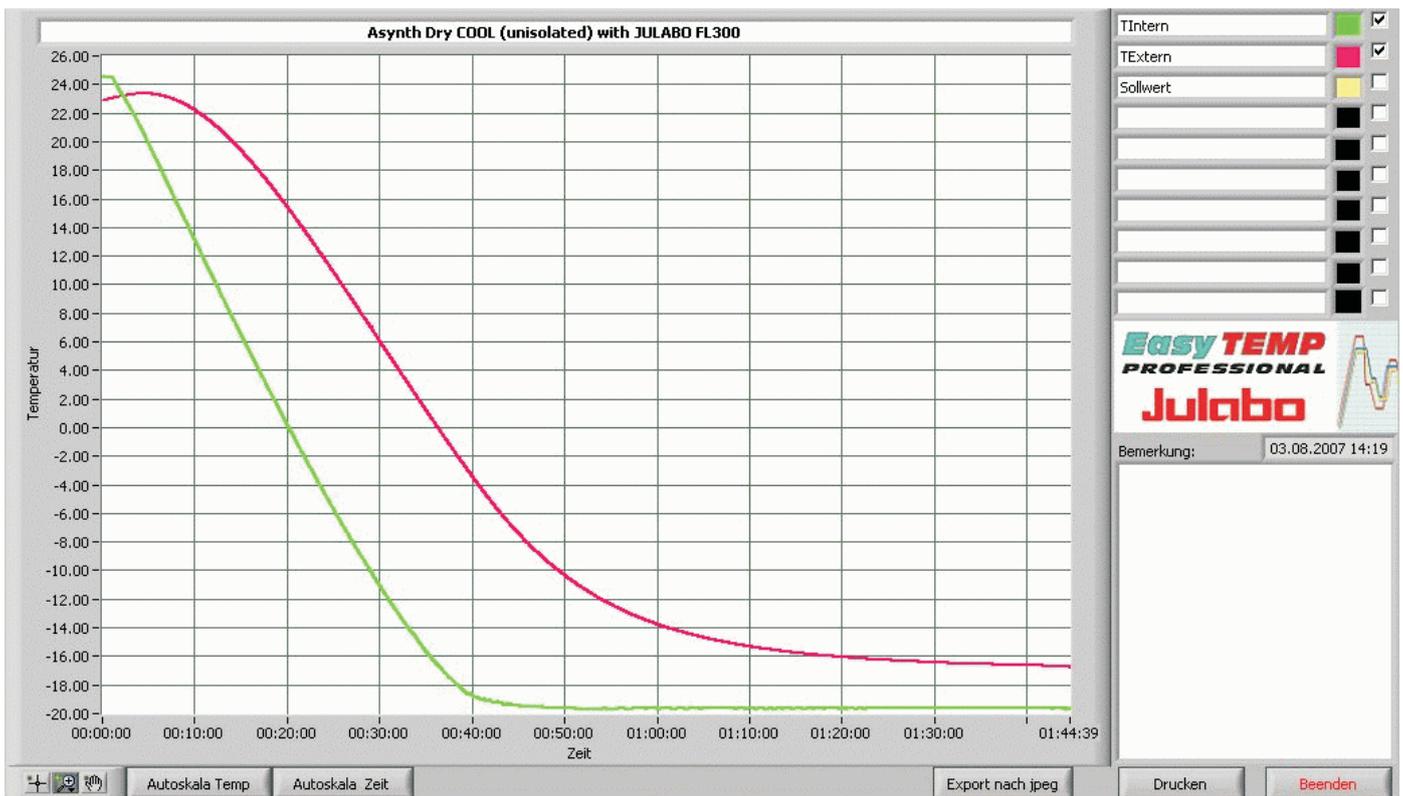

# Case Study

## JULABO CF30, CF41, FL300



Case Study  
03-2008  
Page 7 from 7



### Test 6

#### FL300 Asynth DrySyn COOL, not insulated

Lowest temp. internally:	-20 °C
achieved after:	1h 10min
Lowest temp. externally:	-11.3 °C
achieved after:	2h 10min

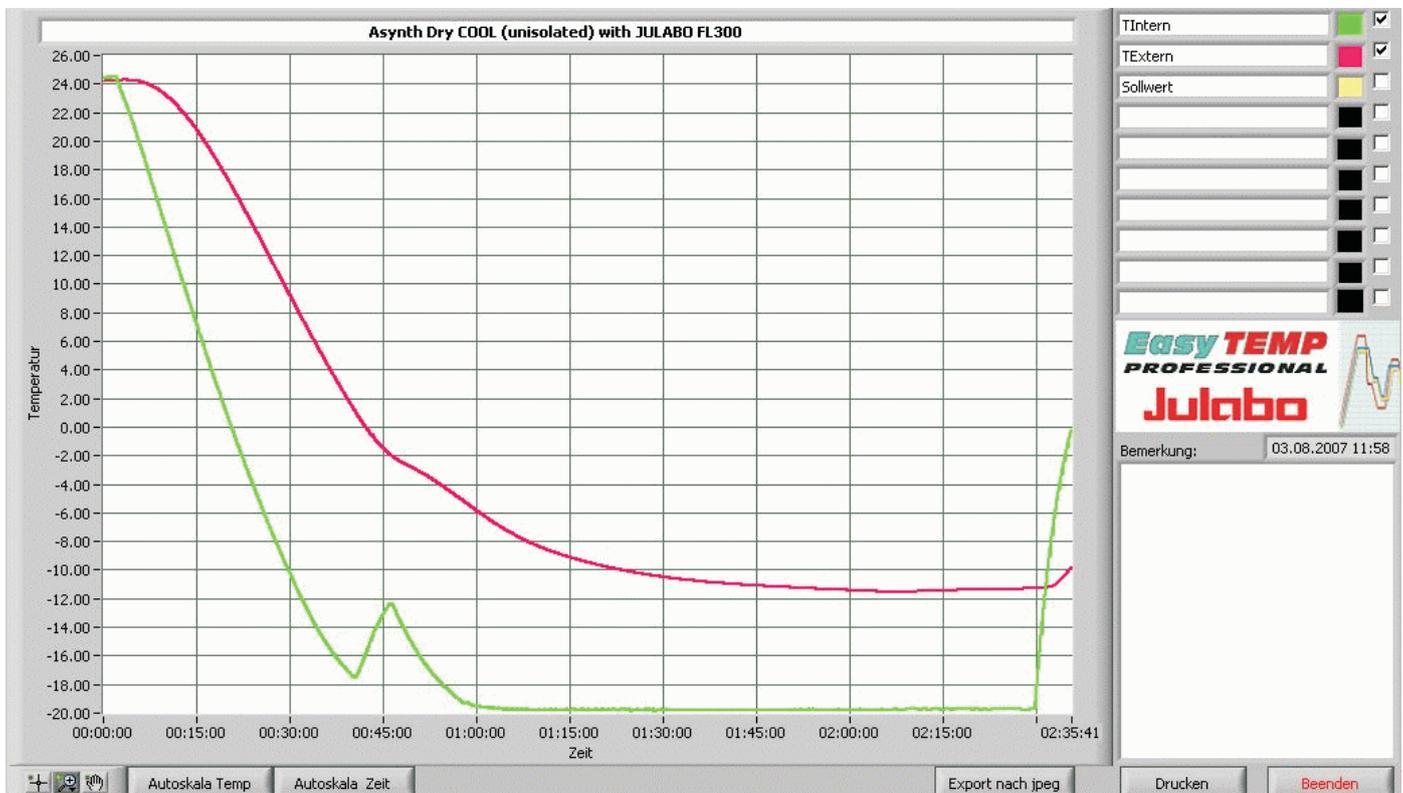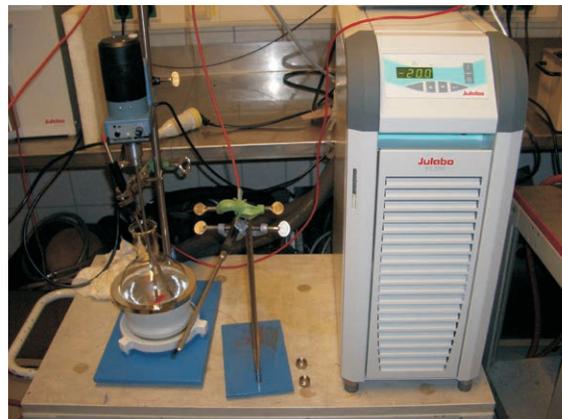

JULABO GmbH  
Eisenbahnstraße 45 • 77960 Seelbach / Germany  
Tel. +49 (0) 7823 51-0 • Fax +49 (0) 7823 24 91  
info@julabo.de • www.julabo.de

**Julabo**  
THE TEMPERATURE CONTROL COMPANY