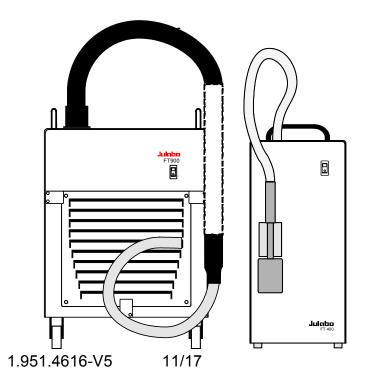
# **English**

# **OPERATING MANUAL**

Immersion Coolers FT200 FT400 FT900

Flow-Through Cooler FD200

JULABO GmbH 77960 Seelbach / Germany Tel. +49 (0) 7823 / 51-0 Fax +49 (0) 7823 / 24 91 info.de@julabo.com

www.julabo.com

#### Congratulations!

You have made an excellent choice.

JULABO thanks you for the trust you have placed in us.

This operating manual has been designed to help you gain an understanding of the operation and possible applications of our immersion coolers. For optimal utilization of all functions, we recommend that you thoroughly study this manual prior to beginning operation.

## The JULABO Quality Management System

Temperature control devices for research and industry are developed, produced, and distributed according to the requirements of ISO 9001 and ISO 14001. Certificate Registration No. 01 100044846

#### **Unpacking and inspecting**

Unpack the immersion cooler and accessories and inspect them for possible transport damage. Damage should be reported to the responsible carrier, railway, or postal authority, and a damage report should be requested. These instructions must be followed fully for us to guarantee our full support of your claim for protecting against loss from concealed damage. The form required for filing such a claim will be provided by the carrier.

Printed in Germany

Changes without prior notification reserved

Important: keep original operating manual for future use

## **TABLE OF CONTENTS**

OP	ERATING MANUAL	5
1.	INTENDED USE	5
	1.1. Description	5
2.	OPERATOR RESPONSIBILITY - SAFETY INSTRUCTIONS	5
	2.1. Disposal	7
	2.2. EC Conformity	8
	2.3. Warranty conditions	12
	2.4. Technical specifications	13
OP	ERATING INSTRUCTIONS	15
3.	OPERATING CONTROLS AND FUNCTIONAL ELEMENTS	15
4.	SAFETY NOTES FOR THE USER	17
	4.1. Explanation of safety notes	17
	4.2. Explanation of other notes	17
	4.3. Safety instructions	18
5.	PREPARATIONS	19
	5.1. Installation	19
	5.2. Immersion Probe	19
	5.3. Tube connection FD200	20
6.	OPERATING PROCEDURES	20
	6.1. Power connection	20
	6.2. Switching On	21
7.	TROUBLESHOOTING	22
8	CLEANING / REPAIRING THE UNIT	22

# **Operating manual**

#### 1. Intended use

JULABO immersion coolers have been designed for temperature application to specific fluids in a bath tank.

For example: Dewar vessels, beakers, or other containers in conjunction

with heating circulators for continuous countercooling

or for dry-ice substitution.

JULABO immersion coolers are not suitable for direct temperature control of foods, semi-luxury foods and tobacco, or pharmaceutical and medical products. Direct temperature control means unprotected contact of the object with the bath medium (bath fluid).

## 1.1. Description

The JULABO immersion coolers FT200, FT 400 and FT900 are employed to cool liquids for working temperatures ranging from +50 °C to -90 °C, such as in:

Dewar vessels, beakers, or other containers in conjunction with heating circulators for continuous countercooling or for dry-ice substitution.

The JULABO FD200 Flow-Through Cooler is employed to cool liquids in closed circuits. This unit is generally installed at the intake of a heating circulator to draw heat away from the circulating bath liquid.

Models FD200, FT200 and FT400 are provided with a handle for portable use.

Model FT900 is equipped with four castors. Two of the castors include locking levers that should be pressed down after setting up the unit to prevent it from moving.

The immersion probe is connected to the instrument with a flexible, specially insulated tube. On model FT900 the immersion probe is also flexible and may be adjusted precisely to different positions within the vessel.

# 2. Operator responsibility – Safety instructions

The products of JULABO ensure safe operation when installed, operated, and maintained according to common safety regulations. This section explains the potential dangers that may arise when operating the circulator and also specifies the most important safety precautions to preclude these dangers as far as possible.

The operator is responsible for the qualification of the personnel operating the units.

- The personnel operating the units should be regularly instructed about the dangers involved with their job activities as well as measures to avert these dangers.
- Make sure all persons tasked with operating, installing, and maintaining the unit have read and understand the safety information and operating instructions.
- When using hazardous materials or materials that could become hazardous, the circulator may be operated only by persons who are absolutely familiar with these materials and the circulator. These persons must be fully aware of possible risks.

If you have any questions concerning the operation of your unit or the information in this manual, please contact us!

**Contact** JULABO GmbH Tel. +49 (0) 7823 / 51-0

Fax +49 (0) 7823 / 24 91

Gerhard-Juchheim-Strasse 1 info.de@julabo.com 77960 Seelbach / Germany www.julabo.com

## Safety instructions for the operator:

- Avoid strikes to the housing, vibrations, damage to the operating-element panel (keypad, display), and contamination.
- ➤ Make sure the product is checked for proper condition regularly (depending on the conditions of use). Regularly check (at least every 2 years) the proper condition of the mandatory, warning, prohibition and safety labels.
- Make sure that the mains power supply has low impedance to avoid any negative effects on the instruments being operated on the same mains.
- ➤ This unit is designed for operation in a controlled electromagnetic environment. This means that transmitting devices (e.g., cellular phones) should not be used in the immediate vicinity.
- Magnetic radiation may affect other devices with components sensitive to magnetic fields (e.g., monitors). We recommend maintaining a minimum distance of 1 m.
- ➤ Permissible ambient temperature: max. 40 °C, min. 5 °C.
- Permissible relative humidity: 50% (40 °C).
- Do not store the unit in an aggressive atmosphere. Protect the unit from contamination.
- > Do not expose the unit to sunlight.

#### Appropriate operation

Only qualified personnel is authorized to configure, install, maintain, or repair the circulator. Persons who operate the circulator must be trained in the particular tasks by qualified personnel. The summarized user guidance (short manual) and the specification table with information on individual parameters are sufficient for this.

#### Use

The bath can be filled with flammable materials. Fire hazard!

There might be chemical dangers depending on the bath medium used.

Observe all warnings for the used materials (bath fluids) and the respective instructions (safety data sheets).

Insufficient ventilation may result in the formation of explosive mixtures. Only use the unit in well ventilated areas.

Only use recommended materials (bath fluids). Only use non-acid and non corroding materials.

When using hazardous materials or materials that could become hazardous, **the operator must** affix the enclosed safety labels **(1 + 2)** to the front of the unit so they are highly visible:

1

Warning label W00: Colors: yellow, black Danger area. Attention! Observe instructions. (operating manual, safety data sheet)

2

Mandatory label M018: Colors: blue, white

Carefully read the user information prior to beginning operation.

Scope: EU

or

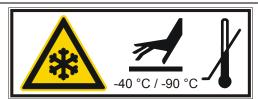
Semi S1-0701 Table A1-2 #9

Carefully read the user information prior to beginning operation.

Scope: USA, NAFTA

Particular care and attention is necessary because of the wide operating range. There are thermal dangers: Touchable parts of the probe can be very cold.

The user must attach the enclosed safety labels to the unit so they are well visible.

Safety label including warning label W017: Colors: yellow, black

Attention: Do not touch cold probe.

#### 2.1. Disposal

The product may be used with oil as bath fluid. These oils fully or partially consist of mineral oil or synthetic oil. For disposal, follow the instructions in the material safety data sheets.

This unit contains the refrigerants R134a R404A, and R-23, which at this time are not considered harmful to the ozone layer. However, over the long operating period of the unit, disposal rules may change. Therefore, only qualified personnel should handle the disposal.

Valid in EU countries

See the current official journal of the European Union – WEEE directive. Directive of the European Parliament and of the Council on waste electrical and electronic equipment (WEEE).

This directive requires electrical and electronic equipment marked with a crossed-out trash can to be disposed of separately in an environmentally friendly manner.

Contact an authorized waste management company in your country. Disposal with household waste (unsorted waste) or similar collections of municipal waste is not permitted!

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The above-named product is in compliance with the following harmonized standards and technical specifications:

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Refrigerating systems and heat pumps - Safety and environmental requirements - Part 1: Basics requirements, definitions, classification and selection criteria

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Refrigerating systems and heat pumps - Safety and environmental requirements - Part 2: Design, construction, testing, marking and documentation

EN 378-3: 2016

Refrigerating systems and heat pumps - Safety and environmental requirements - Part 3: Installation site and personal protection

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2017\_052\_FD200-Durchlaufkühler\_d\_e.docx

## 2.3. Warranty conditions

JULABO GmbH warrants its products against defects in material or in workmanship, when used under appropriate conditions and in accordance with appropriate operating instructions for a period of ONE YEAR.

Extension of the warranty period – free of charge

With the '1PLUS warranty' the user receives a free of charge extension to the warranty of up to 24 months, limited to a maximum of 10 000 working hours.

To apply for this extended warranty the user must register the unit on the JULABO web site <a href="https://www.julabo.com">www.julabo.com</a>, indicating the serial no. The extended warranty will apply from the date of JULABO GmbH's original invoice.

JULABO GmbH reserves the right to decide the validity of any warranty claim. In case of faults arising either due to faulty materials or workmanship, parts will be repaired or replaced free of charge, or a new replacement unit will be supplied.

Any other compensation claims are excluded from this guarantee.

# 2.4. Technical specifications

		ETOOO	ED000
		FT200	FD200
Temperature range	°C	-20 +30	10 +30
Cooling capacity	<u>°С</u>	+20 0 -30	+ <u>20 +10</u>
(medium ethanol)	kW	0.25 0.15 0.04	0.22 0.18
Refrigerant		R134a	R134a
Recommanded flow rate	l/min		2 - 3
Freezing protection	°C		10
Immerson probe (Lxdia.)	cm	9x4	
Connection tubing (L)	cm	120	
Dimensions (WxLxH)	cm	18x27x39	18x27x39
Weight	kg	18	16
Ambient temperature	°C	5 35	5 35
Mains power connection	V/Hz	190-253 /50	230/50
Current draw (at 230 V)	Α	2,0	2,0
Mains power connection	V/Hz	103-127 / 60	115 / 60
Current draw (at 115 V)	А	3,0	3,0

		FT400	FT900	
Temperature range	°C	-40 +30	-90 +30	
Cooling capacity (medium ethanol)	<u>°C</u> kW	+20 +10 -20 -40 0.45 0.36 0.14 0.03	+20 +10 -40 -80 0.3 0.27 0.2 0.07	
Cooling compressor		1-stage	2- stage	
Refrigerant		R404A	R404A/R23	
Immerson probe (Lxdia.)	cm	12x5	65x1.5 (flexible)	
Connection tubing (L)	cm	120	160	
Dimensions (WxLxH)	cm	20x30x43	38x55x60	
Weight	kg	24	50	
Ambient temperature	°C	5 35	5 35	
Mains power connection	V/Hz	230 / 50	230 / 50/60	
Current draw (at 230 V)	Α	3,0	6,0	
Mains power connection	V/Hz	115 / 60	115 / 60	
Current draw (at 115 V)	Α	4,0	7,0	

## Note:

All measurements have been carried out at: rated voltage and frequency; ambient temperature 20 °C;

#### **Environmental conditions according to IEC 61 010-1:**

Use indoors only.

Altitude up to 2000 m - normal zero.

Ambient temperature: see Technical specifications

Humidity:

Max. relative humidity 80% for temperatures up to +31 °C,

linear decrease down to 50% relative humidity at a temperature of +40 °C

Max. mains voltage fluctuations of ±10% are permissible.

Protection class according to IEC 60 529 IP2

The unit corresponds to Class I

Overvoltage category II Pollution degree 2

#### Caution:

The unit is not for use in explosive environment.

#### **EMC** requirements

The device is an ISM device of group 1 per CISPR 11 (uses HF for internal purposes) and is classified in class A (industrial and commercial sector).

#### NOTICE:

- Devices of class A are intended for the use in an industrial electromagnetic environment.
- When operating in other electromagnetic environments, their electromagnetic compatibility may be impacted.

#### Information about the used refrigerants

The Regulation (EU) No. 517/2014 on fluorinated greenhouse gases applies to all systems which contain fluorinated refrigerants and replaces (EC) 842/2006.

The aim of the Regulation is to protect the environment by reducing emissions of fluorinated greenhouse gases.

Among other things it regulates the emission limits, use and recovery of these substances. It also contains requirements for operators of systems which require / contain these substances to function.

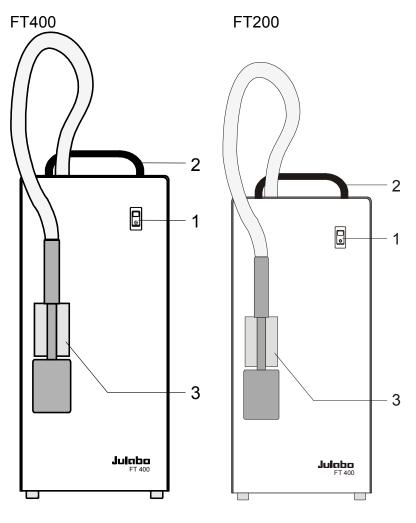
Under Regulation 517/2014, the operator of a system of this nature has the following duties:

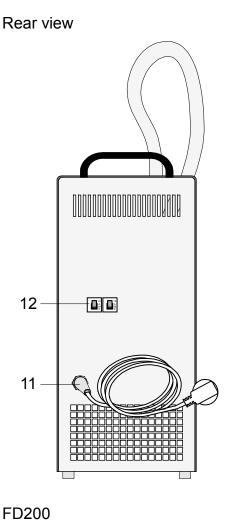
- The operator must ensure that the equipment is checked at regular intervals for leaks.
- These intervals depend on the CO<sub>2</sub> equivalent of the system. This is calculated from the refrigerant fill volume and type of refrigerant. The CO<sub>2</sub> equivalent of your system is shown on the model plate.
- The operator undertakes to have maintenance, repair, service, recovery and recycling work carried out by certified personnel who have been authorized by JULABO.
- All such work must be documented. The operator must keep records and archive them for at least five years. The records must be submitted to the relevant authority on request.

Refer to the text of the Regulation for further information.

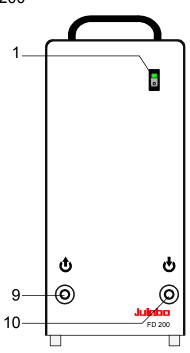
# **Operating instructions**

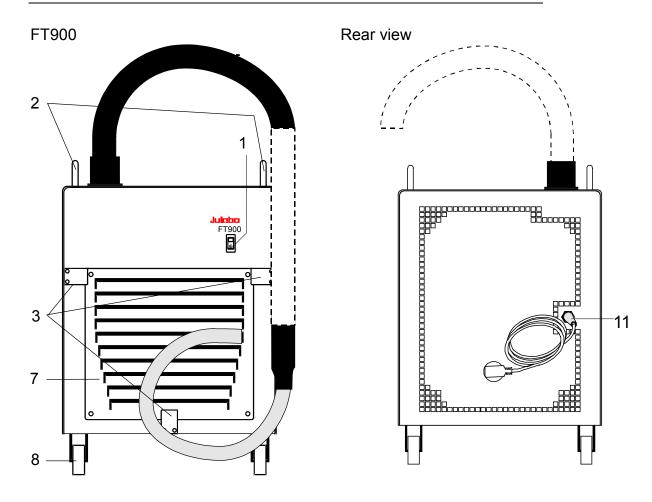
# 3. OPERATING CONTROLS AND FUNCTIONAL ELEMENTS

- 1 Mains switch, illuminated
- 2 Cooling control light
- 3 Clamp for immersion probe
- 7 Removable ventilation grid
- 8 Castor with locking lever
- 9 Tube connection discharge
- 10 Tube connection intake
- 11 Mains power cable with plug
- 12 Safety cutouts: Mains fuses 10 A

# 4. Safety notes for the user

## 4.1. Explanation of safety notes

In addition to the safety warnings listed above, warnings are posted throughout the manual. These warnings are designated by an exclamation mark inside an equilateral triangle. "Warning of a dangerous situation (Attention! Please follow the documentation)."

The danger is classified using a signal word.

Read and follow these important instructions.

## Warning:

Describes a possibly highly dangerous situation. If these instructions are not followed, serious injury and danger to life could result.

#### Caution:

Describes a possibly dangerous situation. If this is not avoided, slight or minor injuries could result. A warning of possible property damage may also be contained in the text.

#### Notice:

Describes a possibly harmful situation. If this is not avoided, the product or anything in its surroundings can be damaged.

### 4.2. Explanation of other notes

## Note!

Draws attention to something special.

#### Important!

Indicates usage tips and other useful information.

#### 4.3. Safety instructions

Follow the safety instructions to avoid personal injury and property damage. Also, the valid safety instructions for workplaces must be followed.

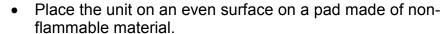
- Only connect the unit to a power socket with an earthing contact (PE protective earth)!
- The power supply plug serves as a safe disconnecting device from the line and must always be easily accessible.
- Place the unit on an even surface on a base made of nonflammable material.
- Do not stay in the area below the unit.
- Make sure you read and understand all instructions and safety precautions listed in this manual before installing or operating your unit.
- Do not touch the immersion probe if it is frosted.
- Do not bend the tube connection of the immersion probe
- Keep the air intake and exhaust grids free of obstructions. (Maintain a sufficient distance from all surrounding surfaces!)
- Do not move the unit from the position where it was set up during operation.
- Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures, or before moving the unit.
- Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit.
- Transport the unit with care.
- Sudden jolts or drops may cause damage in the interior of the unit.
- Observe all warning labels.
- Never remove warning labels.
- Never operate units with damaged mains power cables.
- Repairs are to be carried out only by qualified service personnel.

 There are thermal dangers: Touchable parts of the probe can be very cold. Therefore, exercise particular caution when touching these parts. Use gloves.

# 5. Preparations

#### 5.1. Installation

- Press down the castor levers on model FT900.
- The instrument should be set up at a frost-proof and dry location.
- The place of installation should be large enough and provide sufficient air ventilation to ensure the room does not warm up excessively because of the heat the instrument rejects to the environment. (Max. permissible ambient temperature: 35 °C). For a fault (leakage) in the refrigeration system, the standard EN 378 prescribes a certain room space to be available for each kg of refrigerant.

- > For 0.52 kg of refrigerant R404A, 1 m<sup>3</sup> of space is required.
- > For 0.68 kg of refrigerant R23, 1 m<sup>3</sup> of space is required.
- The ambient temperature must not exceed 35 °C.
- Keep at least 20 cm of open space on the front and rear venting grids.
- Do not set up the unit in the immediate vicinity of heat sources and do not expose to sun light.
- Before operating the unit after transport, <u>wait about one hour after</u>
   <u>setting it up.</u> This will allow any oil that has accumulated laterally
   during transport to flow back down thus ensuring maximum cooling
   performance of the compressor.

#### 5.2. Immersion Probe

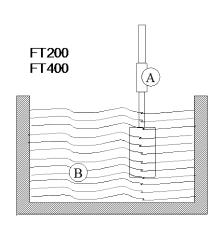
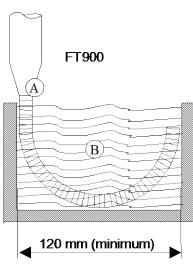
Avoid touching the immersion probe if it is frosted.

DANGER OF INJURY. Use gloves.

Switch the unit on only if the probe is immersed into the bath fluid.

To prevent the immersion probe (A) from icing, it should be completely immersed into the bath liquid (B).

Accessory: Clamp for cooler probe FT200/400 - order no. 8 970 400

#### 5.3. Tube connection FD200

• Connect the tubes and secure with tube clamps. discharge (9)

intake (10)

Recommended flow rate: 2 to 3 l/min

## 6. OPERATING PROCEDURES

#### 6.1. Power connection

#### Caution:

- Only connect the unit to a power socket with earthing contact (PE protective earth)!
- The power supply plug serves as safe disconnecting device from the line and must be always easily accessible.
- Never operate equipment with damaged mains power cables.
- Regularly check the mains power cables for material defects (e.g. for cracks).
- We disclaim all liability for damage caused by incorrect line voltages!

Make sure that the line voltage and frequency match the supply voltage specified on the type plate.

Deviations of ±10 % are permissible.

## 6.2. Switching On

The immersion cooler is turned on and off with the mains switch.
 (1).

The control light in the switch will illuminate.

The immersion probe – as part of the cooling circuit – should not be exposed to bath temperatures above the working temperature of the immersion cooler. This would cause damage to the compressor.

Do not immerse a frosted immersion probe into hot bath oil.

DANGER OF INJURY!

## 7. TROUBLESHOOTING

Malfunction of compressor:

The cooling compressor is equipped with an overload protection device that will be triggered by overheating or excessive current consumption. Possible causes include insufficient ventilation or contamination of the condenser. After a cool-down phase, the motor is automatically switched on again.

• Interruption of the cooling loop (FD200) by a bended tube.

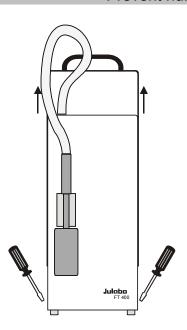
# 8. Cleaning / repairing the unit

#### Caution:

Before cleaning the unit, disconnect the power plug from the mains socket! Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures.

Service and repair work may be performed only by authorized electricians. Prevent humidity from entering into the immersion cooler.

JULABO coolers are designed for continuous operation under normal conditions.

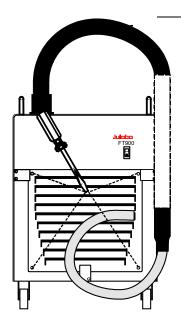
Periodic maintenance is not required.

Regularly check the condensor for dirt contamination. Clean the ribbed condensor, because dust and dirt will reduce cooling performance of the unit.

Cleaning the Cooling Compressor:

Switch off the unit, disconnect mains power cable.

Remove the hood (FD200, FT200, FT400).

- The ventilation grid (7) is detached by unscrewing the four mouting screws (FT900).
- Clean the ribbed condensor with a vacuum cleaner.
- Replace the hood or the ventilation grid.
- Switch on the unit.

Clean the outside of the unit using a wet cloth and low surface tension water.

## Repairs

Before asking for a service technician or returning a JULABO instrument for repair, please contact an authorized JULABO service station.

When returning the unit:

- Clean the unit in order to avoid any harm to the service personnel.
- Attach a short fault description.
- During transport the unit has to stand upright. Mark the packing correspondingly.
- When returning a unit, take care of careful and adequate packing.
- JULABO is not responsible for damages that might occur from insufficient packing.