
# **OPERATING MANUAL**

# ME

Beer Forcing Test Refrigerated/Heating Circulating Bath F38-ME

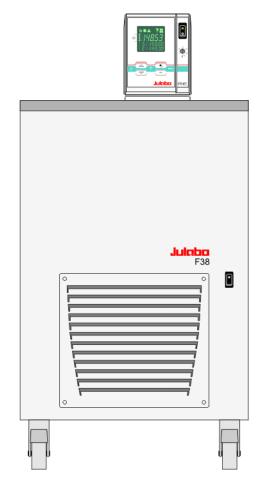



JULABO GmbH 77960 Seelbach / Germany Tel. +49 (0) 7823 / 51-0 Fax +49 (0) 7823 / 24 91 info.de@julabo.com www.julabo.com

1.951.0541-V4

11/17

#### **Congratulations!**

You have made an excellent choice.

JULABO thanks you for the trust you have placed in us.

This operating manual has been designed to help you gain an understanding of the operation and possible applications of our circulators. For optimal utilization of all functions, we recommend that you thoroughly study this manual prior to beginning operation.

### The easy way to the Beer Forcing Test

Notice:

This abbreviated version does not release the user from the duty to instruct the operating personnel on the work with this unit and particularly to point out to possible dangers.

For the general use in the laboratory the refrigerated and heating circulator F38-ME offers a lot of setting possibilities, which are not necessary for doing a "Forcing Test". In the chapter "The easy way to the Forcing Test" the advices for the necessary settings are summarised.

Advices which are particularly important for doing a forcing test, are marked with the symbol of a bottle (see left side).

### The JULABO Quality Management System



Temperature control devices for research and industry are developed, produced, and distributed according to the requirements of ISO 9001 and ISO 14001. Certificate Registration No. 01 100044846

### Unpacking and inspecting

Unpack the circulator and accessories and inspect them for possible transport damage. Damage should be reported to the responsible carrier, railway, or postal authority, and a damage report should be requested. These instructions must be followed to the full extent for us to guarantee our full support of your claim for protection against loss from concealed damage. The form required for filing such a claim will be provided by the carrier

Printed in Germany

Changes without prior notification reserved

page 39

Important: keep operating manual for future use

### TABLE OF CONTENTS


Opera	ting manual	. 5
1. In	tended use	. 5
1.1.	Description	. 5
2. O	perator responsibility – Safety recommendations	6
2.1.	Disposal	. 7
2.2.	EC Conformity	. 8
2.3.	Warranty conditions	10
2.4.	Technical specifications	11
Opera	ting instructions	14
3. Sa	afety notes for the user	14
3.1.	Explanation of safety notes	14
3.2.	Explanation of other notes	14
3.3.	Safety recommendations	15
4. O	perating controls and functional elements	17
4.1.	Circulator	17
4.2.	Cooling Machine	19
5. Pr	reparations	20
5.1.	Installation	20
5.2.	Bath fluids	21
5.3.	Filling / draining	22
6. O	perating procedures	23
6.1.	Power connection	23
6.2.	Switching on / Start - Stop	23
	2.1.Switching on the circulator	
6.	2.2.Switching on the Cooling Machine	
7.	Setting of temperatures	25
7.1.	Using the pre-settings in the <b>T</b> menu	25
7.2.	Direct setting of temperatures	26
8.	Safety installations, warning functions	27
8.1.	Excess temperature protection	27
8.	1.1.Early warning system, low level protection	28
8.2.	Switch-over from warning to shutdown function	
8.3.	Over and Sub temperature warning function	30

9. Menu functions	
9.1. MENU PROGRAM – START	
-	
9.1.2.START Mode Standard	
	ration35
-	
-	
	ıre43
	stics and parameters47
	XTERNAL
9.5.2.Dynamic internal	
-	I-, TV- EXTERNAL49
	NTERNAL
9.6. MENU SERIAL - BAUDRATE, HANDS	HAKE, PARITY51
-	ibration52
	RNAL54
	INT55 rnal control56
	61
	64
12 Remote control	
	dinated data system66
· · · · · ·	•
12.5. Error messages	
13. JULABO Service – Online remote diagnos	iis71
14. Cleaning / repairing the unit	

## **Operating manual**

## 1. Intended use

JULABO circulators have been designed to control the temperature of specific fluids in a bath tank. The JULABO forcing test refrigerated/heating circulating bath in connection with a photometer determines the product life of beer without clouding.

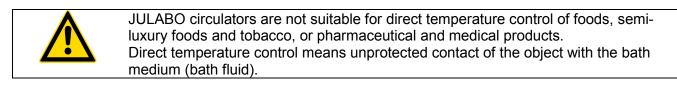

### 1.1. Description

5 15000 E 14530	V	The circulators are operated via the splash-proof keypad. The microprocessor technology allows different values to be set, stored, and displayed on the VFD COMFORT-DISPLAY. Three menu keys facilitate the adjustment of setpoints, warning and safety functions, and menu functions.
	V	Temperature- and time-dependent processes can be stored and executed using the integrated programmer.
	V	The adjustable PID cascade temperature control automatically adjusts the heat supply to the thermal requirements of the bath.
PID3	V	Absolute Temperature Calibration (ATC3) provides high temperature stability at all points in the bath. With the 3-point calibration, an offset is adjusted at three
ATC <sup>3</sup>	<u>ک</u> ا	temperatures to ensure an accurate temperature pattern at the selected spot in the bath over the entire temperature range. Electrical connections:
RS232	V	RS232 interface for modern process technology without an additional interface. Connection for external Pt100 sensor for external temperature measurement and control.
Pt100		Alarm output for external alarm message or control of JULABO refrigerating baths or solenoid valve (cooling water).
40, 100 40, 100 230 c c	V	The excess temperature protection according to IEC 61010-2-010 is a safety device independent from the control circuit. Its safety value can be displayed and adjusted on the VFD COMFORT-DISPLAY.
	V	The early warning system for low level signals that bath fluid needs to be refilled before the low level protection according to IEC 61010-2-010 triggers a safety shutdown of the main functional elements.
smart <b>PUMP</b>	V	Intelligent pump system: the pump capacity (electronically adjustable via the motor speed) can be adapted to different conditions for internal and external temperature-control applications.

## 2. Operator responsibility – Safety recommendations

The products of JULABO ensure safe operation when installed, operated, and maintained according to common safety regulations. This section explains the potential dangers that may arise when operating the circulator and also specifies the most important safety precautions to preclude these dangers as far as possible.

- > The operator is responsible for the qualification of the personnel operating the units.
- The personnel operating the units should be regularly instructed about the dangers involved with their job activities as well as measures to avert these dangers.
- Make sure all persons tasked with operating, installing, and maintaining the unit have read and understand the safety information and operating instructions.
- When using hazardous materials or materials that could become hazardous, the circulator may be operated only by persons who are absolutely familiar with these materials and the circulator. These persons must be fully aware of possible risks.

If you have any questions concerning the operation of your unit or the information in this manual, please contact us!

Contact	JULABO GmbH	Tel. +49 (0) 7823 / 51-0
	Gerhard-Juchheim-Strasse 1	Fax +49 (0) 7823 / 24 91
	77960 Seelbach / Germany	info.de@julabo.com
	-	www.julabo.com

#### Safety instructions for the operator:

- You have received a product designed for industrial use. Nevertheless, avoid strikes to the housing, vibrations, damage to the operating-element panel (keypad, display), and contamination.
- Make sure the product is checked for proper condition regularly (depending on the conditions of use). Regularly check (at least every 2 years) the proper condition of the mandatory, warning, prohibition and safety labels.
- Make sure that the mains power supply has low impedance to avoid any negative effects on instruments being operated on the same mains.
- This unit is designed for operation in a controlled electromagnetic environment. This means that transmitting devices (e.g., cellular phones) should not be used in the immediate vicinity. Magnetic radiation may affect other devices with components sensitive to magnetic fields (e.g., monitors). We recommend maintaining a minimum distance of 1 m.
- Permissible ambient temperature: max. 40 °C, min. 5 °C.
- Permissible relative humidity: 50% (40 °C).
- > Do not store the unit in an aggressive atmosphere.
- > Protect the unit from contamination.
- > Do not expose the unit to sunlight.

#### Appropriate operation

Only qualified personnel is authorized to perform configuration, installation, maintenance and repairs of the circulator.

Routine operation can also be carried out by untrained personnel who should however be instructed by trained personnel.

#### Use:

The bath can be filled with flammable materials. Fire hazard!

There might be chemical dangers depending on the bath medium used.

Observe all warnings for the used materials (bath fluids) and the respective instructions (safety data sheets).

Insufficient ventilation may result in the formation of explosive mixtures. Only use the unit in well ventilated areas.

Only use recommended materials (bath fluids). Only use non-acid and non corroding materials.

When using hazardous materials or materials that could become hazardous, **the operator must** affix the enclosed safety labels (1 + 2) to the front of the unit so they are highly visible:

1		Warning label W00: Colors: yellow, black Danger area. Attention! Observe instructions. (operating manual, safety data sheet)
2 or	Ē	Mandatory label M018: Colors: blue, white Carefully read the user information prior to beginning operation. <b>Scope: EU</b>
2		Semi S1-0701 Table A1-2 #9 Carefully read the user information prior to beginning operation. <b>Scope: USA, NAFTA</b>

Particular care and attention is necessary because of the wide operating range. There are thermal dangers: Burn, scald, hot steam, hot parts and surfaces that can be touched.



Warning label W26: Colors: yellow, black Hot surface warning. (The label is put on by JULABO)

Observe the instructions in the manuals for instruments of a different make that you connect to the circulator, particularly the respective safety recommendations. Also observe the pin assignment of plugs and technical specifications of the products.


### 2.1. Disposal

The circulator contains a back-up battery that supplies voltage to memory chips when the unit is switched off. Do not dispose of the battery with household waste!

Depending on battery regulations in your country, you might be obliged to give back used or defect batteries to gathering places.

The product may be used with oil as bath fluid. These oils fully or partially consist of mineral oil or synthetic oil. For disposal, observe the instructions in the safety data sheets.

These units contains refrigerants– at this time considered not to have any negative effects on the ozone layer. However, during the long operating period of the unit, disposal prescriptions may change. So only qualified personnel should take care of disposal.

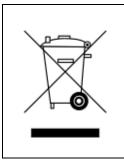

### Valid in EU countries

See the current official journal of the European Union – WEEE directive. Directive of the European Parliament and of the Council on waste electrical and electronic equipment (WEEE).

This directive requires electrical and electronic equipment marked with a crossedout trash can to be disposed of separately in an environmentally friendly manner. Contact an authorized waste management company in your country.

Disposal with household waste (unsorted waste) or similar collections of municipal waste is not permitted!

## 2.2. EC Conformity

#### EG-Konformitätserklärung nach EG Maschinenrichtlinie 2006/42/EG, Anhang II A EC-Declaration of Conformity to EC Machinery Directive 2006/42/EC, Annex II A

Hersteller / Manufa	acturer:	JULABO GmbH Gerhard-Juchheim-Straße 1 77960 Seelbach / Germany Tel: +49(0)7823 / 51 - 0	CE
Hiermit erklären wir We hereby declare, tha	, dass das nachfolgend bezeichne		
	Thermostat / Circulator		
Тур / Туре:	MA, MB, ME	Serien-Nr. / Serial-No.:	siehe Typenschild / see type label
Sicherheits- und Ges due to the design and d	sundheitsanforderungen den nach	ns in Verkehr gebrachten Ausführu nfolgend aufgeführten EG-Richtlinie eted by our Company – complies with f	en entspricht.
EMV-Richtlinie 20 EMV-Richtlinie 20	nie 2006/42/EG; Machinery Dire 004/108/EG; EMC-Directive 2004/ 014/30/EU; EMC-Directive 2014/3 2011/65/EU; RoHS-Directive 2011	/108/EC (bis zum / until 19. April 201 0/EU (vom / from 20. April 2016)	6)
Angewandte harn The above-named pro	nonisierte Normen und tech oduct is in compliance with the follo	n. Spezifikationen: owing harmonized standards and teo	chnical specifications:
EN 50581 : 2012 Technische Dokumentation Technical documentation for	n zur Beurteilung von Elektro- und Elektronikgerät or the assessment of electrical and electronic proc	en hinsichtlich der Beschränkung gefährlicher Stoffe lucts with respect to the restriction of hazardous sub	stances
EN ISO 12100 : 2 Sicherheit von Maschinen Safety of machinery - Gene	010 - Allgemeine Gestaltungsleitsätze - Risikobeurteilu eral principles for design - Risk assessment and ri	ung und Risikominderung (ISO 12100:2010) sk reduction (ISO 12100:2010)	
EN 61010-1 : 201 Sicherheitsbestimmungen Safety requirements for ele	0 für elektrische Mess-, Steuer-, Regel- und Laborg cctrical equiment for measurement, control, and fa	eräte, Teil 1: Allgemeine Anforderungen boratory use, Part 1: General requirements	
EN 61010-2-010 : Sicherheitsbestimungen Safety requirements for ele materials	für elektrische Mess-, Steuer-, Regel- und Laborge	eräte, Teil 2-010: Besondere Anforderungen an Lab boratory use, Part 2-010: Particular requirements foi	orgeräte für das Erhitzen von Stoffen r laboratory equipment for the heating of
EN 61326-1 : 201 Elektrische Mess-, Steuer- Electrical equipment for me	3 , Regel- und Laborgeräte- EMV-Anforderungen- T easurement, control, and laboratory use - EMC rec	eil 1: Allgemeine Anforderungen juirements - Part 1: General requirements	
Authorized represent	für die Zusammenstellung o ative in charge of administering teo e, im Hause / on the manufacturer's	chnical documentation:	
	erklärung wurde ausgestellt nformity was issued and valid of	g,	1.11

Seelbach, 22.02.2016

M 1. M. Juchheim, Geschäftsführer / Managing Director

2016\_007\_MA-MB-ME-Thermostat\_d\_e.docx

#### EG-Konformitätserklärung nach EG Maschinenrichtlinie 2006/42/EG, Anhang II A EC-Declaration of Conformity to EC Machinery Directive 2006/42/EC, Annex II A

Hersteller /	Manufacturer:
--------------	---------------

JULABO GmbH Gerhard-Juchheim-Straße 1 77960 Seelbach / Germany Tel: +49(0)7823 / 51 - 0 CE

Hiermit erklären wir , dass das nachfolgend bezeichnete Produkt We hereby declare, that the following product

Produkt / Product: Kältegerät / Refrigeration Unit

**Тур /** *туре*: F38

Serien-Nr. / Serial-No.: siehe Typenschild / see type label

aufgrund seiner Konzipierung und Bauart in der von uns in Verkehr gebrachten Ausführung den grundlegenden Sicherheits- und Gesundheitsanforderungen den nachfolgend aufgeführten EG-Richtlinien entspricht. due to the design and construction, as assembled and marketed by our Company – complies with fundamental safety and health requirements according to the following EC-Directives.

Maschinenrichtlinie 2006/42/EG; Machinery Directive 2006/42/EC EMV-Richtlinie 2014/30/EU; EMC-Directive 2014/30/EU RoHS-Richtlinie 2011/65/EU; RoHS-Directive 2011/65/EU

#### Angewandte harmonisierte Normen und techn. Spezifikationen: The above-named product is in compliance with the following harmonized standards and technical specifications:

#### EN 50581 : 2012

Technische Dokumentation zur Beurteilung von Elektro- und Elektronikgeräten hinsichtlich der Beschränkung gefährlicher Stoffe Technical documentation for the assessment of electrical and electronic products with respect to the restriction of hazardous substances

EN ISO 12100 : 2010

Sicherheit von Maschinen - Allgemeine Gestaltungsleitsätze - Risikobeurteilung und Risikominderung (ISO 12100:2010) Safety of machinery - General principles for design - Risk assessment and risk reduction (ISO 12100:2010)

EN 61010-1 : 2010

Sicherheitsbestimmungen für elektrische Mess-, Steuer-, Regel- und Laborgeräte, Teil 1: Allgemeine Anforderungen Safety requirements for electrical equiment for measurement, control, and laboratory use, Part 1: General requirements

#### EN 61010-2-010 : 2014

Scherheitsbestimmungen für elektrische Mess-, Steuer-, Regel- und Laborgeräte, Teil 2-010: Besondere Anforderungen an Laborgeräte für das Erhitzen von Stoffen Safety requirements for eletrical equipment for measurement, control, and laboratory use, Part 2-010: Particular requirements for laboratory equipment for the heating of materiale

#### EN 61326-1 : 2013

Elektrische Mess-, Steuer-, Regel- und Laborgeräte- EMV-Anforderungen- Teil 1: Allgemeine Anforderungen Electrical equipment for measurement, control, and laboratory use - EMC requirements - Part 1: General requirements

#### EN 378-1 : 2016

Kälteanlagen und Wärmepumpen – Sicherheitstechnische und umweltrelevante Anforderungen – Teil 1: Grundlegende Anforderungen, Begriffe, Klassifikationen und

Auswahlkriterien Refrigerating systems and heat pumps - Safety and environmental requirements - Part 1: Basics requirements, definitions, classification and selection criteria

#### EN 378-2 : 2016

Kälteanlagen und Wärmepumpen – Sicherheitstechnische und umweltrelevante Anforderungen – Teil 2: Konstruktion, Herstellung, Prüfung, Kennzeichnung und Dokumentation Refrigerating systems and heat pumps - Safety and environmental requirements - Part 2: Design, construction, testing, marking and documentation

#### EN 378-3 · 2016

Kälteanlagen und Wärmepumpen – Sicherheitstechnische und umweltrelevante Anforderungen – Teil 3: Aufstellungsort und Schutz von Personen Refrigerating systems and heat pumps - Safety and environmental requirements - Part 3: Installation site and personal protection

EN 378-4 : 2016

Kälteanlagen und Wärnepumpen – Sicherheitstechnische und umweltrelevante Anforderungen – Teil 4: Betrieb, Instandhaltung, Instandsetzung und Rückgewinnung Refrigerating systems and heat pumps - Safety and environmental requirements - Part 4: Operation, maintenance, repair and recovery

#### Bevollmächtigter für die Zusammenstellung der techn. Unterlagen: Authorized representative in charge of administering technical documentation:

Hr. Torsten Kauschke, im Hause / on the manufacturer's premises as defined above

#### Die Konformitätserklärung wurde ausgestellt The declaration of conformity was issued and valid of

Seelbach, 05.10.2017

M. Juchheim, Geschäftsführer / Managing Director

2017\_015\_F38-Kältegerät\_d\_e.docx

### 2.3. Warranty conditions

JULABO GmbH warrants its products against defects in material or in workmanship, when used under appropriate conditions and in accordance with appropriate operating instructions

#### for a period of ONE YEAR.

Extension of the warranty period - free of charge



With the '1PLUS warranty' the user receives a free of charge extension to the warranty of up to 24 months, limited to a maximum of 10 000 working hours.

To apply for this extended warranty the user must register the unit on the JULABO web site www.julabo.com, indicating the serial no. The extended warranty will apply from the date of JULABO GmbH's original invoice.

JULABO GmbH reserves the right to decide the validity of any warranty claim. In case of faults arising either due to faulty materials or workmanship, parts will be repaired or replaced free of charge, or a new replacement unit will be supplied.

Any other compensation claims are excluded from this guarantee.

		-	F38-ME
Working temperature range		°C	-38 80
Temperature stability		°C	±0,05
		•	
Temperature selection			digital
via keypad			VFD COMFORT-Display
remote control via personal comp	outer		indication on monitor
Temperature indication			VFD COMFORT-Display
Resolution		°C	0.01
Absolute Temperature Calibration	INT/EXT	°C	±3 / ±9
Temperature control			PID3 cascade temperature control
Heater wattage (at 230 V)		kW	2,0
Cooling capacity		°C	<u>+20 0 -20</u>
Medium ethanol		kW	0.92 0.66 0.32
Refrigerant			R404A
Electronically adj. pump capacity	stages		1 4
Flow rate	at 0 bar	l/min	11 16
Max. pressure	at 0 liters	bar	0.23 0.45
Electrical connections:			04.03/41.7
External alarm output			24-0 Vdc / max. 25 mA
Computer interface			RS232
External temperature sensor			Pt100
Overall dimensions	(WxDxH)	cm	44x64x82
Bath opening	(WxDxII) (WxL)	cm	35x41
Bath depth		cm	27
•	from to	liters	45
Weight		kg	80
Ambient temperature		°C	5 40
		0	5 10
Mains power connection		V/ Hz	230 / 50
Current draw (	(at 230 V)	А	14
Current draw GB, CH (	(at 230 V)	А	<9+5>
Mains power connection		V/ Hz	230 / 60
Current draw (	(at 230 V)	А	14

All measurements have been carried out at:

rated voltage and frequency ambient temperature: 20 °C Technical changes without prior notification reserved.

Safety installations according to IEC 61010-2-010: Excess temperature protection Low liquid level protection Classification according to DIN 12876-1	adjustable from 0 °C 230 °C float switch class III
Supplementary safety installations Early warning system for low level High temperature warning function Low temperature warning function Supervision of working sensor Reciprocal sensor monitoring between working and safety sensors Alarm message Warning message	float switch optical + audible (in intervals) optical + audible (in intervals) plausibility control difference >35 K optical + audible (permanent) optical + audible (in intervals)
Environmental conditions according to IEC 61 010-	1:

Use only indoor. Altitude up to 2000 m - normal zero. Ambient temperature: +5 ... +40 °C Air humidity: Max. rel. humidity 80 % for temperatures up to +31 °C, linear decrease down to 50 % relative humidity at a temperature of +40 °C Max. mains fluctuations of ±10 % are permissible.

Protection class according to IEC 60 529	IP21
The unit corresponds to Class I	
Overvoltage category	II
Pollution degree	2



#### Caution:

The unit is not for use in explosive environment

#### **EMC requirements**

The device is an ISM device of group 1 per CISPR 11 (uses HF for internal purposes) and is classified in class A (industrial and commercial sector).

#### NOTICE

- Devices of class A are intended for the use in an industrial electromagnetic environment.
- When operating in other electromagnetic environments, their electromagnetic compatibility may be impacted.

#### Information about the used refrigerants

The **Regulation (EU) No. 517/2014 on fluorinated greenhouse gases** applies to all systems which contain fluorinated refrigerants and replaces (EC) 842/2006.

The aim of the Regulation is to protect the environment by reducing emissions of fluorinated greenhouse gases.

Among other things it regulates the emission limits, use and recovery of these substances. It also contains requirements for operators of systems which require / contain these substances to function.

Under Regulation 517/2014, the operator of a system of this nature has the following duties:

- The operator must ensure that the equipment is checked at regular intervals for leaks.
- These intervals depend on the CO<sub>2</sub> equivalent of the system. This is calculated from the refrigerant fill volume and type of refrigerant. The CO<sub>2</sub> equivalent of your system is shown on the model plate.
- The operator undertakes to have maintenance, repair, service, recovery and recycling work carried out by certified personnel who have been authorized by JULABO.
- All such work must be documented. The operator must keep records and archive them for at least five years. The records must be submitted to the relevant authority on request.

Refer to the text of the Regulation for further information.

## **Operating instructions**

#### 3. Safety notes for the user

#### Explanation of safety notes 3.1.

In addition to the safety warnings listed, warnings are posted throughout the operating manual. These warnings are designated by an exclamation mark inside an equilateral triangle. "Warning of a dangerous situation (Attention! Please follow the documentation)." The danger is classified using a signal word. Read and follow these important instructions for averting dangers.
Warning: Describes a <b>possibly</b> highly dangerous situation. If these instructions are not followed, serious injury and danger to life could result.
<b>Caution:</b> Describes a <b>possibly</b> dangerous situation. If this is not avoided, slight or minor injuries could result. A warning of possible property damage may also be contained in the text.
Nation

Notice: Describes a **possibly** harmful situation. If this is not avoided, the product or anything in its surroundings can be damaged.

#### 3.2. Explanation of other notes

	<b>Note!</b> Draws attention to something special.
í	Important! Indicates usage tips and other useful information.
) <u> </u>	This icon is used in the operating instructions to indicate flashing values or parameters which have to be set or confirmed.

### 3.3. Safety recommendations

Follow the safety instructions to avoid personal injury and property damage. Also, the valid safety instructions for workplaces must be followed.



- Only connect the unit to a power socket with an earthing contact (PE protective earth)!
- The power supply plug serves as a safe disconnecting device from the line and must always be easily accessible.
- Place the unit on an even surface on a base made of nonflammable material.
- Do not stay in the area below the unit.
- Make sure you read and understand all instructions and safety precautions listed in this manual before installing or operating your unit.
- Adjust excess-temperature safety device below the flash point of the bath fluid.
- Observe the limited working temperature range when using plastic bath tanks.
- Never operate the unit without bath fluid in the bath.
- Pay attention to the thermal expansion of bath oil during heating to avoid overflowing of the fluid.
- Prevent water from entering the hot bath oil.
- Do not drain the bath fluid while it is hot! Check the temperature of the bath fluid prior to draining (e.g., by switching the unit on for a short moment).
- Use suitable connecting tubing.
- Avoid sharp bends in the tubing, and maintain a sufficient distance from surrounding walls.
- Make sure that the tubing is securely attached.
- Regularly check the tubing for material defects (e.g., for cracks).
- Never operate damaged or leaking units.
- Always turn off the unit and disconnect the mains cable from the power source before performing any service or maintenance procedures, or before moving the unit.
- Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit.
- Always empty the bath before moving the unit.
- Transport the unit with care.
- Sudden jolts or drops may cause damage in the interior of the unit.
- Observe all warning labels.
- Never remove warning labels.
- Never operate units with damaged mains power cables.
- Repairs are to be carried out only by qualified service personnel.



Some parts of the bath tank and the pump connections may become extremely hot during continuous operation. Therefore, exercise particular caution when touching these parts.



#### Caution:

The temperature controlling i.e. of fluids in a reactor constitutes normal circulator practice.

We do not know which substances are contained within these vessels. Many substances are:

- inflammable, easily ignited or explosive
- hazardous to health
- environmentally unsafe

#### i.e.: dangerous

The user alone is responsible for the handling of these substances! The following questions shall help to recognize possible dangers and to reduce the risks to a minimum.

Are all tubes and electrical cables connected and installed?
 Note:

sharp edges, hot surfaces in operation, moving machine parts, etc.

- Do dangerous steams or gases arise when heating? Is an exhaust needed when working?
- What to do when a dangerous substance was spilled on or in the unit? Before starting to work, obtain information concerning the substance and determine the method of decontamination.




Notice: Check the safety installations at least twice a year!

- Excess temperature protection according to IEC 61010-2-010.
   With a screwdriver turn back the adjustable excess temperature protection until the shut-down point (actual temperature).
- Low level protection according to IEC 61010-2-010. To check the function of the float, it can be manually lowered with a screwdriver for example.

#### **Operating controls and functional elements** 4.


#### 4.1. Circulator

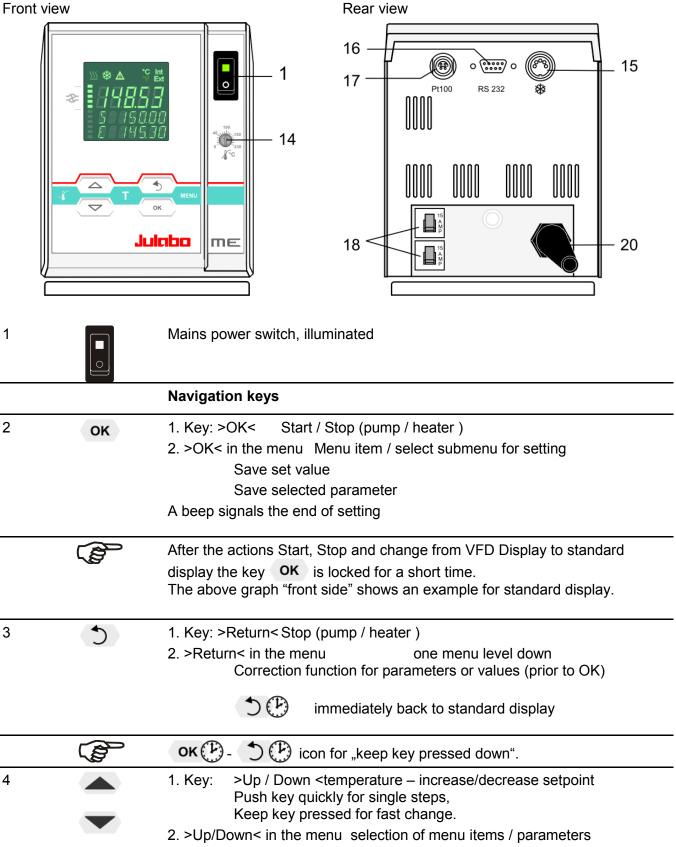

		Menu keys
5		Key: start the menu > warning and safety values<
6		Key: start the menu >temperature setpoints<
7	MENU	Key: display of MENU structure
10	S 15000 E 145.30	VFD COMFORT-DISPLAY         Header: Control indicators see sections 11 and 12         Line 1:       Actual value internal or external         The display is depending on the selected control mode in the menu > Control < (internal or external).
11	<u>₩</u>	<u>Control indicators in the header:</u> Heating / Cooling / Alarm / <b>R</b> emote control
12	°C Int °F Ext	Control indicators in the header: Temperature indication <b>Int</b> ernal or <b>Ext</b> ernal actual value Temperature indication in °C (°F not possible on this unit)
13		Display of set pump pressure stage Four stages, can be set via the key . , under >MENU - PUMP<.
14	100 40 0 230 230	Adjustable excess temperature protection according to IEC 61010-2-010
15	© *	Socket: control cable of JULABO refrigerated circulator or output for alarm messages
16	ہ (ﷺ) ہ RS232	Interface RS232: remote control via personal computer
17	()) Pt100	Socket for external measurement and control sensor or external setpoint programming
18		Mains circuit breakers (resettable) 15 A
20		Mains power cable with plug

#### 4.2. **Cooling Machine**

Front view

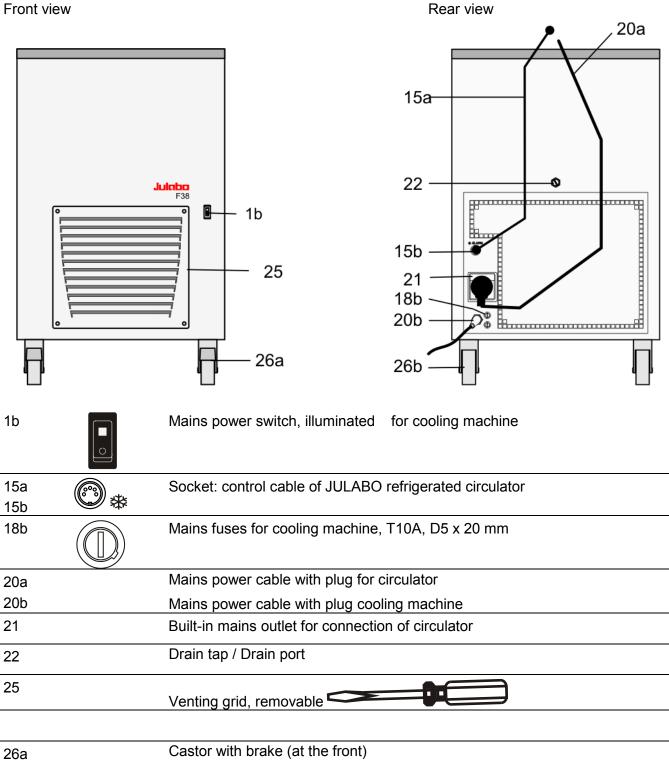

26b


Castor without brake (at the back)

## 5. Preparations

#### 5.1. Installation





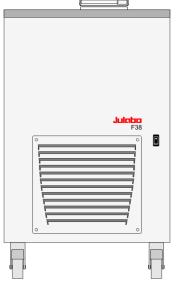

- Place the unit on an even surface on a pad made of **non-flammable** material.
  - Using the castors move the unit to the intended location.
  - The place of installation should be large enough and provide sufficient air ventilation to ensure the room does not warm up excessively because of the heat the instrument radiates to the environment. (Max. permissible ambient temperature: 40 °C). With regard to a disturbance in the cooling loop (leakage), the guideline EN 378 prescribes a certain room space to be available for each kg of refrigerant. The necessary amount of refrigerant is specified on the type plate

> For 0.52 kg of refrigerant R404A, a room space of 1 m<sup>3</sup> is required.

- F38-ME: The circulator fitted with a stainless steel bridge is placed on on the back of the bath tank leaving the bath open on the front side.
- Keep at least 20 cm of open space on the front and rear venting grids.
- Do not set up the unit in the immediate vicinity of heat sources and do not expose to sun light
- Before operating the unit after transport, <u>wait about one hour after</u> <u>setting it up.</u> This will allow any oil that has accumulated laterally during transport to flow back down thus ensuring maximum cooling performance of the compressor.



#### Caution:

Carefully read the safety data sheet of the bath fluid used, particularly with regard to the fire point!

If a bath fluid with a fire point of  $\leq$ 65 °C is used, only supervised operation is possible.

Water: The quality of water depends on local conditions.

- Due to the high concentration of lime, hard water is not suitable for temperature control because it leads to calcification in the bath.
- Ferrous water can cause corrosion even on stainless steel.
- Chloric water can cause pitting corrosion.
- Distilled and deionized water is unsuitable. Their special properties cause corrosion in the bath, even in stainless steel.

#### **Recommended bath fluids:**

Bath fluid	Temperature range
soft/decalcified water	5 °C to 80 °C
mixture water/glycol, mixture 1:1	-20°C to 50°C

#### JULABO bath fluids

JULABO	-	Thermal	Thermal	Thermal
Description		G	HY	H5
Order Number	10 liters	8 940 124	8 940 104	8 940 106
	5 liters	8 940 125	8 940 105	8 940 107
Temperature range	°C	-30 80	-80 55	-50105
Flash point	°C		78	124
Fire point	°C		80	142
Color		light yellow	clear	clear



See website for list of recommended bath fluids.

#### ATTENTION:

The maximum permissible viscosity is 50 mm<sup>2</sup> /s·



### Caution:

#### Fire or other dangers when using bath fluids that are not recommended:

Use only nonacidic and noncorrosive bath fluids. JULABO assumes no liability for damage caused by the selection of an unsuitable bath liquid.

Unsuitable bath fluids are fluids which, e.g.,

- are highly viscous (much higher than recommended at the respective working temperature)
- have a low viscosity and have creep characteristics
- have corrosive characteristics or
- tend to crack.
- No liability for use of other bath fluids!

### 5.3. Filling / draining




### Notice:

• Pay attention to the thermal expansion of bath oil during heating to avoid overflowing of the liquid.

Do not drain the bath fluid while it is hot! Check the temperature of the bath fluid prior to draining (by switching the unit on for a short moment, for example).

- Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit, or before moving the unit.
- Store and dispose the used bath fluid according to the laws for environmental protection.

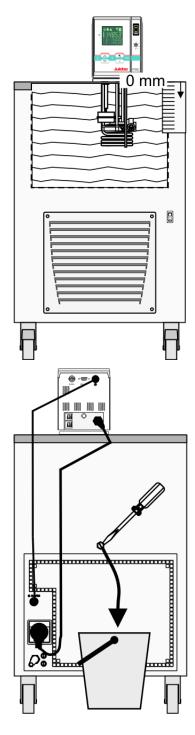

### Filling

Take care that no liquid enters the interior of the circulator.

(1) By inserting the goods there is a displacement of the bath liquid. This should be considered when filling the bath tank.

Recommendation for the use as forcing test with 20 beer bottles:

Filling level below					
the tank rim:	46	63	80	97	[mm]
when using	5	10	15	20	bottles
				<del></del> .	

Fill the bath tank according to the table (e.g. Thermal G). Put the stainless steel insert cage with the beer bottles into the bath tank. Then the filling height should be approx. 25 mm below the bath rim.

Limit value:

Recommended filling level with water as bath fluid:

max. 25 mm below the tank rim

- min. 70 mm below the tank rim (WARNING >CODE 40<)
- (1) The circulator provides an early warning system for low level that may be triggered when changing samples in the bath.

### Draining


- Turn off the circulator and cooling machine.
- Unscrew the drain tap (22) and empty the unit completely.
- Tighten the drain tap.

## 6. Operating procedures

#### 6.1. Power connection

#### Caution:

- Only connect the unit to a power socket with earthing contact (PE protective earth)!
- The power supply plug serves as safe disconnecting device from the line and must be always easily accessible.
- Never operate equipment with damaged mains power cables.
- Regularly check the mains power cables for material defects (e.g. for cracks).
- We disclaim all liability for damage caused by incorrect line voltages!

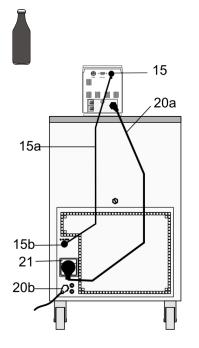

Check to make sure that the line voltage matches the supply voltage specified on the identification plate. Deviations of  $\pm 10$  % are permissible.

- Connect the circulator with mains power cable (20a) to the mains outlet (21).
- Connect the control cable (15a) between the connectors & (15, 15b).
- Connect the refrigerated circulator with mains power cable (20b) to the mains socket.

6.2. Switching on / Start - Stop

#### 6.2.1. Switching on the circulator









- Switching on:
- Turn on the mains power switch (1).
- (i) The unit performs a self-test.

Then the software version (example: V 1.xx) appears. The display **"OFF**" or **"R OFF**" indicates the unit is ready to operate.

 The circulator enters the operating mode activated before switching the circulator off: keypad control mode (manual operation)

or

remote control mode (operation via personal computer).

#### Operating procedures



#### Start:

Press **OK** key. The actual bath temperature is displayed on the VFD COMFORT-DISPLAY. The circulating pump starts with a slight delay.

#### Stop:

 Press OK key. or
 Keep D (P) key pressed.
 The VFD COMFORT-DISPLAY indicates the message "OFF".

### 6.2.2. Switching on the Cooling Machine



Switching on:

• Switch on the cooling machine using the switch (1b).



### (i) Control of the cooling machine:

With the mains switch (1b) turned on, the circulator automatically switches the cooling machine off and on.

It is switched off, if:
 the actual working temperature is increased by >30 °C (cooling is not required).

- the heater operates at full power (>800 W) for longer than 5 minutes.

• It is switched on, if:

- cooling is necessary for maintaining the bath temperature. After switch-off, the cooling machine automatically switches on only after a delay of 5 minutes for protecting the cooling compressor.

## 7. **T** Setting of temperatures

## 7.1. Using the pre-settings in the **T** menu

Pr

Ĵ	7	
		Ì

	_							
ress the `		key to	call up	the	menu	for	temperature	selection.

3 different working temperatures can be adjusted. Their values are freely selectable within the operating temperature range.

- (i) The temperatures can be set in start or stop mode.
- Press key if a value is to be retained

#### Setting of working temperature in the

1. Press the key  $\frown$  . The value flashes  $\parallel$ 

Werkseinstellungen:				
SETPNT 1	25 °C			
SETPNT 2	37 °C			
SETPNT 3	70 °C			

- 2. Select SETPOINT 1 or 2 or 3 using the key 💙 or 🔺
- 3. Confirm by pressing the **OK** key.
- (i) The circulator uses the new working temperature value for temperature control.

# Example: Adjustment/modification of the pre-setting of "SETPOINT 3"

- 1. Press the **T** key.
- 2. Select SETPOINT 3 by pressing the key. Example: SETPNT 3 / 70.00 °C
- 4. Adjust value by pressing the key and the key to 85.00 °C and confirm by pressing the OK key. The decimal digits flash '' and can be adjusted if desired. Confirm once more by pressing the OK key. Example on the left: SETPNT 3 / 85.00.
- If the active setpoint (SETPNT) is changed, the new value is immediately used for the control of the working temperature. The heater control indicator flashes.
- If the other two setpoints (not activated for control) are changed the
   MENU has to be left by pressing the key after the decimal digits have been confirmed

Notice: Refer to SETPOINT MAX / MIN in chapter 9.8. MENU LIMITS

88.8.88
SETPNT
з 10.00 ж





### 7.2. Direct setting of temperatures



SETPNT 1 2500 XK SETPNT 5 5000 The circulator uses the setpoint of SETPNT 1 or 2 or 3 for temperature control

The indicated setpoint temperature can be changed directly any time. Example: change 25.00 °C to 50.00 °C

- 1. By pressing the key the circulator switches to the active SETPOINT< example on the left: >SETPNT / 1 25.00°C<. The integer digits flash \\ (example: <25>).
- 2. By pressing the keys and the value is changed to 50.00 °C and

is confirmed by pressing the **OK** key. The decimal digits flash and can be adjusted if desired. Confirm once more by pressing the **OK** key.

- (1) The circulator uses the new working temperature value for temperature control.
- ① The temperatures can be set in start or stop mode.

## Safety installations, warning functions



Check the safety installations at least twice a year! Refer to (page 16)

SECVAL (Security Values)

> SAFETMP

8.

- > AL-TYPE
- OVERTMP
- > SUBTEMP

Settings for the excess temperature protection > **SAFETMP**< and for the warning functions for high > **OVERTMP**< and low >

**SUBTEMP**< temperature are made in a menu which is called up by pressing the key

- Menu item > **AL-TYPE**< allows choosing between a warning and
  - an alarm cut-off for the menu items > **OVERTMP**< and > **SUBTEMP**<.

#### 8.1. Excess temperature protection



### Warning:

Adjust excess-temperature safety device below the flash point of the bath fluid. In case of wrong setting there is a fire hazard! We disclaim all liability for damage caused by wrong settings!



### Notice:

On the forcing test avoid an alarm CODE 14. The program is cancelled. On the restart the count of the cycles (CYCLE) restarts from zero. >**SAFETMP**< einstellen auf 70 °C





This excess temperature protection is independent of the control circuit. When activated heater and circulating pump are completely shut down. The alarm is indicated by optical and audible signals (continuous tone) and the error message "ALARM-CODE **14**" appears on the VFD COMFORT-DISPLAY together with the ticker: > EXCESS TEMPERATURE PROTECTOR ALARM-CHECK ADJUSTMENT <

Setting range: 20 °C ... 230 °C





SRFEIME

(i) Rough setting can be effected by using the temperature scale.

### Exact setting:

- Press the key to display menu >SAFETMP<.</li>
- 2. Press the **OK** key and the set shutdown value is indicated.

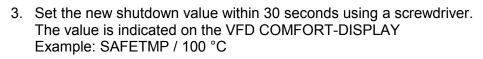

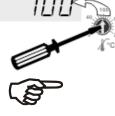

Recommendation:


Set the excess temperature protection at 5  $^\circ\text{C}$  to 10  $^\circ\text{C}$  above the working temperature setpoint.

### 8.1.1. Early warning system, low level protection



This low level protection is independent of the control circuit and is divided into two sections:

 Switch in stage 1 recognizes a defined fluid level ⊕. An audible warning sounds (interval tone) and together with the ticker: > LOW LEVEL WARNING-FILL MEDIUM < a message appears on the VFD COMFORT-DISPLAY:

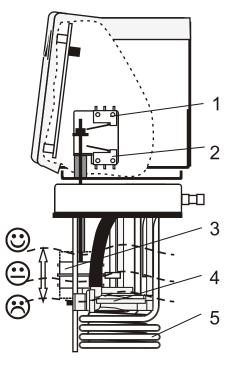

**HARNING** URRNING EDJE 40 Refill the bath fluid!

 Switch in stage 2 recognizes a low fluid level .
 If stage 2 of the low level protection according to IEC 61010-2-010 is triggered, a complete, all-pole shutdown of heater and circulating pump is effected
 A continuous alarm sounds and together with the ticker: > LOW LEVEL ALARM-FILL MEDIUM < a message appears on the VFD COMFORT-DISPLAY:

RI RRM ΓΠΤΓΓ

Turn off the unit with the mains switch, refill bath fluid and turn the unit on again!

- 3. Float
- 4. Circulating pump
- 5. Heater



### Notice:

On the forcing test avoid a low level alarm CODE 1. The program is cancelled. On the restart the count of the cycles (CYCLE) restarts from zero. Refill the bath liquid **in time**!



### Warning:

When adding bath fluid, always us the type of fluid which is identical with the fluid in the bath.

Bath oils must not contain any water and should be pre-heated approximately to the current bath temperature! Explosion hazard at high temperatures!

### 8.2. Switch-over from warning to shutdown function

ЪE	Ľ	ľ	RL
RL		T	ΥΡΕ

If a shutdown of functional elements (e.g. heater, circulating pump) is required when the limit values are exceeded or undercut the circulator can be changed over from warning function >WARNING< to shutdown function >ALARM<.

Factory setting: >WARNING<

AL - TYPE

AL - TYPE

AI ARM

WARNING XX

1. Press the key

- 2. Select the menu >SECVAL -AL-TYPE< by pressing the **v** key.
- 3. Press the **OK** key and the set parameter will flash  $\sum_{i=1}^{M}$  (Example: WARNING)
- 4. Change the parameter by pressing the very key and confirm by pressing the **OK** key.

or

press the 💙 key if the parameter is to retained.

#### • Setting >WARNING<

A mere warning function with optical and audible warning signal (interval tone) A message appears on the VFD COMFORT-DISPLAY:

888.88	
WARNING EDIE D3 OVERTMP	or

WRRNING EDIE DY SUBTEMP

Setting >ALARM<</p>

Temperature limit with shutdown of heater and circulating pump. An audible alarm sounds (continuous tone) and a message appears on the VFD COMFORT-DISPLAY:

- **[]FF-**RLRRM EDIE D3 or OVERTMP

- **[]FF**-RLRRM EDJE D4 SUBTEMP

#### 8.3. Over and Sub temperature warning function

(See chapter 8.2.

Over temperature OVERTMP חחחק

Sub temperature SUBTEMP - 99.00

test

°C 150 SETPNT OVERIMP 8500 8 7.DD 87 85 83 Recommendation: SUBIEMP Settings of the forcing 8300 20 OVERTEMP 65,0 °C SUBTEMP -5.0 °C

)

- 1. Press the key
  - 2. By pressing the or key select the menu >OVERTMP< or</p> >SUBTEMP<.

If the observance of a working temperature value >SETP< has to be supervised for a sensitive temperature application, then set over and sub

temperature warning values. In the example below the SETPOINT 85 °C is

electronics immediately register if the actual temperature breaches one of the set limit values. The resulting reaction is defined in a further menu item.

surrounded by the values OVERTMP 87 °C and SUBTEMP 83 °C. The

- 3. Press the **OK** key. The integer digits flash
- 4. Change the values to 87. °C and/or 83. °C by pressing the 💙 and **A** key and confirm with the **OK** key. The decimal digits flash and can be adjusted if desired. Confirm once more by pressing the **OK** key. See above examples.
- (1) The warning functions are only activated if the actual bath temperature remains within the set limit values for 3 seconds after switch-on.



#### **Recommendation:**

Set the over temperature warning value >OVERTMP< 5 °C to 10 °C above the working temperature setpoint.

Set the sub temperature warning value >SUBTMP< 5 °C to 10 °C below the working temperature setpoint.



## 9. MENU Menu functions

## MENU 🗸

The term "Menu functions" refers to settings such as

Menu	level	1
INICI IU	10,001	

MENU P-STRRT		Page 32	
MENU PROGRAM	Administration and creation of programs Mode Forcing test Mode Standard	Page 35	
MENU PUMP	Electronically adjustable pump capacity	Page 38	
MENU EONFIG	Configuration of the unit REMOTE – on / off (remote control via RS232) AUTOST – AUTOSTART on / off OFF-MODE – pump on / off TIME / DATE – setting time and date RESET – factory settings		
MENU EONTROL	Control characteristics and parameters C-TYPE – Internal or external control DYNAMICS - internal Control parameter - XP-, TN-, TV- INTERNAL Control parameter - XP-, TN-, TV- XPU-, EXTERNAL	Page 47	
MENU SERIAL	Adjustable interface parameters BAUD RATE, PARITY, HANDSHAKE	Page 51	
MENU RTE	ATC - Absolute Temperature Calibration, Page 52 Sensor calibration INTERNAL SENSOR, Sensor calibration EXTERNAL SENSOR 3-point calibration		
MENU LIMITS	Limitations of temperature and capacity Page 58 SETPOINT MAX / MIN - Maximum and minimum setpoint HEAT MAX – Set maximum heating COOLING MAX – Set maximum cooling INTERN MAX / MIN – Limitation of the temperature range BAND HIGH / LOW – Band limit		
	1. Open the menu by pressing the <b>MENU</b> key.		
	2. Use the keys to scroll in menu level 1.		
	3. Press the <b>OK</b> key to change to menu level 2.		
	Press the Constraints are to be retained.		

### 9.1. MENU PROGRAM – START

#### 9.1.1. START Mode Forcing test

	This menu will start a	previously set program.	
Start-Menu	2. Return to the	ram. (refer to chapter >MENU PROGRAM<) Start-MENU and confirm the desired setting of each ith the key <b>OK</b>	
		key if a parameter is to be retained. on for parameters or values (prior to OK)	
Menu level 1	> RUNS < Number of repetitions of the warm/cool phase 1 10		
MENLI	> GO < Start the forcing test (NOW)		
P-STRRT			
ок 🕹			
Level 2	Parameter level		
P-START RUNS or →	RUNS	<ul> <li>Set number of runs with  and  oκ and  oκ</li> <li>example: 1 run</li> </ul>	
	' /IN		
Р-START 60 ок →	60 NDW 米	<ul> <li>Confirm &gt;NOW&lt; with the ok key and the program will start immediately</li> </ul>	
	l		

### 9.1.2. START Mode Standard

This menu will start a previously set program.

(i) Requirements:

3. Create a program. (refer to next chapter)

Start-Menu

- Return to the Start-MENU and confirm the desired setting of each MENU item with the key Οκ
- 5. Set a start time (>TIME< >DATE< >YEAR<) if the program is to be started by the internal timer.

Menu level 1	> STEP<	Program start at section 1 10
MENLI	> RUNS <	Number of repetitions 1 99
P-START	> END<	Status at end of program (STDBY/SETPNT) Standby or last setpoint
ок 🕹	> GO <	Time of start (NOW/TIMER)

Level 2	Parameter level	<ul> <li>Press the  key if a parameter is to be retained.</li> <li>Correction function for parameters or values (prior to OK)</li> </ul>
Р-5ТАРТ 5ТЕР ок →	STEP / Ж	• Set program step with and ok example: STEP 1
P-START RUNS or →	RUNS / XK	• Set number of runs with and ok example: 1 run
P-START ENII ok →	ENI STILY or ENI SETPNT	<ul> <li>Set desired parameters with and ok.</li> <li>(STDBY / SETPNT) Parameter STanDBY: the circulator switches to – OFF Parameter SETPoiNT: the circulator constantly keeps the temperature at the value of the last step.</li> </ul>
Р-STЯRТ 60 ок →	GD NDW XK or GD TIMER	<ul> <li>Confirm &gt;NOW&lt; with the OK key and the program will start immediately</li> <li>Or start at the set time under parameter (TIMER ). Set time in the example below: 09. August 2009, 11:15 hrs</li> </ul>
GD TIMER OK ↓		③ set the time for the start of the program in the submenu >TIMER<.
Submenu TIMER	Parameter level	
TIMER TIME oK →	TIME ILIS Ж	<ul> <li>&gt;TIME&lt; hours/minutes (hh:mm), set both values one after the other and confirm</li> <li>hours flash, set by pressing         <ul> <li>+ οκ</li> <li>minutes flash, set by pressing</li> <li>+ οκ</li> </ul> </li> </ul>
TIMER DRTE οκ→	JRTE 09.08 XK	<ul> <li>&gt;DATE&lt; day/months (TT/MM), set both values one after the other and confirm.</li> <li>day flashes, set by pressing <ul> <li>+ ok</li> <li>month flashes, set by pressing</li> <li>+ ok</li> </ul> </li> </ul>

#### Menu functions

TIMER YERR οκ→	YERR year 2009 XX Set the year with ▲ ▲ ▲ ■ and ○ĸ.
TIMER START ok →	$STRRT$ $\square K$ <
<b>- []FF -</b> S 2500 TIMER	<ul> <li>Display of time until start: In line 3 the notice &gt;TIMER&lt; and the set values for "TIME" and "DATE/YEAR" are alternately indicated</li> <li>Check the correct setting of the internal real time clock if required (see MENU CONFIG)</li> </ul>
$ \begin{array}{c} \mathbf{P} \mathbf{P} \mathbf{P} \mathbf{P} \mathbf{P} \mathbf{P} \mathbf{P} \mathbf{P} $	The started program         After the start the program will indicate the currently calculated setpoint in line         2         S XX.XX. The value increases within the time period >TSLICE< until the target temperature >SETPNT< of the section is reached.
<b>- []F F -</b> 5 x x x x 1 x x x x	<ul> <li>Termination / Interruption of a program</li> <li>The program can be terminated any time by pressing the key</li> <li>In case of power failure the program is interrupted. The circulator switches to –OFF</li> </ul>

- The circulator switches to -OFF- .
- (i) If the AUTOSTART-function is activated the programmer starts again at the point in time approx. 5 minutes prior to the interruption. However, an uncontrolled change of the bath temperature has occurred.

## 9.2. MENU PROGRAM – creation, administration

	PRDGRAM MDDE Progr Set this menu item fir	ram Mode r <b>st!</b>
	and the "Forcing Test-F	which submenu items are available in >PROGRAM
Menu level 1		
MENU PROGRAM	<ul> <li>Press  key, if a parameter is to be retained.</li> <li>Correction function for parameters or values (prior to OK)</li> </ul>	
ок 🕁		
Level 2	Level 3	
PROGRAM E DI T		
PROGRAM Delete		
Р <i>R</i> ОБ <i>R</i> ЯМ МО∏Е ок →	MD IE STANIRI XX	<ul> <li>Set program mode with and ок</li> <li>FORCING-Test Program page 35</li> </ul>
	MD DE FOREIER	STANDARD-Program page 37
9.2.1. FORCING-1	Fest Program	
	>EDIT< PROGF	RAM – creation, administration
	> W-TE	MP< Temperature of the warm phase Adjustable from –40 °C up to +80 °C
	>W-TIN	
	> C-TE	-
	>C-TIM	
	> DELETE< without function when forcing test program is set	
	>MODE< Program	n Mode

#### Menu functions

Menu level 1 MENU PRDGRAM		a parameter is to be retained. for parameters or values (prior to OK)
Level 2	Level 3	Parameter level
PR⊡GRAM E III ▼	E IIII W-TEMP ok →	<ul> <li>SETPNT EDDD ;;;</li> <li>Set temperature setpoint of the warm phase (°C).</li> <li>Integer digits flash, set by pressing</li> <li>Decimal digits flash, set by pressing</li> </ul>
		+ ок
	ЕЛІІ И-ТІМЕ ок →	ISLICE 24.00 X (time slice)
		Set duration of the warm phase (h.Min) <ul> <li>Set duration by pressing</li> <li>аnd ок</li> </ul>
	Е ∏ I I E - TEMP ок →	Set temperature setpoint of the cool phase (°C). Integer digits flash, set by pressing Decimal digits flash, set by pressing + OK
	ЕШІІ Е-ТІМЕ ок →	TSLIEE 2400 X (time slice)
		Set duration of the cool phase (h.Min) <ul> <li>Set duration by pressing</li> <li>and ок</li> </ul>
PRDGRRM DELETE oK →	IELETE STEP	Delete program     without function when forcing test program is set
РКОБКАМ МОЛЕ ок 🔸	MDIE FOREIER <mark>X</mark> S	


#### 9.2.2. STANDARD Program

Menu level 1 MENU PRD5RRM

1 program

10 sections

The integrated programmer permits fast and easy programming of setpoint temperature sequences. This temperature sequence is called program. A program is composed of individual sections (STEP). The sections are defined by duration (TSLICE) and target temperature. The target temperature is the setpoint (SETPNT), which is achieved at the end of a section. The programmer calculates the temperature ramp from the difference in time and temperature.

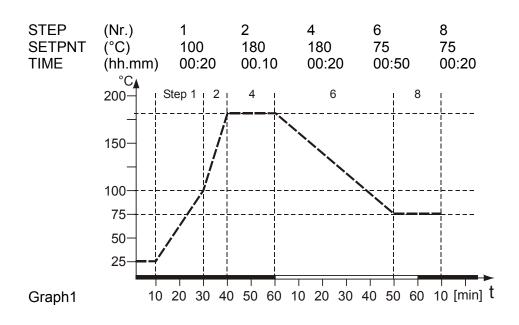

③ Sections without set value and time are skipped. They can be defined retroactively and the integrated into the program.

Menu level 1 MENU PREGRAM		Create, administer program > STEP< Program step (1 10) >SETPNT < Temperature setpoint of step >TSLICE< Duration of step delete program step (01 10, ALL) a parameter is to be retained. for parameters or values (prior to OK)
Level 2	Level 3	Parameter level
РПБЛЯМ ЕЛІТ •	ЕЛІІ STEP ок >	<ul> <li>Set program step with</li> </ul>
	EDIT DI STEP 🗲	and OK
	STEP	<ul> <li>← (Example: EDIT STEP 01)</li> <li>④ For STEP 01 the values for SETPOINT 01 and</li> </ul>
		TSLICE 01 are set one after the other
	EDIT DI SETPNT OK →	SETPNT IDDDD ): (values within working temp. range)
		<ul> <li>Integer digits flash, set by pressing</li> <li>+ οκ</li> </ul>
		<ul> <li>Decimal digits flash, set by pressing</li> <li>+ οκ</li> </ul>
	EIII II ISLIEE OK >	TSLIEE DD. ID XX
	(time slice)	<ul> <li>Set duration by pressing</li> </ul>
PROGRAM DELETE OK -	IELETE STEP Ж	<ul> <li>Delete program Program steps can be deleted individually or entirely. (STEP 01, 02, 10, ALL).</li> <li>Set parameters by pressing and ок</li> </ul>
РКОБКАМ МОЛЕ ок →	MDIE STRNIRI <mark></mark> ‰⊅	

#### 9.2.3. The easy way to the forcing test



#### Notice:

On the forcing test avoid low level alarm CODE 1. The program is cancelled. On a restart the count of the cycles (CYCLE) restarts from zero. Refill the bath liquid in time!

- 1. **Installation** see page 20.
- 2. **Power connection** see picture "Rear view" page 23
- 3. Filling / Bath fluid Filling: see page 22
- **4. Insert the the product** Put the stainless steel insert cage with beer bottles into the bath tank and refill the fluid. Maximum filling height approx. 25 mm under the bath edge.
- Put on the bath cover In order to reduce the loss of fluid due to evaporation of water, in case of long-term operation.
  - (i) The circulator ME is provided with an early-warning system for low level, which activates at a filling height of 65-70 mm below the bath edge. For measuring cycles, which last for several days the liquid level should be peered once a day, in order to not to endanger the course of the forcing test.

	6.	<b>Switching on:</b> Circulator and cooling machine may be turned on and off with separate mains switches (1a, 1b).
SAFETMP 700-00-00-00-00-00-00-00-00-00-00-00-00-	7.	Safety installations, warning functions Set the following value in the MENU Excess temperature protection see page 28 Set >SAFETMP< to 70 °C
		Over and Sub temperature warning functions see page 30DVERTMP 500SUBTEMP - 500Set > SUBTMP< to -5 °C
- []FF- SETPNT 1 2000	8.	Setting the test temperature Select e.g. SETPNT 1 in the MENU and set 20 °C. see page 25

Menu functions	
мели МЕЛЦ РКПБКЯМ ок →	9. Set MENU Program to >Mode Forcier< See page 35 PROGRRM MODE → MODE FOREIER ;;; ок
	10. Start the forcing test
MODE	As factory-setting the "Forcing test 60 °C" is stored
FORTIER	Measuring cycle:RUNS1 flow(48 h)Warm phaseW-TEMP60 °CDurationW-TIME24 hCool phaseC-TEMP0 °CDurationC-TIME24 h
	Setting of other values see page 35 "FORCING-Test Program "
Start-Menu MENU ↓ Menu level 1 MENU P STHRT ok ↓	<ul> <li>Press the  key if a parameter is to be retained. Correction function for parameters or values (prior to OK)</li> <li>RUNS &lt; Number of repetitions of the warm/cool phase 1 10</li> <li>GO &lt; Start the forcing test (NOW)</li> </ul>
Level 2	Parameter level Level 2
P-STRRT RUNS or →	RUNS P-START RUNS OK >
Р-STRRT БО ок →	60 NDW;;; 80 ок →

11. **Program sequence** A program sequence is described from page 41 on.

#### Program sequence:

#### Example forcing test 60 °C

As factory-setting the "Forcing Test 60 °C" is stored

RUNS	1 repetition
W-TEMP	60 °C
W-TIME	24 h
C-TEMP	0 °C
C-TIME	24 h
	W-TEMP W-TIME C-TEMP

In the course of the program the VFD-DISPLAY shows messages concerning the process of the forcing test in line 2 and 3.


In line 3 two messages are shown alternately.

Message 1 always indicates the program status (phase of the forcing test), message 2 indicates the status >RUN< (Test runs).

Message 2 can be changed by the editing keys . The remaining time of a phase or >E -----< can still be chosen.

(Message >E ------<, temperature value of the external temperature sensor is not important for the forcing test.)

Action diagram:

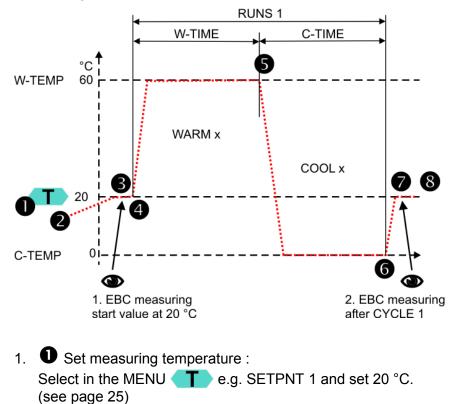


2. Start forcing test (see page 32)

Р-START 60 60 NDW Жок

#### Menu functions

5 2000 STARTUP 5 2000 8 2000 RUN	Beginning of the STARTUP-phase First the circulators tempers to the measuring ten Line 2: S 20.00 (SETPNT 1) is shown (Example Line 3: Message 1 >STARTUP<, Message 2 >RU or >00h00.00<. (The time of the start phase is not defined).	: 20.0 °C)
S 2000 STARTUP S 2000 MEASURE	The measuring temperature is reached. Measurin Line 3: Message 1 >STARTUP<, Message 2 >MI Additionally the signal tone sounds, which asks for in order to enter the start value.	EASURE<
S 6000 WRRM I S 6000 S 6000 S 6000 S 6000	Der Thermostat heizt. Die Zeit, >W-TIME< wird a Press the key ok . Beginning of the warm phas Line 2: S 60.00 (W-TEMP) is shown (Example: Line 3: Message 1 >WARM 1<, Message 2 >RUI or >23h58.23<. The circulator is heating. The time >W-TIME< is a	se. 60 °C) N<
2001 2001 2001 2000 2000 2000 2000	The time >W-TIME< is over. Direct beginning of t Line 2: S 0.00 (C-TEMP) is shown (Example: 0 Line 3: Message 1 >COOL 1<, Message 2 >RUN or >21h18.06<. The circulator is cooling. The time, >C-TIME< is o	).0 °C) I<
5 2000 WRRMUP 5 2000 RUN	The time >C-TIME< is over. Direct beginning of the WARMUP-phase Line 2: S 20.00 (SETPNT 1) is shown (Example Line 3: Message 1 >WARMUP <, Message 2 >R or >00h00.00<. (Time of the WARMUP -Phase is not defined.)	
S 2000 EYELE I S 2000 MERSURE	The measuring temperature is reached. Line 3: Message 1 >CYCLE 1<, Message 2 >ME Additionally a signal tone sounds, which asks for The number of repetitions (RUNS) form <b>one</b> (CYCLE). In the example the number of >RU >1<,due to which already after a warm- cool preached and the demand for measuring is eff	measuring measuring cycle NS< has been set to phase >CYCLE 1< is
5 6000 WARM 2 5 000 COOL 2 5 2000 CYCLE 2	In order start the second measuring cycle for the ok has to be pressed. The program skips to position 4. Direct beginning With each cycle warm phase and cool phase are	of the warm phase.

#### 9.3. MENU PUMP – Setting of pump pressure

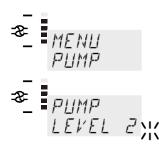

The capacity of the circulating pump is set by adjusting the motor speed Settings: stage / LEVEL 1 ... 4

Display: 🍣 with illuminated indicator

 Flow rate:
 11 ... 16 l/m

 Pump pressure:
 0,22 ... 0,45 bar

Factory setting: stage 2

РЦМР	
LEVEL	1

- 1. Press the **MENU** key.
- Select the menu >PUMP< pressing the key and confirm by pressing the OK key The set parameter flashes (example: >LEVEL 2<)</li>
- Change the parameter by pressing and confirm by pressing the OK key.


or

2

Press the 🕥 key if the parameter is to be retained.

The pump flow is pre-adjusted in the factory and can be modified to suit user requirements.

- Using a screwdriver turn the screw (1) anti-clockwise by 360 °.
- Using flat pliers turn the marking of the slide (2) to the desired position.
- Tighten the screw.

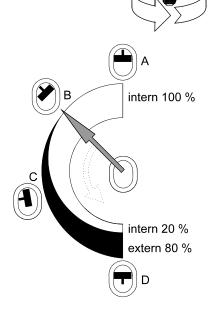

Examples:

Internal applications in the bath

- A 100 % internal bath circulation (for large bath tanks)
- B Reduced internal bath circulation (for smooth surface of bath fluid)

#### External/internal applications

- C 40 % external discharge,60 % internal circulation(for large bath tanks)
- D 80 % external discharge,20 % internal circulation(for small bath tanks)

## 9.4. MENU CONFIG – Configuration of unit

Menu level 1 MENU E DNF I 5 ok		ffected only in the >OFF< mode. $\overrightarrow{RESET}$ to by pressing the <b>OK</b> key and call up the menu
Level 2	Parameter level	<ul> <li>Press the key if a parameter is to be retained. Correction function for parameters and values (prior to OK).</li> </ul>
EONFIG	REMOTE OFF XX or REMOTE ON RUTOST	<ul> <li>Switch on and off remote control by pressing and ок</li> <li>Control display in the topline</li></ul>
	or RUTOST DN	AUTOSTART on = on AUTOSTART off = off See WARNING page 45
E DNFIG DFFMD]]E ok →	DFFMDJE PMP DNX or DFFMDJE PMP DFF	<ul> <li>Switch on and off OFFMODE by pressing and ok</li> <li>PUMP ON continuous operation of circulating pump</li> <li>PUMP OFF circulating pump is linked to Start/Stop</li> </ul>
	Level 3	Parameter level
E⊡NFIG TIME/∐T ок →	TIME / IT TIME OK >	TIME TIME X
		<ul> <li>Hours flash, set by pressing + οκ</li> <li>Minutes flash, set by pressing + οκ</li> </ul>
	ТІМЕ/ПТ ПЯТЕ ок →	<ul> <li>IRTE 08.09 \;</li> <li>Day flashes, set by pressing ▲ ▼ + ок</li> </ul>
		<ul> <li>Month flashes, set by pressing</li></ul>

-		
	Level 3	Parameter level
	TIME∕∐T YERR οκ→	<ul> <li>YERR</li> <li>• Year flashes, set by pressing + ок</li> </ul>
E ⊡NFIG RESET ок →	- []FF- RESET DK ]:: - []FF- RESET - RUN-	<ul> <li>Return to factory settings by pressing ок RESET returns all set values to the factory setting except for date and time.</li> <li>A RESET can be effected only in the –OFF- mode.</li> <li>During the message –RUN- all parameters are reset to factory settings.</li> </ul>

#### 9.4.1. REMOTE

Factory setting: OFF	<ul> <li>The control electronics offer two ways of adjusting a setpoint.</li> <li>Adjustment of setpoint using the keypad or the integrated programmer.</li> <li>Adjustment of setpoint via the serial interface RS232 using a PC or a superordinated process control system.</li> <li>The topline of the VFD-DISPLAY shows a bright "<b>R</b>"FOR remote control; - remote control discontinued.</li> </ul>
o  o RS232	IMPORTANT: additional measures for remote control  Connect the circulator to the PC using an interface cable.
	<ul> <li>Check the interface parameters of both interfaces (circulator and PC) and make sure they match. (refer to 12.1. Setup for remote control page 65)</li> </ul>

#### 9.4.2. AUTOSTART



#### Warning

For supervised or unsupervised operation with the "AUTOSTART" function avoid any hazardous situation to persons or property

Take care to fully observe the safety and warning functions of the circulator.

Factory settings: OFFNotice:<br/>The circulator has been configured and delivered by JULABO in<br/>accordance with the NAMUR recommendations. This means for the start<br/>mode that the unit must enter a safe operating status after a power<br/>failure. This safe operating status is indicated by the message "OFF" or<br/>"R OFF" on the VFD COMFORT-DISPLAY.<br/>A complete, all-pole shutdown of the main functional elements such as<br/>heater and pump motor is effected.<br/>The values set on the circulator remain saved and the unit is restarted by<br/>pressing the start/stop key in manual control.<br/>In remote control mode the values need to be resent by the PC via the

interface.

If such a safety standard is not required, the NAMUR recommendations can be bypassed with the AUTOSTART function thus allowing a direct start of the circulator by pressing the mains switch or using a timer.

#### 9.4.3. OFF-MODE

Factory setting: Usually the circulating pump is controlled with the key **OK** or the start/stop command. If the circulating pump is to work in the –OFF-mode, the adjustment can be set in a sub-menu.

(1) The pump motor will be shutdown in case of alarm anyhow.

#### 9.4.4. Setting of clock and date



TIME

1115

The internal real time clock allows starting a program any time. The clock is set to the local mean time (MEZ) at the factory.

- (i) If the unit is operated in a different time zone, the clock can be adjusted in this menu.
- (i) Change summer/winter time in this menu

#### 9.4.5. RESET – Factory settings

A Reset will return all values to factory setting except for date and time.



A RESET can be effected in the >OFF< mode only.</li>
 Switch off the circulator by pressing the key OK and call up the menu CONFIGURATION.

Menu level 1		ed for internal and external temperature control ed out in the menu >C-TYPE< .(INT or EXT).
MENU EDNTROL	For external temperature control and measurement connect a Pt100 external sensor to the socket at the rear of the circulator.	
ок ↓	Press the ke for parameters or v	ey if a parameter is to be retained. Correction function alues (prior to OK)
Level 2	Parameter level	
EENTROL E-TYPE or →	E-TYPE INT XX	<ul> <li>Switchover of control type by pressing</li> <li>and οκ</li> </ul>
	or	① The control type can be adjusted in the -OFF-
	Ε-ΤΥΡΕ	mode only.
	E×T	Depending on the adjustment only the active parameters are displayed.
C-TYPE INTERNAL		
EONTROL DYNRMIE ok →	IYNAMIE APER Ж	<ul> <li>The parameter flashes, switch by pressing</li> <li>and οκ</li> </ul>
	or IIYNRMIE NORM	① This parameter affects the temperature sequence in case of internal control.
EENTREL XP INT OK >	×ア INT 15 );; 0.1 99.9	<ul> <li>The parameter flashes, set by pressing</li> <li>+ οκ</li> </ul>
EDNIRDL IN INT OK >	TN INT 100 XX 39999	<ul> <li>The parameter flashes, set by pressing</li> <li>+ οκ</li> </ul>
EENTROL TV INT OK ->	TV INT 5 0999	<ul> <li>The parameter flashes, set by pressing</li> <li>+ οκ</li> </ul>
C-TYPE EXTERNAL		
EDNTROL XPEXT OK >	×ア E×T ロフ);; 0.1 99.9	<ul> <li>The parameter flashes, set by pressing</li> <li>+ οκ</li> </ul>
EDNTROL TN EXT OK >	TN E×T 72日)::( 3 9999	<ul> <li>The parameter flashes, set by pressing</li> <li>+ οκ</li> </ul>

#### 9.5. MENU CONTROL – Control characteristics and parameters

Menu functions

Level 2	Parameter level	
EDNTROL TV EXT OK >	<i>TV E×T</i> 55 0999	<ul> <li>The parameter flashes, set by pressing</li> <li>         + οκ     </li> </ul>
Е ПИТ Я П ХРЦ ОК →	× <i>PU</i> ヨロ ╎╎ 0.1 99.9	<ul> <li>The parameter flashes, set by pressing</li> <li>         + οκ     </li> </ul>

#### 9.5.1. **CONTROL – Control INTERNAL / EXTERNAL**

	Switchover can only be effected if a Pt100 external sensor is connected.
Pt100	Factory setting: INT
(B)	
	<b>IMPORTANT:</b> Additional measures for external temperature control
	Suggested settings for external temperature control: BAND HIGH / LOW and INTERN MAX / MIN see chapter >LIMITS< page 58.
	Sensor calibration of the Pt100 external sensor is carried out in the menu >ADJUST<, submenu >ATC SENOR - EXT<; set ATC STATUS< to >OFF< (See page 52).



Attention: Place the external sensor into the temperature-controlled medium and securely fix the sensor.



#### Accessory: Pt100 external sensor


Order No.	Description	Material	Cable
8981003	200x6 mm Ø,	stainless steel	1.5 m
8981005	200x6 mm Ø,	glass	1.5 m
8981006	20x2 mm Ø,	stainless steel	1.5 m
8981010	300x6 mm Ø,	stainless steel	1.5 m
8981015	300x6 mm Ø,	stainless steel / PTFE coated	3 m
8981013	600x6 mm Ø,	stainless steel / PTFE coated	3 m
8981016	900x6 mm Ø,	stainless steel / PTFE coated	3 m
8981014	1200x6 mm Ø,	stainless steel / PTFE coated	3 m
8981103	Extension cable fo	r Pt100 sensor	3.5 m
8981020	M+R in-line Pt100	sensor	

Pt100

M+R

The M+R in-line Pt100 sensor is a flow sensor and can be installed loop circuit

#### 9.5.2. Dynamic internal

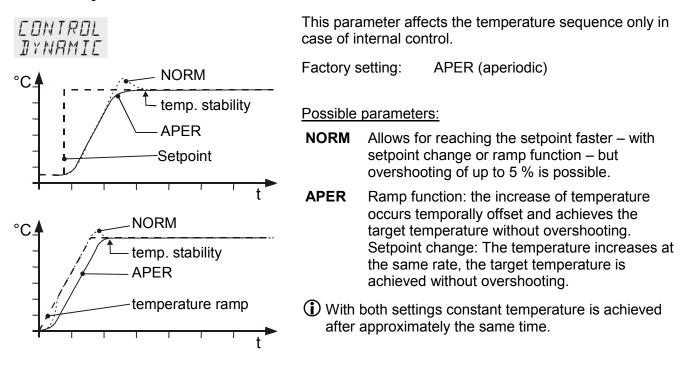

#### 9.5.3. Control parameters – XPU-, XP-, TN-, TV- EXTERNAL

In most cases the control parameters preset in the factors are adequate XP EXT for achieving an optimum temperature sequence. 0.7 The control parameters allow adjustment to special control processes. Setting range: 0.1 ...99.9 TN EXT 720 Setting range: 3 ... 9999 TV EXT 55 Setting range: 0 ... 999 Proportional range >Xpu< XPH The proportional range Xpu of the cascaded controller is only needed for 30 external control. Setting range: 0.1 ... 99.9

#### 9.5.4. Control parameters- XP-, TN-, TV- INTERNAL

In most cases the control parameters preset in the factory are adequate for achieving an optimum temperature sequence. The control parameters allow adjustment to special control processes.

# XP TNT

Proportional range >Xp<

The proportional range is the range below the setpoint in which the control circuit reduces the heating capacity from 100% to 0 %

Setting range: 0.1 ... 99.9

Reset time >Tn< (Integral component)

Compensation of the remaining control deviation due to proportional regulation. An insufficient reset time may cause instabilities. Excessive reset times will result in unnecessary prolongation of compensation of the control difference.

#### ΤĽ INT


Setting range: 0 ... 999

Setting range: 3 ... 9999

Lead time >Tv< (Differential component)

The differential component reduces the transient time. An insufficient lead time will prolong the time required for compensation of disturbance effects and cause high overshooting during run-up. An excessive lead time could cause instabilities (oscillations)

#### **Optimization instructions for the PID control parameters**

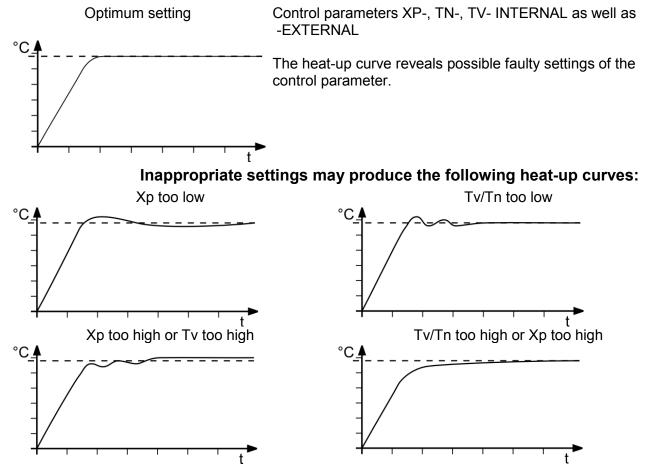

### 9.6. MENU SERIAL - BAUDRATE, HANDSHAKE, PARITY

Menu level 1 MENU SERIAL		
Level 2	Parameter level	Press the  key if a parameter is to be retained.
SERIAL BAUBRAT OK -	BAU DRAT 4800 XX BAU DRAT 9600	<ul> <li>The parameter flashes, switch by pressing</li> <li>and οκ</li> </ul>
	BRUDRAT 19200 BRUDRAT 38400	
SERIAL PARITY OK >	PARITY EVEN PARITY	<ul> <li>The parameter flashes, switch by pressing</li> <li>and οκ</li> </ul>
	ע ע	even: Data bits = 7; Stop bits = 1
	PARITY ND	odd: Data bits = 7; Stop bits = 1 no: Data bits = 8; Stop bits = 1
SERIAL HSHAKE OK →	н Shrke Нвр ј Ж	<ul> <li>The parameter flashes, switch by pressing</li> <li>and οκ</li> </ul>
	HSHRKE SOFT	Xon/Xoff-protocol (Software handshake) Protocol RTS/CTS (Hardware handshake)

#### 9.7. MENU ATC - Absolut Temperature Calibration



ATC serves to compensate a temperature difference that might occur between circulator and a defined measuring point in the bath tank because of physical properties.

Principle: For ATC calibration, in steady state the bath temperature at the location Example: of the temperature sensor (CT) is determined at the respective adjusted 1-point calibration working temperature. This value is then set on the circulator in the menu °C >ATCalibration< under menu item > CTEMP X <. This can be a 1-point, 2-point or 3-point calibration. Μ B (INT) Т T<sub>⊤</sub>1 88888 °C 3-point calibration °C CT  $\mathsf{T}\mathsf{T}$ B (EXT) T<sub>⊤</sub>1 °C  $T_T 1 = Original curve$ Pt100 ext 88888 °C CT TΤ M M = Temperature measuring instrument with temperature sensor B = Bath tank (INTernal or EXTernal) T = circulator CT = Temperature on measuring point TT = Temperature on circulator

Menu level 1		
MENLI		
ATE		
ок 🔸		
Level 2	Parameter level	<ul> <li>Press the  key if parameter is to be retained. Correction function for parameters or values (prior to OK).</li> </ul>
RTE SENSOR or 🚽	SENSOR INTERN Ж	<ul> <li>The parameter flashes, switch by pressing</li> <li>and οκ</li> </ul>
	or SENSOR	<ol> <li>On level 2 a (I) is indicated for internal or an (E) for external.</li> </ol>
	EXTERN	Example: RTE (I) RTE (E)
RIE (I) STRTUS ok →	STATUS YES Ж	<ul> <li>The parameter flashes, switch by pressing</li> <li>and οκ</li> </ul>
	or	I >NO< Carry out an ATC calibration
	STRTUS ND	<ul> <li>YES&lt; return to standard operation after calibration.</li> </ul>
<i>ПТЕ (I)</i> ТҮРЕ ок →	IYPE I-POINT Ж	• The parameter flashes, switch by pressing $\mathbf{a}_{nd}$ $\mathbf{o}_{\kappa}$
	TYPE 2-PDINT	A >1-point<, >2-point< or >3-point< calibration can be carried out.
	TYPE 3-POINT	The selected calibration is indicated on level 2 by 1 or 2 or 3.
RIE (I)	TMPVALI	The value >TMPVAL< is only indicated
IMPVALI ok →	80.00	<ul> <li>In addition the measured temperature value</li> <li>CALVAL X&lt; is saved during the next step.</li> </ul>
RIE (I) ERLVRLI ok →	ERLVALI 79.70 <b>Ж</b>	<ul> <li>Integer digits flash, set by pressing</li> <li>+ οκ</li> </ul>
		<ul> <li>Decimal digits flash, set by pressing</li> <li>+ οκ</li> </ul>
	<ul> <li>If only a 1-point cal indicated anymore</li> </ul>	ibration is carried out, the following menu items are not

Menu functions

RIE (I) TMPVRL2 οκ→	TMPV AL 2 120.00 🕤	The value is only indicated
RTE (I) ERLVRL2 ok →	EALVAL2 1 19.50 X	<ul> <li>Integer digits flash, set by pressing         <ul> <li>+ ok</li> </ul> </li> <li>Decimal digits flash, set by pressing         <ul> <li>+ ok</li> </ul> </li> </ul>
	If only a 2-point cal indicated anymore	ibration is carried out, the following menu items are not
ATE ⟨I⟩ TMPĽAL∃ ok →	TMPV AL 3 160.00	The value is only indicated
RIE (I) ERLVRL∃ οκ →	EALVAL3 159.30 <b>X</b>	<ul> <li>Integer digits flash, set by pressing         <ul> <li>• oκ</li> </ul> </li> <li>Decimal digits flash, set by pressing         <ul> <li>• oκ</li> </ul> </li> </ul>

#### 9.7.1. ATC SENSOR - INTERNAL / EXTERNAL

RTE SENSOR

SENSOR	
INTERN	
SENSOR	
EXTERN	

In the first submenu the ATC function is set for the **>INTERN**< internal or the **>EXTERN**< external temperature sensor.

Calibration can be carried out for the internal temperature sensor and for the external temperature sensor connected to the socket "ext. Pt100". The circulator is able to save both parameter sets. However only the one which has been set under menu item >ATC SENSOR < is displayed.

#### 9.7.2. ATC STATUS - YES / NO

RIE (I)	In the second submenu the ATC function for the temperature sensor selected above is activated >YES< or deactivated >NO<.
STATUS	>YES< (factory setting) The controller of the circulator uses the original
STRTUS	curve of the temperature sensor or the new curve measured during the ATC calibration.
YES	Important: Set to >NO< during the calibration process
STRTUS	
ND	>NO< An ATC calibration is to be carried out.
	Important: Set to >YES< after calibration.
	(i) In the > ATC STATUS < >YES< the ATC calibration always affects the

In the > ATC STATUS < >YES< the ATC calibration always affects the current working temperature; also the one set via interface.</p>

#### 9.7.3. CALIBRATION TYPE: 1 -/ 2 -/ 3 POINT

RIE (I)	A >1-point<, >2-point< or >3-point< calibration can be carried out.
TYPE	First geometrically define the location for calibration (measuring point CT),
TYPE I-POINT	then determine the temperature values of the calibration points. The type of calibrations also determines the number of the following pairs of values indicated on the LCD DIALOG-DISPLAY.
TYPE 2-POINT	
TYPE 3-PDINT	

TMPVALI	EALVALI
80.00	19.70
TMPV AL 2	CALVAL2
120.00	1 19.50
TMPV RL 3	CALVAL 3
160.00	159.30

#### Pairs of values:

**TMPVAL X:** Circulator temperature 1 or 2 or 3 (actual value TT) The actual temperature of the bath is simultaneously saved with the "calibration value" >CALVAL< and can be indicated for control purposes (value does not flash).

**CALVAL X:** Calibration temperature 1 or 2 or 3 (actual value CT) The "calibration value" is determined with a temperature measuring device and saved under menu item >CALVAL<. (value flashes /1)

#### 9.7.4. Example: 3-point calibration for internal control

In the temperature range from 80 °C to 160 °C the calibration curve of the temperature sensor (TT) is to be adjusted to the actual temperatures at measuring point (CT).

#### 1. Set circulator to internal control: **MENU CONTROL page 47**

Menu level 1 The type of control can be set only in the -OFF- mode. MENLI  $\Gamma - TYPF$ EONTROL EONTROL OK \_ INT X OK E - TYPEок 🔺 **)** ('F) 2. Set working temperature setpoint – SETPNT: Refer to "Direct temperature setting " page 26 By pressing the key *(* **he** circulator switches to the active 5 >SETPOINT< see example on the left: >SETPNT / 1 25.00°C<. The integer digits flash  $\cancel{1}$  (Example: <25>). Change the value to 80.00 °C by pressing the keys 💌 and 📥 and SETPN confirm by pressing the key **OK**. 8000 XX The decimal digits flash. 80.00 °C OK Confirm once more by pressing the key 120.00 °C The bath is heated up. 160.00 °C Wait for approx. 5 minutes until the temperature is constant. 3. Reading of temperature measuring device Pt100 Read the value of measuring point CT on the device and enter under menu ורפר item >CALVAL X< by using the keypad. >CALVAL 1< (79.70 °C) >CALVAL 2< (119.50 °C) >CALVAL 3< (159.30 °C) 4. Calibration Menu level 1 Press the key if parameter is to be retained. Correction function for MENLI parameters or values (prior to OK). RTE Setting is required only for the first calibration point. OK J Level 2 Parameter level SENSOR Set SENSOR INTERN: RTE INTERNX

The parameter flashes, switch by pressing

and ok

SENSOR

ок 🔺

RIE (I) STRIUS οκ→	STATUS ND \\( O	<ul> <li>An ATC calibration is to be carried out.</li> <li>Set to &gt;NO</li> <li>The parameter flashes, switch by pressing</li> </ul>
HTE (I) TYPE OK ->	TYPE 3-POINT XX	<ul> <li>and οκ</li> <li>The parameter flashes, switch by pressing</li> <li>and οκ</li> </ul>
		A >3-point< calibration is carried out.
RIE (I) IMPVRLI οκ→	TMPVALI 80.00 S	The value >TMPVAL< is only indicated In addition the measured value >CALVAL X< is saved during the following step
ΠΤΕ ⟨I⟩ ERLVRLI οκ →	EALVALI 79.70 <b>%</b>	<ul> <li>Integer digits flash, set by pressing</li> <li>(79) + οκ</li> </ul>
		<ul> <li>Decimal digits flash, set by pressing</li> <li>(70) + οκ</li> </ul>
		The first of 3 points is calibrated.
	Return to 2. Set workin	g temperature value SETPNT: 120.00 °C
RTE (I) TMPVRL2 ok →	TMPV AL 2 120.00	The value is only indicated
	ERLVRL2 1 19.50 <b>%</b>	<ul> <li>Integer digits flash, set by pressing         <ul> <li>(119) + ок</li> </ul> </li> <li>Decimal digits flash, set by pressing         <ul> <li>(50) + ок</li> <li>The second of 3 points is calibrated.</li> </ul> </li> </ul>
	Deturn to 2 pot working	a temperature velue SETENT: 160.00 °C
RTE (I) TMPI/RL∃ ok →	TMPVAL 3 IE0.00 S	g temperature value SETPNT: 160.00 °C The value is only indicated.
ATE (I) ERLVAL∃ ok →	ERLVRL3 159.30 <b>X</b>	<ul> <li>Integer digits flash, set by pressing         <ul> <li>(159) + ок</li> </ul> </li> <li>Decimal digits flash, set by pressing         <ul> <li>(30) + ок</li> </ul> </li> <li>The 3-point calibration is completed</li> </ul>
	5. Return to standar	d operation
RTE (I) STRTUS (ok →	STRTUS YESX OK	<ul> <li>Set &gt;YES&lt; after calibration. (Standard operation)</li> </ul>

#### 9.8. MENU LIMITS

Menu level 1		
MENLI		
LIMITS		
ок 🕁		
Level 2	Parameter level	Press the two key if parameter is to be retained. Correction function for parameters or values (prior to OK).
LIMITS SETMAX or →	SETMA× 20000 <b>)</b> ;;	<ul> <li>Integer digits flash, set by pressing</li> <li>+ οκ</li> </ul>
		<ul> <li>Decimal digits flash, set by pressing</li> <li>+ οκ</li> </ul>
LIMITS SETMIN or 🔺	SETMIN - 94.99 \\\	<ul> <li>Integer digits flash, set by pressing</li> <li>+ οκ</li> </ul>
		<ul> <li>Decimal digits flash, set by pressing</li> <li>+ οκ</li> </ul>
LIMITS НЕПТМП× ок →	HERTMR× IDD\X	<ul> <li>The value flashes, set by pressing</li> <li>+ οκ</li> </ul>
	0 100 %	
LIMITS ЕППLМЯ× ок →	EOOLMA× D\X	<ul> <li>The value flashes, set by pressing mit</li> <li>→ → → → → → → → → → → → → → → → → → →</li></ul>
	0 100 %	
	In case of external cont	trol these menu items are additionally indicated.
LIMITS INTMR× or →	INTMA× 20000;X	<ul> <li>Integer digits flash, set by pressing</li> <li>+ οκ</li> </ul>
		<ul> <li>Decimal digits flash, set by pressing</li> <li>+ οκ</li> </ul>
LIMITS INTMIN OK 🔺	INTMIN -9499 XX	<ul> <li>Integer digits flash, set by pressing</li> <li>+ οκ</li> </ul>
		<ul> <li>Decimal digits flash, set by pressing</li> <li>+ οκ</li> </ul>
LIMITS ∄ЯN]-Н ок⇒	BAND-H 200 XX	<ul> <li>The value flashes, set by pressing</li> <li>+ οκ</li> </ul>
LIMITS BAND-L ok →	BAND-L 200 XX	<ul> <li>The value flashes, set by pressing</li> <li>+ οκ</li> </ul>
	/ \	<u> </u>

#### 9.8.1. Limits for internal control

	Ε-ΤΥΡΕ	SETPOINT MAX / MIN – Maximum and minimum setpoint
	INT	Restriction of the adjustable temperature range
		The limitation of the operating temperature range effects the temperature
	SEIMR×	setting in the menu with the key <b>C</b> .
	200.00	Only setting of working temperatures which lie within the determined limits is possible
	5E TMIN - 50.00	Existing settings for SETPNT 1, -2, -3, as well as those for >OVERTMP< and > SUBTMP < (refer to page 30), are automatically deferred into the limit range.
		<u>Setting range:</u> -94,90 °C +200,0 °C
		SET MAX > SET MIN Interchange of values is not possible.
		Set maximum heating / cooling
	HERTMRX	The heating and cooling capacity of the unit are adjustable. 100 % corresponds to the technical specification of the equipment.
m	100	Setting range:
	EOOLMA× D	HEAT MAX – 0 to 100 % in 1 % steps COOLING MAX – 0 to 100 % in 1 % steps

#### 9.8.2. Limits for external control

:	INTERN MAX / MIN Restriction for the temperature range of the internal bath.
INTM8× 200.00 INTMIN - 50.00	<u>Setting range:</u> -94,9 °C +200,0 °C The limits INT MAX and INT MIN are only active in external control. INT MAX and INT MIN determine fixed limits for the temperature within the internal bath. The temperature controller cannot exceed these limits even if it would be necessary for achieving the temperature in an external system. Therefore it is possible that the external setpoint cannot be achieved.
	Sense of limit setting:
> SAFETMP <	Protects the bath fluid from overheating.
	Prevents an undesired alarm shutdown by the excess temperature protection - >ALARM CODE 14<. Set the value of > INT MAX at least 5 °C below the value of >SAFETMP<.
	Protects the pump motor from high viscosity of the bath fluid at low temperatures.
	For refrigerated circulators. Freezing protection when using water as bath fluid.


#### BAND HIGH / LOW – Band limitation



The band limitation is active during external control. Varied, practiceoriented setting are feasible for heat-up and cool-down phases.

Setting range: 0 °C ... 200 °C

**BAND HIGH** and **BAND LOW** allow for the limitation of the difference between the temperatures in the internal bath and the external system to any maximum value for the heat-up and cool-down phase. During the heat-up phase this difference value is always added to the actual external temperature. During the cool-down phase the difference value is subtracted.

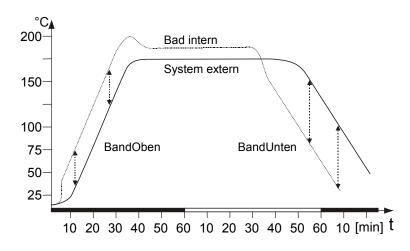

Sense of a band limitation:

- Protection of objects and samples by gentle temperature control
- Protection of e.g. glass reactors from thermal shock.

## 10. Troubleshooting guide / error messages





**BBBBB** WARNING CODE 40

#### Alarm with complete shutdown:

If one of the following failures occur a complete, all-pole shutdown of the heater and circulating pump is effected.

"A" lights up and a continuous signal sounds. The code for the cause of alarm is indicated on the VFD COMFORT-DISPLAY.

#### Alarm without shutdown:

The code for the cause of alarm is indicated on the VFD COMFORT-DISPLAY. The warning signal sounds in regular intervals. The messages appear every 10 seconds.

	Press the key ok to stop the signal
ALARM EDIE DI	Error message with ticker: >LOW LEVEL ALARM-FILL MEDIUM < Low level alarm
	The circulator is operated without or insufficient bath fluid. Switch the unit off with the mains switch, refill bath fluid and switch on!
	Tube breakage has occurred (insufficient filling level of bath fluid caused by pumping-out) Replace the tubing and refill bath liquid.
	The float is defect (e.g. transport damage). Repair by authorized JULABO service personnel.
ALARM EDIE D2	Error message with ticker: > REFRIGERATOR ALARM-CHECK CONNECTION < During the self-test after switch-on a short –circuit is detected between pin 2 and pin 4 of the control line or the control line was disconnected during operation. Reconnect the control line or repair short-circuit.
WRRNING EDIE D3 RLRRM EDIE D3	Error message with ticker: > EXCESS TEMPERATURE WARNING-CHECK LIMITS < Excess temperature warning or Excess temperature alarm <b>Type of warning:</b> set to >warning< or >alarm< (refer to page 29)
WARNING EDIE OY ALARM EDIE OY	Error message with ticker: > LOW TEMPERATURE WARNING-CHECK LIMITS < Low temperature warning or Low temperature alarm. <b>Typ of warning:</b> set to >warning< or >alarm< (refer to page 29)
ALARM EDIE OS	Error message with ticker: > WORKING SENSOR ALARM-CALL SERVICE < Cable of working temperature sensor is disconnected or short-circuited.

#### Troubleshooting guide / error messages

ALARM EDIE 06	Error message with ticker: >SENSOR DIFFERENCE ALARM-CHECK VISCOSITY AND PUMP STAGE< Defect of working or excess temperature protector. Working temperature sensor and excess temperature protector report a temperature difference of more than 35 K.
RLARM EQJE OT	Error message with ticker: > INTERNAL HARDWARE ERROR-CALL SERVICE < Other errors
ALARM EDIE 12	Error in A/D converter
ALARM EDIE IY	Error message with ticker: > EXCESS TEMPERATURE PROTECTOR ALARM-CHECK ADJUSTMENT <
	Excess temperature protector defect.
	The protection temperature is below the set working temperature setpoint. Set the protection temperature to a higher value.
ALARM	Error message with ticker: > EXTERNAL SENSOR ALARM-CHECK EXTERNAL SENSOR <
EDDE IS	External control was set but the Pt100 external senor was not connected or is defect.
WRRNING EDIE 20	Error message with ticker: > CLEAN CONDENSER OR CHECK COOLING WATER <
	Insufficient cooling of condenser. Clean the air-cooled condenser. Check the flow and the temperature of the cooling water of a water-cooled condenser.
HRRNING EDIE 21	Error message with ticker: > COMPRESSOR FAILURE-CHECK REFRIGERATOR <
	Stage 1 of the compressors does not work. Automatic restart after short cool-down, message E 21 goes off.
WARNING	Stage 2 of the compressor does not work. Cooling machine – overload protection
CO11E 22	The driving motor of the cooling compressor is equipped with an overload protection which is triggered by increased internal_temperatures or excessive current consumption.
	Shutdown can be caused by - insufficient ventilation, - insufficient wall distance,
	- soiled condenser,
	- high room temperature
	- switching off and on in short sequence
WARNING EDIE 23	Excess temperature in stage 1 of the compressor.
WARNING EDIE 24	Excess temperature in stage 2 of the compressor.
WARNING EDIE 25	Short circuit of control line to cooling machine during self-test.

ALARM COIE 33	Error message with ticker: > SAFETY SENSOR ALARM-CALL SERVICE < The cable of the excess temperature protector has been disconnected or short-circuited
WARNING EDIE YD	Error message with ticker: > LOW LEVEL WARNING-FILL MEDIUM <
	The early warning system for low level reports a critical fluid level. Refill bath fluid.
	By quickly switching off and restarting the unit the alarm is cancelled. If the error occurs once more after the restart, a remote diagnosis is required.
E-Err PRESS	Error message with ticker: > CONFIGURATION ERROR-PRESS OK<
PRESS OK	The configuration of the circulator does not correspond with its current application.
	Press the <b>OK</b> key for a non-recurring, automatic change of the configuration.
	In this case please call the JULABO Technical Service or an authorized dealer.
	Disturbances that are not indicated.
	$-\mathbf{T}$ is the state of the st

The electronic pump motor is overload-protected by an electronic current limiter. If viscosity of the bath fluid is or becomes too high, the motor stops running.



Mains circuit breakers (resettable) 15 A



Cooling machine: Fuse T 10.0 A, dia.5 x 20 mm The mains fuses (8b) on the rear of the unit may easily be exchanged as shown on the left.



#### Warning:

Before exchanging the fuses, turn off the mains power switch and disconnect the power plug from the mains socket! Only use fine fuses with a nominal value as specified.

Example:

Manufacturer	Supplier	Туре	Order No.
Wickmann	Wickmann	G- fuse insert T10,0A 5x20 mm	No. 19195


#### 11. Electrical connections



## Notice:

Use shielded cables only.

The shield of the connecting cable is electrically connected to the plug housing. The unit ensures safe operation if connecting cables with a maximum length of 3 m are used. The use of longer cables does not affect proper performance of the unit, however external interferences may have a negative impact on safe operation (e.g. cellular phones).

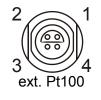

# Pt100 Shield Plug

Pin assignment: <u>Pin Signal</u>

Socket for external Pt100 sensor

Pin	Sign
1	+
2	U+
3	U-
4	I-

The shield of the connecting cable is electrically connected to the plug housing and the sensor tube.

#### **RS232 serial interface**

This port can be used to connect a computer with an RS232 cable for remote control of the circulator.

5	
9	0
RS2	32C

#### Pin assignments RS232:

Pin 2	RxD
Pin 3	TxD
Pin 5	0 V
Pin 7	RTS
Pin 8	CTS

Receive Data Transmit Data Signal GND Request to send Clear to send

Pin 1; 4; 6, 9 Reserved - do not use!

#### RS232 interface cable


Circulator (9-pol)		PC (9-pol)
Pin 2 RxD	$\Leftrightarrow$	Pin 3 TxD
Pin 3 TxD	$\Leftrightarrow$	Pin 2 RxD
Pin 5 GND	$\Leftrightarrow$	Pin 5 GND
Pin 7 RTS	$\Leftrightarrow$	Pin 8 CTS
Pin 8 CTS	$\Leftrightarrow$	Pin 7 RTS

Accessories:	Order No.	Description
	8 980 073	RS232 interface cable 9-pol./9-pol., 2,5 m
	8 900 110	USB interface adapter cable

## ₩ / Control output

The onnector may be used for control of JULABO refrigerated circulators or as output for alarm messages.

Circuit:	Operation	= relay powered
	Alarm	= relay not powered

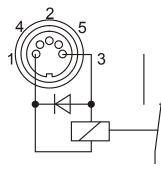

Pin assignme	ent:
Pin	Signal
1	+24 V (I max. current 25 mA)
2	0 V
3	Alarm relay
4	Reserved - do not use!
5	Cooling pulse

#### 12. Remote control

#### 12.1. Setup for remote control

5 19 RS232C 00 RS320C 00 RS3200C 00 RS	<ol> <li>Check the interface parameters for both interfaces (on circulator and PC) and make sure they match. (Serial interface refer to page 51)</li> </ol>
<b>- []FF-</b> REMOTE DN	<ol> <li>In the menu &gt; MENU CONFIG &lt; set the menu item &gt; REMOTE &lt; to &gt; ON &lt; .</li> <li>Connect both units with an interface cable</li> </ol>

Like all parameters which can be entered through the keypad, interface parameters are stored in memory even after the circulator is turned off.

#### 12.2. Communication with a PC or a superordinated data system

If the circulator is put into remote control mode via the configuration level, the VFD COMFORT-DISPLAY will read "R -OFF-" = REMOTE STOP. The circulator is now operated via the computer. In general, the computer (master) sends commands to the circulator (slave). The circulator sends data (including error messages) only when the

computer sends a query.



In remote control mode, the start command and all values to be set must be resent by the PC via the interface in case of a power interruption. AUTOSTART is not possible.

A transfer sequence consists of:

- command
- space (⇔; Hex: 20)
- parameter (decimal separation with a period)
- end of file (↓; Hex: 0D)

The commands are divided into **in** and **out** commands. **in** commands: retrieve parameters **out** commands: set parameters

#### Important times for a command transmission:

To ensure a safe data transfer, the time gap between two commands should be at least 250 ms.

The circulator automatically responds to an **in** command with a data string followed by a LF (Line Feed). The next command should only be sent after 10 ms.

The out commands are valid only in remote control mode.

Command to set the working temperature > SETPNT 1< to 55.5 °C  $out\_sp_00 \Leftrightarrow 55.5 \downarrow$ 

Command to retrieve the working temperature > SETPNT 1< in\_sp\_00, J

Response from the circulator:

55.5.⊣





out commands: Setting temperature values or parameters.			values or parameters.
	Command	Parameter	Response of circulator
	out_mode_01	0	Use working temperature > SETPNT 1<
	out_mode_01	1	Use working temperature > SETPNT 2<

out_mode_01	0	Use working temperature > SETPNT 1<
out_mode_01	1	Use working temperature > SETPNT 2<
out_mode_01	2	Use working temperature > SETPNT 3<
out_mode_04	0	Temperature control of internal bath.
out_mode_04	1	External control with Pt100 sensor.
out_mode_05	0	Stop the unit = R –OFF
out_mode_05	1	Start the unit.
out_mode_08	0	Set the control dynamics - aperiodic
out_mode_08	1	Set the control dynamics - standard
out_sp_00	XXX.XX	Set working temperature. "SETPNT 1"
out_sp_01	XXX.XX	Set working temperature. "SETPNT 2"
out_sp_02	XXX.XX	Set working temperature. "SETPNT 3"
out_sp_03	XXX.XX	Set high temperature warning limit "OVERTMP"
out_sp_04	XXX.XX	Set low temperature warning limit "SUBTMP"
out_sp_07	x	Set the pump pressure stage. (1 4)
out_par_04	X.X	CoSpeed 0 5.0 Band limit during external control.
		Setting the maximum difference between the temperatures
		in the internal bath and external system.
out par 06		Xp control parameter of the internal controller. 0.1 99.9
out_par_06	XXX	
out_par_07	XXX	Tn control parameter of the internal controller.3 9999Tv control parameter of the internal controller.0 999
out_par_08	XXX	
out_par_09		
out_par_10	XXX	Proportional portion of the cascade controller. 1 99.9
out_par_11	XXX	The control parameter of the cascade controller. 3 9999
out_par_12	XXX	Tv control parameter of the cascade controller. 0 999
out_par_13	XXX	Maximum internal temperature of the cascade controller.
out_par_14	XXX	Minimum internal temperature of the cascade controller.
out_par_15	XXX	Band limit (upper) 0 200 °C
out_par_16	XXX	Band limit (lower) 0 200 °C

Command	Parameter	Response of circulator
version	none	Number of software version (V X.xx)
status	none	Status message, error message (see page 69)
in_pv_00	none	Actual bath temperature.
in_pv_01	none	Heating power being used (%).
in_pv_02	none	Temperature value registered by the external Pt100 sensor.
in_pv_03	none	Temperature value registered by the safety sensor.
in_pv_04	none	Setpoint temperature of the excess temperature protection
in_sp_00	none	Working temperature "SETPNT 1"
in_sp_01	none	Working temperature "SETPNT 2"
in_sp_02	none	Working temperature "SETPNT 3"
in_sp_03	none	High temperature warning limit "OVERTEMP"
in_sp_04	none	Low temperature warning limit "SUBTEMP"
in_sp_07	none	Pump pressure stage
in_par_01	none	Te - Time constant of the external bath.
in_par_02	none	Si - Internal slope
in_par_03	none	Ti - Time constant of the internal bath.
in_par_04	none	CoSpeed - Band limit (max. difference between the temperatures in the internal bath and external system).
in_par_05	none	Factor pk/ph0:
		Ratio of max. cooling capacity versus max. heating capacity
in_par_06	none	Xp control parameter of the internal controller.
in_par_07	none	Tn control parameter of the internal controller.
in_par_08	none	Tv control parameter of the internal controller.
in_par_09	none	Xp control parameter of the cascade controller.
in_par_10	none	Proportional portion of the cascade controller.
in_par_11	none	Tn control parameter of the cascade controller.
in_par_12	none	Tv control parameter of the cascade controller.
in_par_13	none	Adjusted maximum internal temperature of the cascade controller.
in_par_14	none	Adjusted minimum internal temperature of the cascade controller.
in_par_15	none	Band limit (upper)
 in_par_16	none	Band limit (lower)

Command	Parameter	Response of circulator
in_mode_01	none	Selected setpoint:
		0 = SETPNT 1
		1 = SETPNT 2
		2 = SETPNT 3
		3 = Last setpoint setting was carried out through an external programmer
in_mode_04	none	Internal/external temperature control:
		0 = Temperature control with internal sensor.
		1 = Temperature control with external Pt100 sensor.
in_mode_05	none	Circulator in Stop/Start condition:
		0 = Stop
		1 = Start
in_mode_08	none	Adjusted control dynamics
		0 = aperiodic
		1 = standard

#### 12.4. Status messages

Status messages	Description
00 MANUAL STOP	Circulator in "OFF" state.
01 MANUAL START	Circulator in keypad control mode.
02 REMOTE STOP	Circulator in "r OFF" state.
03 REMOTE START	Circulator in remote control mode.

## 12.5. Error messages

Error messages	Description
-01 LOW LEVEL ALARM	Low liquid level alarm.
-02 REFRIGERATOR ALARM	Control cable of the refrigerated circulator or MVS solenoid valve controller short-circuited or interrupted.
-03 EXCESS TEMPERATURE WARNING	High temperature warning.
-04 LOW TEMPERATURE WARNING	Low temperature warning.
-05 WORKING SENSOR ALARM	Working temperature sensor short-circuited or interrupted.

Remote control

Error messages	Description
-06 SENSOR DIFFERENCE ALARM	Sensor difference alarm. Working temperature and safety sensors report a temperature difference of more than 35 K.
-07 I <sup>2</sup> C-BUS ERROR	Internal error when reading or writing the I <sup>2</sup> C bus.
-08 INVALID COMMAND	Invalid command.
-09 COMMAND NOT ALLOWED IN CURRENT OPERATING MODE	Invalid command in current operating mode.
-10 VALUE TOO SMALL	Entered value too small.
-11 VALUE TOO LARGE	Entered value too large.
-12 TEMPERATURE MEASUREMENT ALARM	Error in A/D converter.
-13 WARNING : VALUE EXCEEDS TEMPERATURE LIMITS	Value lies outside the adjusted range for the high and low temperature warning limits. But value is stored.
-14 EXCESS TEMPERATURE PROTECTOR ALARM	Excess temperature protector alarm
-15 EXTERNAL SENSOR ALARM	External control selected, but external Pt100 sensor not connected.
-20 WARNING: CLEAN CONDENSOR OR CHECK COOLING WATER CIRCUIT OF REFRIGERATOR	Cooling of the condenser is affected. Clean air-cooled condenser. Check the flow rate and cooling water temperature on water- cooled condenser.
-21 WARNING: COMPRESSOR STAGE 1 DOES NOT WORK	Compressor stage 1 does not work.
-22 WARNING: COMPRESSOR STAGE 2 DOES NOT WORK	Compressor stage 2 does not work.
-23 WARNING: HIGH TEMPERATURE ON COMPRESSOR STAGE 1	Excess temperature on compressor stage 1.
-24 WARNING: HIGH TEMPERATURE ON COMPRESSOR STAGE 2	Excess temperature on compressor stage 2.
-25 REFRIGERATOR WARNING	Error in the cooling machine.
-30 CONFIGURATION ERROR: CONFIRM BY PRESSING <ok> ON CIRCULATOR</ok>	The configuration of the circulator does not conform to its present use. Press OK to automatically perform a single modification of the configuration.
-33 SAFETY SENSOR ALARM	Excess temperature sensor short-circuited or interrupted.
-40 NIVEAU LEVEL WARNUNG	Low liquid level warning in the internal reservoir.

#### 13. JULABO Service – Online remote diagnosis

×

 $\nabla$ 

JULABO circulators of the HighTech series are equipped with a black box. This box is implemented in the controller and records all significant data for the last 30 minutes. In case of a failure, this data can be read out from the unit by using special software. This software is available as a **free** download from www.julabo.com \ EasyBlackBox.

- Installation is easy and is performed step by step. Please observe the instructions.
- Data read-out is possible in the conditions "OFF", "R OFF" or "ALARM".
- Connect the circulator to the computer using an interface cable.
- Start the EasyBlackBox program. The program asks for the port used (COM1, .....) and the baud rate of the unit.

You do not have this information on hand? Simply try it out! The program continues to send the request until the correct settings are made.

EasyBlackBox.vi	
Julaba	EasyBlackBox
Einstellungen/Settings	Alarmspeicher/Alarms stored Blackbox
JULABO TopTech Series ME Software Version 1.0 Voltage Supply 230 Volt Barcode: 4294967295 Adjust Offset: 0.00	
Bath: 4 Start Mode: Normal *** Pump Stage *** Pump Stage 1	
*** Serial Interface R5232 Baudrate: 4800 Parity: Even Handshake: Hardware	***
*** TEMPERATURE SETPOI Topical Setpoint/Setpoint3: Setpoint T1: 15.00 C Setp *** TEMPERATURE LIMITS	61.00 C oint T2: 37.00 C Setpoint T3: 61.00 C
Working Temperature Rang Select Temperature Range SavePoti Temperature: 90	ge: -94,90 C to 200.00 C -94,90 C to 200.00 C C it: -99.90 C , High Limit: 105.00 C
Speichern/Save	Hilfe/Help Beenden/Quit

4800 Baud

PortDef.vi

Mit OK bestätigen!

the used baud rate! Confirm with OK!

COM1

 $\nabla$ 

Bitte den verwendeten COM Port und die Übertragungsrate auswählen!

Please choose desired COM port and

OK

 Data is read out and shown on the monitor divided into the sections
 >Einstellungen/Settings<,</li>
 >Alarmspeicher/Alarms stored<,</li>
 >Blackbox<</li>

← see example

- After pressing >Speichern/Save<, a text file is created. The program suggests a filename ->C:\model description and barcode no.<. Modifications are possible.
- E-mail this file to <u>service@julabo.de</u>, JULABO's service department. JULABO is thus able to provide rapid support.

## 14. Cleaning / repairing the unit



#### Caution:

- Always turn off the unit and disconnect the mains cable from the power source before cleaning the unit.
- Prevent humidity from entering into the circulator.
- Electrical connections and any other work must be performed by qualified personnel only.



#### Maintaining the cooling performance

To maintain the full cooling performance, clean the condenser from time to time.

- Switch off the unit, disconnect mains power cable.
- Remove the venting grid.
- Clean the ribbed condenser with a vacuum cleaner.
- Replace the venting grid.
- Switch on the unit.

#### **Cleaning:**

For cleaning the bath tank and the immersed parts of the circulator, use low surface tension water (e.g., soap suds).

Clean the outside of the unit using a wet cloth and low surface tension water.

The circulator is designed for continuous operation under normal conditions. Periodic maintenance is not required.

The tank should be filled only with a bath fluid recommended by JULABO. To avoid contamination, it is essential to change the bath fluid from time to time.

#### Repairs

# Before asking for a service technician or returning a JULABO instrument for repair, please contact an authorized JULABO service station.

When returning the unit:

- Clean the unit in order to avoid any harm to the service personnel.
- Attach a short fault description.
- During transport the unit has to stand upright. Mark the packing correspondingly.
- When returning a unit, take care of careful and adequate packing.
- JULABO is not responsible for damages that might occur from insufficient packing.

JULABO reserves the right to carry out technical modifications with repairs for providing improved performance of a unit.